

Nov 19, 2019

Version 1

# Plasmid Modification by Annealed Oligo Cloning V.1

DOI

[dx.doi.org/10.17504/protocols.io.47tgznn](https://dx.doi.org/10.17504/protocols.io.47tgznn)



Addgene The Nonprofit Plasmid Repository<sup>1</sup>

<sup>1</sup>Addgene



Addgene The Nonprofit Plasmid Repository

Addgene

## Create & collaborate more with a free account

Edit and publish protocols, collaborate in communities, share insights through comments, and track progress with run records.

Create free account

OPEN  ACCESS



DOI: <https://dx.doi.org/10.17504/protocols.io.47tgznn>

External link: <https://www.addgene.org/protocols/annealed-oligo-cloning/>

**Protocol Citation:** Addgene The Nonprofit Plasmid Repository 2019. Plasmid Modification by Annealed Oligo Cloning. protocols.io <https://dx.doi.org/10.17504/protocols.io.47tgznn>

**License:** This is an open access protocol distributed under the terms of the [Creative Commons Attribution License](https://creativecommons.org/licenses/by/4.0/), which permits unrestricted use, distribution, and reproduction in any medium, provided the original author and source are credited



**Protocol status:** Working

**We use this protocol and it's working**

**Created:** July 05, 2019

**Last Modified:** November 19, 2019

**Protocol Integer ID:** 25555

**Keywords:** plasmid modification by annealed oligo cloning, plasmid modification, annealed oligo cloning, oligo cloning, plasmid, cloning, addgene protocol page


## Abstract

This protocol describes plasmid modification by annealed oligo cloning. To see the full abstract and additional resources, please visit the [Addgene protocol page](#).

## Troubleshooting

## Designing overlapping oligos

- 1 Briefly, we will design overlapping oligos that once annealed can be cloned directly into the overhangs generated by restriction digest of existing sites in the original vector.

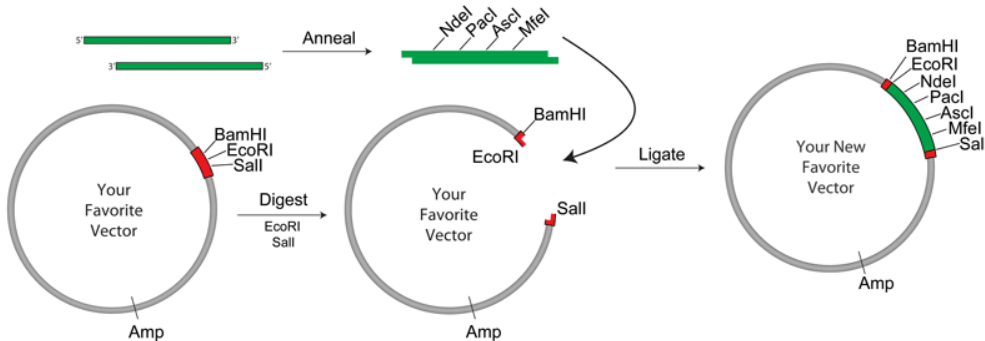

To add NdeI, PaeI, Ascl and MfeI sites between the EcoRI and Sall sites of the vector, we design a top oligo with each of the additional sites in tandem (NdeI - CATATG, PaeI - TTAATTAA, Ascl - GGCGCGCC, MfeI - CAATTG). The bottom oligo will be the reverse complement so that they can anneal.

Top oligo: 5' - CATATGTTAATTAAAGGCGCGCCCAATTG - 3' = 28 bp

Bottom oligo: 3' - GTATACAATTAATTCCGCGCGGGTTAAC - 5' = 28 bp

We also need to include additional bases to complement the overhangs generated when digesting the vector with EcoRI and Sall (see [diagram](#)). To do this, we add 5' - AATTC and G - 3' to the top oligo and 3' - G and CAGCT - 5' to the bottom oligo, making our final oligos 34 bp each:

Top oligo: 5' - AATTCATATGTTAATTAAAGGCGCGCCCAATTGG - 3'

Bottom oligo: 3' - GGTATACAATTAATTCCGCGCGGGTTAACAGCT - 5'

### Note

Note, we could leave off the 3' G on each oligo (and the complementary C of the other oligo), but this would destroy the EcoRI and Sall sites in the final vector.




Order the following oligos from your favorite oligo synthesis company:

Top oligo:	5' - AATTCCATATGTTAATTAAGGCGCGCCCAATTGG - 3'
Bottom oligo:	5' - TCGACCAATTGGGCGCGCCTTAATTAACATATGG - 3'

#### Note

Note, if you plan to phosphatase treat your cut vector it is necessary to use 5'-phosphorylated oligos. This is an option that can be added when ordering them or can be performed enzymatically later.


## Digest and Purify Vector

- 2 While waiting for your oligos to arrive, conduct a restriction digest of  1  $\mu\text{g}$  of vector with EcoRI and Sall
- 3 Run an agarose gel and cut out the band containing your vector DNA
- 4 Gel purify your DNA away from the agarose using a commercially available kit or standard protocol.

## Anneal Oligos

### 5 Anneal oligos

The oligos should be resuspended in annealing buffer (10mM Tris, pH7.5-8.0, 50mM NaCL, 1mM EDTA) and mixed in equimolar concentrations. We recommend mixing

 2  $\mu\text{g}$

each in a total volume of

 50  $\mu\text{L}$

- add additional annealing buffer if necessary to get to 50 $\mu\text{L}$ . Efficient annealing can be achieved by one of two methods:



## 5.1 Method 1

- Place the mixed oligos in a 1.5mL microfuge tube.
- Place tube in  $90^{\circ}\text{C}$  -  $95^{\circ}\text{C}$  hot block and leave for 00:03:00 - 00:05:00 .
- Remove the hot block from the heat source (turn off or move block to bench top) allowing for slow cooling to Room temperature (~ 00:45:00 ).

## 5.2 Method 2

- Place mixed oligos in a PCR tube.
- Place tube in a thermocycler programmed to start at  $95^{\circ}\text{C}$  for 00:02:00 .
- Then, gradually cool to  $25^{\circ}\text{C}$  over 00:45:00 .

## Ligation

- Dilute 5  $\mu\text{L}$  of annealed oligos with 45  $\mu\text{L}$  nuclease-free water and quantify the concentration (should be about 8ng/ $\mu\text{L}$ ).
- Mix the annealed oligos with cut vector in molar ratios (vector:insert) between 4:3 and 1:6 in a standard ligation reaction (ex. to ligate an annealed oligo insert of 50bp in length into a 5kb vector, mix 100ng of the vector with 750  $\mu\text{g}$  - 6 ng of annealed oligos).
- Transform** 2  $\mu\text{L}$  - 3  $\mu\text{L}$  into your favorite competent bacteria and plate.
- Be sure to pick multiple colonies for **mini-prepping** and verify insert by sequencing.

