

Nov 17, 2025

Pole Test Protocol for Motor Function Assessment in Mice

DOI

<https://dx.doi.org/10.17504/protocols.io.5qpvodozdg4o/v1>

Roberto Garcia Swinburn¹, Ernest Arenas¹

¹Karolinska Institute Stockholm

SOX6 mDA differentiation



Roberto Garcia Swinburn

Karolinska Institute Stockholm

Create & collaborate more with a free account

Edit and publish protocols, collaborate in communities, share insights through comments, and track progress with run records.

[Create free account](#)

OPEN  ACCESS



DOI: <https://dx.doi.org/10.17504/protocols.io.5qpvodozdg4o/v1>

Protocol Citation: Roberto Garcia Swinburn, Ernest Arenas 2025. Pole Test Protocol for Motor Function Assessment in Mice. **protocols.io** <https://dx.doi.org/10.17504/protocols.io.5qpvodozdg4o/v1>

License: This is an open access protocol distributed under the terms of the **Creative Commons Attribution License**, which permits unrestricted use, distribution, and reproduction in any medium, provided the original author and source are credited

Protocol status: Working

We use this protocol and it's working

Created: September 09, 2025

Last Modified: November 17, 2025

Protocol Integer ID: 226864

Keywords: pole test protocol for motor function assessment, motor function assessment, general motor skill, mice this protocol, pole test protocol, mice, grafted animal

Abstract

This protocol was used to assess general motor skills and balance on grafted animals.

Guidelines

- Ensure consistent environmental conditions (noise, light, temperature).
- Use the **same pole and orientation** throughout the experiment.
- Observe for abnormal behaviors or jumping, which may invalidate a trial.

Troubleshooting



Training Phase (Minimum 2 Days Prior to Testing)

2d

- 1 Train animals for **at least 2 consecutive days** before the actual test. Each day, perform **3 mock trials per animal** using the same procedure as the test. Ensure mice are acclimated and capable of descending the pole.

2d

Test Day Procedure

- 2 Place the mouse **facing upwards** (head toward the top) on **top of the 20 cm vertical pole** (preferably inside the home cage).
- 3 Begin timing once the mouse rotates and orients downward (i.e., head pointing toward the ground).
- 4
 - Allow the mouse to climb down naturally and stop time when mouse touches the floor. **If the mouse does not descend within 1 minute**, stop the trial and allow the animal to rest. *For incomplete or timed-out trials, repeat after a **minimum rest period of 10 minutes**.*
- 5 Perform a total of **3 successful trials per animal**.

Data Analysis

- 6 Calculate the **mean descent time** (in seconds) across the 3 valid trials for each animal.