

Jan 09, 2026

Postnatal primary hippocampal neuronal cultures

DOI

<https://dx.doi.org/10.17504/protocols.io.kxygx4d4dl8j/v1>

Akio Mori¹, Robert Edwards¹

¹University of California San Francisco



Akio Mori

Create & collaborate more with a free account

Edit and publish protocols, collaborate in communities, share insights through comments, and track progress with run records.

Create free account

OPEN  ACCESS



DOI: <https://dx.doi.org/10.17504/protocols.io.kxygx4d4dl8j/v1>

Protocol Citation: Akio Mori, Robert Edwards 2026. Postnatal primary hippocampal neuronal cultures. **protocols.io**
<https://dx.doi.org/10.17504/protocols.io.kxygx4d4dl8j/v1>

License: This is an open access protocol distributed under the terms of the **[Creative Commons Attribution License](#)**, which permits unrestricted use, distribution, and reproduction in any medium, provided the original author and source are credited

Protocol status: Working

We use this protocol and it's working

Created: November 21, 2025



Last Modified: January 12, 2026

Protocol Integer ID: 233108

Keywords: postnatal primary hippocampal neuronal cultures this protocol, preparation of postnatal primary hippocampal, postnatal primary hippocampal, neuronal culture, preparation

Abstract

This protocol describes the preparation of postnatal primary hippocampal neuronal cultures.

Materials

■ Dissection solution (HBSS++) 500 mL

	A	B
	491 mL	Hank's Balanced Salt Solution (HBSS) (UCSF Cell Culture Facility)
	5 mL	10 mM HEPES (UCSF Cell Culture Facility)
	4 mL	20 mM D-(+)-Glucose solution (G8769, Sigma)

■ Serum Neuronal Media (SNM) 25 mL

	A	B
	22.8 mL	Minimal Essential Medium (MEM) (11090081, Gibco)
	213 µL	45% D-(+)-Glucose solution (G8769, Sigma)
	250 µL	Glutamax (3505006, Gibco)
		(Sterilize with a 0.22 µm syringe filter, and add the following reagents)
	1.25 mL	FBS (SH30070.02, Cytiva)
	500 µL	B27 (17504-044, Gibco)

■ Dissociation

Trypsin 0.25% (15090-046, Gibco)

■ Neurobasal Media (NBM) 25 mL

	A	B
	24.175 mL	Neurobasal Media (21103-049, Gibco)
	325 µL	Glutamax (35050061, Gibco)
		(Sterilize with a 0.22 µm syringe filter, and add the following reagents)
	500 µL	B27 (17504-044, Gibco)

■ Other Supplies

Warner Instruments 1.5 Glass coverslip 25 mm round, 54-0715

Cloning cylinders, 10×10 mm (CLS-1777-03, Chemglass Life Sciences)

Poly-L-lysine for coating (P2636, Sigma-Aldrich)

Plate and coverslip preparation and coating

- 1 Sterilize coverslips and clonal rings by submerging them in 70% (v/v) ethanol and briefly flaming them over a Bunsen burner.
- 2 Place sterile coverslips into each well of a 6-well plate.
- 3 Prepare clonal rings with grease applied to the bottom, and place each ring in the center of a coverslip.
- 4 Add poly-L-lysine (PLL) solution to cover the surface of each well (with coverslip and clonal ring in place). Incubate for 3 hours at 37°C in a 5% CO₂ incubator.
- 5 Remove PLL solution and wash the wells three times with sterile PBS.



Dissection and Dissociation

- 6 Dissect hippocampi from postnatal day 0–1 (P0–P1) mouse pups (appropriate genotypes) in HBSS++ solution.
- 7 Transfer tissue to 4.5 mL of HBSS++ containing 500 µL of 0.25% trypsin (15090-046, Gibco) in a 15 mL tube.
- 8 Incubate tissue in a 37°C water bath for 17 minutes for enzymatic digestion.
- 9 Wash tissue three times with SNM medium to remove trypsin.

Mechanical Dissociation

- 10 Add 1 mL of SNM to the tissue.
- 11 Using a glass Pasteur pipette of original diameter, triturate 3–4 times.



- 12 Flame the pipette to reduce its diameter by half. Triturate the tissue 5 times with the 1/2-diameter pipette.
- 13 Flame the pipette again to reduce its diameter to 1/4 of the original size. Triturate 5 times with the 1/4-diameter pipette.
- 14 Collect the 1 mL suspension of the resulting single cells and transfer to a 1.5 mL tube.

Cell counting and plating

- 15 Take 10 μ L of the cell suspension and transfer to a 0.5 mL tube. Add Trypan Blue solution and count viable cells using a hemocytometer.
- 16 Plate dissociated neurons onto PLL-coated coverslips inside the 6-well plate at a density of 30,000–40,000 viable cells per well.
- 17 Incubate at 37°C, 5% CO₂.
- 18 On the day after plating, carefully replace 75% of the SNM with fresh, pre-warmed Neurobasal Medium (NBM) in each well using a 1 mL pipette.