

Sep 03, 2024

Preferential Lysis of *S. rosetta* for Total RNA (VERSION 2)

 [mSphere](#)

DOI

<https://dx.doi.org/10.17504/protocols.io.261ge4bojv47/v1>

David Booth¹

¹University of California, San Francisco



Vicki Deng

UT Austin

Create & collaborate more with a free account

Edit and publish protocols, collaborate in communities, share insights through comments, and track progress with run records.

Create free account

OPEN  ACCESS



DOI: <https://dx.doi.org/10.17504/protocols.io.261ge4bojv47/v1>

External link: <https://doi.org/10.1128/msphere.00917-24>

Protocol Citation: David Booth 2024. Preferential Lysis of *S. rosetta* for Total RNA (VERSION 2). **protocols.io**
<https://dx.doi.org/10.17504/protocols.io.261ge4bojv47/v1>

Manuscript citation:

Leon F, Espinoza-Esparza JM, Deng V, Coyle MC, Espinoza S, Booth DS Cell differentiation controls iron assimilation in the choanoflagellate *Salpingoeca rosetta*. *mSphere* 10(3). doi: [10.1128/msphere.00917-24](https://doi.org/10.1128/msphere.00917-24)



License: This is an open access protocol distributed under the terms of the **[Creative Commons Attribution License](#)**, which permits unrestricted use, distribution, and reproduction in any medium, provided the original author and source are credited

Protocol status: Working

We use this protocol and it's working

Created: August 19, 2021

Last Modified: September 03, 2024

Protocol Integer ID: 52478

Keywords: choanoflagellate, mRNA, RNA-seq, rna purification, rnas for rna purification, eukaryotic membranes of the choanoflagellate species, membranes with sterol, rna, rnase inhibitor, eukaryotic membrane, protein extraction, protocol suitable for protein extraction, lysis buffer, membrane, different membrane compositions of eukaryote, total rna, purification, choanoflagellate species, based detergent, different membrane composition, sterol, protein, cdna synthesis, preferential lysi, bacteria, protease inhibitor, included protease inhibitor

Abstract

Sterol-based detergents, like digitonin more effectively disrupt membranes with sterols, like those of eukaryotes. This protocol leverages the different membrane compositions of eukaryotes and bacteria to preferentially lyses the eukaryotic membranes of the choanoflagellate species, *S. rosetta*, from a co-culture with bacteria. The lysis buffer includes RNase inhibitors to preserve RNAs for RNA purification or cDNA synthesis, and the included protease inhibitors make this protocol suitable for protein extractions for western blots.

Prepare Lysis Buffer

- 1 Combine the following components for the lysis buffer:

A	B	C	D	E
Chemical	[Final]	[Stock]	Final Vol	Chemical Vol
Water			10 ml	0.41 ml
Tris-HCl, pH 8.0	20 mM	1 M		200 µl
KCl	150 mM	2 M		750 µl
MgCl ₂	5 mM	1 M		50 µl
Sucrose	250 mM	1.75 M (60% w/v)		1ml 420 uL
Cycloheximide	100 ug/ml	100 mg/ml		10 ul
Protease inhibitor tablet	2 mini tablet / 10 ml			1 ml
Digitonin	10 mM	20 mM		5 ml
Sodium Heparin	1 mg/ml	100 mg/ml		100 µl
Pefabloc SC	1 mM	200 mM		50 µl
<i>DTT</i>	<i>1 mM</i>	<i>1 M</i>		<i>10 µl</i>
<i>Turbo DNase</i>	<i>0.1 U/ml</i>	<i>2 U/µl</i>		<i>500 µl</i>
<i>SUPERaseIN</i>	<i>1 U/ml</i>	<i>20 U/µl</i>		<i>500 µl</i>

Notes:

1. Prepare the buffer ahead of time by combining all but the italicized reagents (DTT, Turbo DNase, and SUPERaseIN), splitting into 449.5 µl aliquots, and storing at -20°C.
2. Just before use, thaw the prepared lysis buffer on ice and then add 0.5 µl of 1 M DTT, 25 µl of Turbo DNase, and 25 µl SUPERaseIN to the 449.5 µl aliquot for a total volume of 500 µl.

Count Cells

- 2 Determine the total number of cells in the culture that will be harvested using a hemocytometer.



- 2.1 Fix 200 μ l of cells with 2 μ l of 37% formaldehyde and vortex well.
- 2.2 Pipet up and down to homogenize cells, and pipet 12 μ l of cells into the chamber of a hemocytometer.
- 2.3 Count the number of cells (N) in the four corner quadrants of a Neubauer, bright-line hemocytometer.
- 2.4 Calculate the cell concentration ($[Cells]$ cells/ml) according to this equation:
$$[Cells] = N/4 \bullet 10000 \text{ cells/ml}$$
- 2.5 Determine the volume of the culture (Vol) and then calculate the total number of cells in the culture ($Total$) according to this equation:
$$Total = [cells] \bullet Vol \text{ cells}$$
- 2.6 Calculate the volume of lysis buffer ($LysisBuffer$ μ l) to add to the cells for lysis:
$$LysisBuffer = Total/100000 \mu\text{l}$$

Harvest Cells

- 3 Harvest cells of *S. rosetta*.
- 3.1 Centrifuge the cells in 50 ml conical tubes at  2400 x g, 4°C, 00:05:00
- 3.2 Gently remove supernatant with a serological pipette, leaving a small amount of liquid of the pellet.
- 3.3 Gently remove the remaining supernatant with a fine-tip transfer pipet.

Lyse Cells

- 4 Lyse cells in preferential lysis buffer.



- 4.1 Resuspend the cell pellet ( [go to step #3.3](#)) in the calculated volume of *Lysis Buffer* ( [go to step #2.6](#)).
- 4.2 Pipet the cells gently up and down and then incubate  On ice for  00:10:00 .
- 4.3 Homogenize the lysate by passing 5 times through a 30G needle attached to a luer lock syringe.
- 4.4 Clarify the lysate by centrifugation  6000 x g, 4°C, 00:10:00 10m
- 4.5 Separate the *S. rosetta* lysate from the bacterial pellet by using a gel-loading tip to gently transfer the supernatant into a new tube. Pay attention not to disturb the pellet.

Storage

- 5 Flash freeze lysate in liquid nitrogen and store at -80°C for long term storage.