

Jun 11, 2019

Prep for FACS on blood samples

DOI

<https://dx.doi.org/10.17504/protocols.io.32dgqa6>

Grace Burgin¹, Inbal Benhar¹

¹Broad Institute



Grace Burgin

Create & collaborate more with a free account

Edit and publish protocols, collaborate in communities, share insights through comments, and track progress with run records.

Create free account

OPEN  ACCESS



DOI: <https://dx.doi.org/10.17504/protocols.io.32dgqa6>

Protocol Citation: Grace Burgin, Inbal Benhar 2019. Prep for FACS on blood samples. **protocols.io**

<https://dx.doi.org/10.17504/protocols.io.32dgqa6>

License: This is an open access protocol distributed under the terms of the [Creative Commons Attribution License](#), which permits unrestricted use, distribution, and reproduction in any medium, provided the original author and source are credited

Protocol status: Working

We use this protocol and it's working

Created: June 11, 2019



Last Modified: June 11, 2019

Protocol Integer ID: 24357

Keywords: blood samples for flow, blood sample, prep for fac, fac, sample, prep, blood, flow

Abstract

To prepare blood samples for flow



- 1 Collect blood into EDTA coated tubes
- 2 Transfer 50ul into clean tube - include 1 unstained
***mix from several mice to have 50ul for unstained ctrl tube
- 3 Add 1ml ACK, incubate at RT for 7 min, centrifuge 5' 350g
- 4 Resuspend pellet in FACS buffer + antibody mix
- 5 Incubate in dark for 15' on ice
- 6 Wash (centrifuge) in FACS buffer, 5' 350g
- 7 Resuspend in 300ul FACS Buffer + calcein blue