

Aug 24, 2022

## Preparation and imaging of enriched Golgi from GolgiTAG-IP using Transmission Electron Microscopy

 Proceedings of the National Academy of Sciences of the United States of America

DOI

<https://dx.doi.org/10.17504/protocols.io.x54v9y9nqg3e/v1>

Rotimi Fasimoye<sup>1</sup>, Alan Prescott<sup>2</sup>, Dario R Alessi<sup>1</sup>

<sup>1</sup>Medical Research Council Protein Phosphorylation and Ubiquitylation Unit, School of Life Sciences, University of Dundee, Dow Street, Dundee DD1 5EH, UK;

<sup>2</sup>Dundee Imaging Facility, School of Life Sciences, University of Dundee, Dow Street, Dundee DD1 5EH, UK

asap



Dario R Alessi

Dundee

### Create & collaborate more with a free account

Edit and publish protocols, collaborate in communities, share insights through comments, and track progress with run records.

Create free account

OPEN  ACCESS



**DOI:** <https://dx.doi.org/10.17504/protocols.io.x54v9y9nqg3e/v1>

**External link:** <https://doi.org/10.1073/pnas.2219953120>

**Protocol Citation:** Rotimi Fasimoye, Alan Prescott, Dario R Alessi 2022. Preparation and imaging of enriched Golgi from GolgiTAG-IP using Transmission Electron Microscopy. protocols.io <https://dx.doi.org/10.17504/protocols.io.x54v9y9nqg3e/v1>

**Manuscript citation:**

Fasimoye R, Dong W, Nirujogi RS, Rawat ES, Iguchi M, Nyame K, Phung TK, Bagnoli E, Prescott AR, Alessi DR, Abu-Remaileh M, Golgi-IP, a tool for multimodal analysis of Golgi molecular content. Proceedings of the National Academy of Sciences of the United States of America 120(20). doi: [10.1073/pnas.2219953120](https://doi.org/10.1073/pnas.2219953120)

**License:** This is an open access protocol distributed under the terms of the [Creative Commons Attribution License](#), which permits unrestricted use, distribution, and reproduction in any medium, provided the original author and source are credited

**Protocol status:** Working

**We use this protocol and it's working**

**Created:** August 22, 2022

**Last Modified:** May 31, 2024

**Protocol Integer ID:** 69022

**Keywords:** Enriched Golgi imaging, GolgiTAG-IP, Transmission Electron Microscopy, ASAPCRN, imaging of enriched golgi, cells by golgitag immunoprecipitation, golgitag immunoprecipitation, using transmission electron microscopy transmission electron microscopy, transmission electron microscopy transmission electron microscopy, golgi, organelle, enriched golgi, structure of organelle, using various organelle, various organelle, golgitag, intact cell, cell, fixed intact cell, imaging

**Funders Acknowledgements:**

Aligning Science Across Parkinson's (ASAP) initiative

Grant ID: ASAP-000463

UK Medical Research Council

Grant ID: MC\_UU\_00018/1

## Abstract

Transmission electron microscopy (TEM) is a tool used to image, in good resolution, the structure of organelles within the cell. Available protocols are designed to image structures within fixed intact cells. Here, we described a protocol where Golgi, isolated from cells by GolgiTAG immunoprecipitation (IP) (as described in [ddx.doi.org/10.17504/protocols.io.6qpvrjrogm/v1](https://dx.doi.org/10.17504/protocols.io.6qpvrjrogm/v1)), can be prepared and imaged using TEM. This protocol can also be used to image any organelles isolated using various organelle-IP protocols that are available.

## Attachments



[ivgnbgrdf.docx](#)

804KB



## Materials

### Materials:

-  Paraformaldehyde **Merck MilliporeSigma (Sigma-Aldrich) Catalog #158127**
-  Glutaraldehyde EM Grade 25% **Merck MilliporeSigma (Sigma-Aldrich) Catalog #G5882-50ML**
-  Sodium cacodylate trihydrate **Merck MilliporeSigma (Sigma-Aldrich) Catalog #C0250**
- Osmium tetroxide (Agar. Catalog # R1023)
-  Sodium ferrocyanide decahydrate **Merck MilliporeSigma (Sigma-Aldrich) Catalog #13425**
- Tannic acid (BDH. Catalog # 30337)
- Uranyl acetate (Agar. Catalog # R1260A)
-  Durcupan™ ACM **Merck MilliporeSigma (Sigma-Aldrich) Catalog #44611**
-  Lead(II) citrate tribasic trihydrate **Merck MilliporeSigma (Sigma-Aldrich) Catalog #15326**
-  Sodium citrate monobasic **Merck MilliporeSigma (Sigma-Aldrich) Catalog #71497**
-  Sodium hydroxide (NaOH) **Merck MilliporeSigma (Sigma-Aldrich) Catalog #S5881**
- Potassium chloride (KCl) (VWR. Catalog# 267264)
-  Potassium phosphate monobasic **Merck MilliporeSigma (Sigma-Aldrich) Catalog #P5379**
- Calcium chloride (Calbiochem. Catalog# 208291)
-  (R)-( )-Propylene oxide **Merck MilliporeSigma (Sigma-Aldrich) Catalog #540048**
-  Ethyl alcohol, 200 proof, anhydrous, ≥99.5% **Merck MilliporeSigma (Sigma-Aldrich) Catalog #459836**

### Buffer:

- KPBS (  136 millimolar (mM) KCl,  10 millimolar (mM) KH<sub>2</sub>PO<sub>4</sub>. Adjust to pH 7.25 with KOH).
-  0.4 Mass Percent sodium cacodylate buffer (  17.12 g sodium cacodylate dissolved in water. Adjust to pH7.4 with HCl).
- Reynold lead citrate (  1.33 g lead citrate,  1.76 g sodium citrate;  8 mL 1N NaOH in  50 mL water).

### Equipment:

-  Pierce&trade; Anti-HA Magnetic Beads **Thermo Fisher Catalog #88837**
-  DynaMag&trade;- Spin Magnet **Thermo Fisher Catalog #12320D**
- Microcentrifuge with thermostat (VWR Micro Star 17R. S/N 42209232)
- Scalpel
- P1000 pipette tips
- Pasteur pipette



- Oven (TAAB. I064)
- Leica UCT ultramicrotome (Leica)
-  Clear glass with solid top black phenolic 14B rubber lined cap **VWR International (Avantor) Catalog #215-3571**
- JEOL 1200EX TEM using SIS III camera



## Method

2d 5h 4m

- 1 Perform GolgiTAG-IP as previously described in [dx.doi.org/10.17504/protocols.io.6qpvr djro gmk/v1](https://doi.org/10.17504/protocols.io.6qpvr djro gmk/v1)
- 2 After last wash of beads with KPBS, fix the beads in  1 mL solution containing 4% (w/v) paraformaldehyde and 2.5% (v/v) glutaraldehyde in  0.1 Mass Percent sodium cacodylate buffer dilute in water for  01:00:00 at  Room temperature .  
- 3 Pellet beads with magnet DynaMag™ Spin Magnet (or equivalent) and remove the supernatant.
- 4 Wash beads pellet three times in  1 mL  0.1 Mass Percent sodium cacodylate buffer, pelleting the beads with magnet after each wash. 
- 5 After third wash, spin down beads in tabletop centrifuge at  1000 x g for  00:01:00 . Pellet should be tightly packed.  
- 6 Use a needle to gently dislodge the pellet and use a Pasteur pipette to remove the pellet to a clean glass vial. (Note: All steps are carried out in a glass vial). 
- 7 To ensure rapid dehydration and embedding, cut pellet with scalpel into approximately 1mm<sup>3</sup> pieces, then post-fix in  1 mL 1% (w/v) Osmium tetroxide with 1.5% (w/v) sodium ferrocyanide in  0.1 Mass Percent sodium cacodylate buffer for  01:00:00 at  Room temperature .  
- 8 Wash pellets three times in  0.1 Mass Percent sodium cacodylate buffer, by adding the buffer, allowing the pellets to settle and gently taking out supernatant with P1000 tip. 
- 8.1 Wash pellets three times in  0.1 Mass Percent sodium cacodylate buffer for  00:01:00 (1/3).   
- 8.2 Wash pellets three times in  0.1 Mass Percent sodium cacodylate buffer for  00:01:00 (2/3). 



8.3 Wash pellets three times in [M] 0.1 Mass Percent sodium cacodylate buffer for 00:01:00 (3/3).

1m



9 After the third wash, wash with water three times (as in step 8), then resuspend in 1% w/v tannic acid and 1% w/v uranyl acetate diluted in water for 01:00:00. Remove supernatant.

1h

10 Dehydrate pellet through alcohol series:

50% ethanol for 00:10:00

70% ethanol for 00:10:00

80% ethanol for 00:10:00

90% ethanol for 00:10:00

95% ethanol for 00:10:00

then twice in 100% ethanol for 00:10:00 (1/2)

100% ethanol for 00:10:00 (2/2)

1h 10m



#### Note

Use a P1000 pipette to gently remove ethanol after each series.

11 Further dehydrate in 100% propylene oxide twice for 10 min.

11.1 Further dehydrate in 100% propylene oxide twice for 00:10:00 (1/2).

10m

11.2 Further dehydrate in 100% propylene oxide twice for 00:10:00 (2/2).

10m

12 Resuspend pellet into 50% v/v propylene oxide and 50% v/v Durcupan resin Overnight.

10m





13 Open glass vial to allow propylene oxide to evaporate (This will take approximately  01:00:00 ).

1h

14 To embed pellet in resin, resuspend in 100% Durcupan resin.

15 Embed polymerise resin in an oven at  60 °C for  48:00:00 .

2d

16 Cut sections of polymerised embedded resin with Leica UCT ultramicrotome.

**Note**

Sections should be 70-100nm thick

17 Stain sections with 3% aqueous uranyl acetate for  00:20:00 .

20m


18 Stain with Reynold lead citrate for  00:20:00 .

20m

19 Image stained sections on JEOL 1200EX TEM using SIS III camera.

**Note**

Note: Due to the cutting and many centrifugation steps, it is expected that the magnetic particles dissociate from the plastic bead core.)

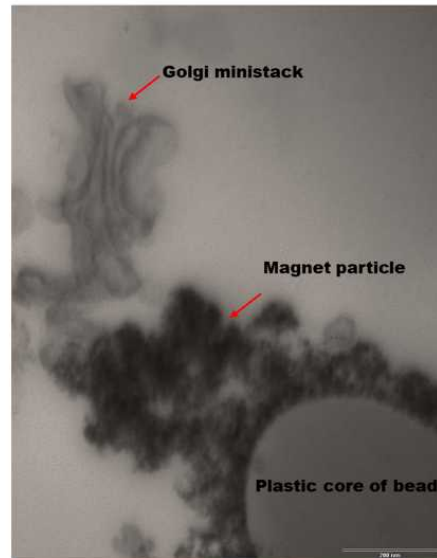

**Figure 1: Transmission electron micrograph of enriched Golgi.** Intact Golgi ministack is shown to be captured by GolgiTAG-IP. Scale bar 200nm