



May 09, 2024

Preparation and imaging of lipid bilayer-coated silica microspheres

 [Nature Methods](#)

DOI

<https://dx.doi.org/10.17504/protocols.io.rm7vzjy2rlx1/v1>

Ezra Bruggeman¹

¹University of Cambridge



Ezra Bruggeman

University of Cambridge

Create & collaborate more with a free account

Edit and publish protocols, collaborate in communities, share insights through comments, and track progress with run records.

Create free account

OPEN  ACCESS



DOI: <https://dx.doi.org/10.17504/protocols.io.rm7vzjy2rlx1/v1>

External link: <https://doi.org/10.1101/2023.02.07.527479>

Protocol Citation: Ezra Bruggeman 2024. Preparation and imaging of lipid bilayer-coated silica microspheres. **protocols.io**
<https://dx.doi.org/10.17504/protocols.io.rm7vzjy2rlx1/v1>

Manuscript citation:

Bruggeman et al., POLCAM: Instant molecular orientation microscopy for the life sciences. bioRxiv 2023.02.07.527479 (Feb 2023), doi: <https://doi.org/10.1101/2023.02.07.527479>



License: This is an open access protocol distributed under the terms of the **Creative Commons Attribution License**, which permits unrestricted use, distribution, and reproduction in any medium, provided the original author and source are credited

Protocol status: Working

We use this protocol and it's working

Created: May 08, 2024

Last Modified: May 31, 2024

Protocol Integer ID: 99426

Keywords: ASAPCRN, imaging of lipid bilayer, lipid bilayer, instant molecular orientation microscopy for the life science, lipid vesicle, microscopy, lipid extrusion, scale glass bead, instant molecular orientation, coated micrometer, 3d localization of dipole, tip probe sonicator

Funders Acknowledgements:

Aligning Science Across Parkinson's (ASAP)

Grant ID: ASAP-000509

Abstract

This protocol describes the preparation and imaging of lipid bilayer-coated micrometer-scale glass beads. This protocol was used to generate the data presented in **Figure 2d-g** in the following publication:

- Bruggeman et al., POLCAM: Instant molecular orientation microscopy for the life sciences. bioRxiv 2023.02.07.527479 (Feb 2023), doi: <https://doi.org/10.1101/2023.02.07.527479>

The described protocol is based on a previously published protocol that uses lipid extrusion instead of a tip probe sonicator to generate lipid vesicles:

- Tingting Wu, Jin Lu and Matthew D. Lew. *pixOL: pixel-wise dipole-spread function engineering for simultaneously measuring the 3D orientation and 3D localization of dipole-like emitters*. bioRxiv (2022). <https://doi.org/10.1101/2021.12.30.474544>

Guidelines

Chloroform is toxic (**SDS**).



Materials

Reagents:

- DPPC in chloroform (**850355C**, Avanti Polar Lipids, sold by Merck). *Note: solid at -20 °C.*
- Cholesterol (**C8667-5G**, Sigma-Aldrich). *Note: will need to be diluted in chloroform, after which it is liquid at -20 °C.*
- Chloroform (**366927**, Sigma-Aldrich). *Note: chloroform is toxic (**SDS**).*
- Tris buffer: 100 mM NaCl, 10 mM **Tris base**, pH 7.4
- Tris-Ca²⁺ buffer: 100 mM NaCl, 3 mM CaCl₂, 10 mM **Tris base**, pH 7.4
- 5 µm diameter silicon dioxide (SiO₂, i.e. silica) microspheres (**44054-5ml-F**, **Sigma-Aldrich**)
- 0.01% poly-L-lysine solution (P4707, Sigma-Aldrich).

Equipment:

- Tip sonicator.
- Plasma cleaner.
- Small centrifuge for Eppendorfs (**Cat. No. 5453000015**, **MiniSpin plus**, **Eppendorf**).
- Heated water bath that can reach 65 °C

Other:

- 0.02 µm syringe filters (6809-1102, Whatman) with compatible syringes.
- Glass vials with PTFE (Teflon) cap liners (**14-955-327**, Thermo Fisher Scientific, or **600460**, Avanti Polar Lipids, sold by Merck). If you don't have vials with PTFE caps, you can use PTFE sealing tape (**Z104388-1PAK**, Merck) before screwing a regular cap on a vial.
- Parafilm (**P7543-1EA**, Sigma-Aldrich)
- Cover glass.
- Frame-seal slide chamber (9×9 mm, SLF0201, Bio-rad).



Preparation of reagents

- 1 Prepare Tris buffer: 100 mM NaCl, 10 mM **Tris base**, pH 7.4
- 2 Prepare Tris-Ca²⁺ buffer: 100 mM NaCl, 3 mM CaCl₂, 10 mM **Tris base**, pH 7.4
- 3 Filter 50 ml of Tris buffer and Tris-Ca²⁺ buffer using a 0.02 µm syringe filter (6809-1102, Whatman).

Preparation of DPPC + 40% cholesterol lipid vesicles

1h 22m

- 4 Dissolve DPPC (850355C, Avanti Polar Lipids) in chloroform (366927, Sigma-Aldrich) to a concentration of **10** 25 mg/mL in a glass vial with a PTFE (Teflon) lined cap (14-955-327, Fisher Scientific). You will need **23** 23 µL per sample.
- 5 Dissolve cholesterol (C8667-5G, Sigma-Aldrich) in chloroform (366927, Sigma-Aldrich) to a concentration of **10** 10 mg/mL in a glass vial with a PTFE (Teflon) lined cap (14-955-327, Fisher Scientific). You will need **20** 20 µL per sample.
- 6 Prepare a DPPC + 40% cholesterol mixture by combining **23** 23 µL **10** 25 mg/mL DPPC with **20** 20 µL **10** 10 mg/mL cholesterol in a new glass vial with a PTFE (Teflon) lined cap (14-955-327, Fisher Scientific).
- 7 Evaporate the solvent **Overnight** under vacuum.
- 8 Re-hydrate the lipids/cholesterol mixture using **1** 1 mL of Tris-Ca²⁺ buffer (**100** 100 millimolar (mM) NaCl, **3** 3 millimolar (mM) CaCl₂, **10** 10 millimolar (mM) Tris base, **pH** 7.4).
- 9 **00:40:00** Vortex for **00:00:30** .

40m 30s



10 Sonicate the solution using a tip sonicator for  00:40:00 (cycles of 45 seconds on and 15 seconds off, 60% amplitude) until the solution runs clear.

40m

11 Centrifuge the solution for  00:01:30 at  14000 rcf to remove residu from the sonicator probe. Keep the supernatant.

1m 30s

Coating silica microspheres with the DPPC/40% cholesterol mixture

35m

12 Dilute the glass beads (5 μ m diameter, 44054-5ML-F, Sigma-Aldrich) to approximately  2.8 mg/mL .

13 Clean the beads by centrifuging and replacing the supernatant with Tris-Ca²⁺ buffer.

14 Heat the glass beads and the lipid vesicle solutions to  65 °C using a heated water bath.

15 Once heated to  65 °C , mix the glass bead and lipid vesicle solution together in a 1:1 ratio. Leave the mixture at  65 °C and wait for  00:30:00 .

30m

16 Turn the heating element of the water bath off, and let the glass bead/lipid vesicle mixture slowly cool down to room temperature inside the water bath. The lipid vesicles will attach to the glass beads, open and form a lipid-bilayer on the glass surface of the beads.

17 Gradually replace the buffer by Tris (100 mM NaCl, 10 mM Tris base, pH 7.4):

17.1 Centrifuge for  00:05:00 at  0.3 rcf .

5m

17.2 Gently replace two thirds of the supernatant with Tris.

17.3 Repeat the last two steps (14.1 and 14.2) a total of 6 times.



- 18 Store the lipid-coated glass beads at  4 °C and use as soon as possible, or within a week of preparation.

Imaging

50m

- 19 Argon plasma clean cover glass (VWR collection, 631-0124) for  00:30:00 in a plasma cleaner (Expanded Plasma Cleaner, PDC-002, Harrick Plasma). 30m
- 20 Create a sample well on the cleaned cover glass by sticking a frame-seal slide chamber (9×9 mm, SLF0201, Bio-rad) on the cover glass.
- 21 Pipet  70 µL of 0.01% PLL (0.01% poly-L-lysine solution, P4707, Sigma-Aldrich) into the well and wait for  00:15:00. The PLL will coat the surface of the cover glass. 15m
- 22 Use a pipet to remove the excess PLL from the well and immediately replace it with  70 µL of filtered PBS.
- 23 Use a pipet to remove the excess filtered PBS from the well and immediately replace with  70 µL filtered PBS. Gently pipet up and down in the corners of the well. Repeat this step 2 more times.
- 24 Use a pipet to remove the excess PBS from the well and immediately replace with  50 µL of the lipid bilayer-coated beads. Wait for  00:05:00. 5m
- 25 Use a pipet to gently remove the excess PBS from the well and immediately replace with  50 µL of a dye of choice, e.g.  1 nanomolar (nM) Nile red for PAINT of the lipid membrane.
- 26 Image the sample the same day.