



Mar 09, 2023

## Preparation of 1L of Foraging Medium

DOI

[dx.doi.org/10.17504/protocols.io.bp2l69n3rlqe/v1](https://dx.doi.org/10.17504/protocols.io.bp2l69n3rlqe/v1)

Alfonso Pérez Escudero<sup>1</sup>, Alid Al-Asmar<sup>1</sup>, gabrielmadirolas<sup>1</sup>

<sup>1</sup>Research Centre on Animal Cognition (CRCA), Centre for Integrative Biology (CBI), Toulouse III University, CNRS, Toulouse, France



APE Lab

CNRS

### Create & collaborate more with a free account

Edit and publish protocols, collaborate in communities, share insights through comments, and track progress with run records.

Create free account

OPEN  ACCESS



DOI: <https://dx.doi.org/10.17504/protocols.io.bp2l69n3rlqe/v1>

**Protocol Citation:** Alfonso Pérez Escudero, Alid Al-Asmar, gabrielmadirolas 2023. Preparation of 1L of Foraging Medium. [protocols.io https://dx.doi.org/10.17504/protocols.io.bp2l69n3rlqe/v1](https://dx.doi.org/10.17504/protocols.io.bp2l69n3rlqe/v1)

**License:** This is an open access protocol distributed under the terms of the [Creative Commons Attribution License](https://creativecommons.org/licenses/by/4.0/), which permits unrestricted use, distribution, and reproduction in any medium, provided the original author and source are credited

**Protocol status:** Working

We use this protocol and it's working



**Created:** September 27, 2022

**Last Modified:** March 09, 2023

**Protocol Integer ID:** 70548

**Keywords:** foraging medium protocol, medium preparation, nematode growth medium, experiments with the worm *Caenorhabditis elegans*, standard nematode growth medium, foraging experiment, making medium, worm *Caenorhabditis elegans*, same basis as the standard nematode growth medium, term worm growth, preparation of 1l, preparation, pouring, bacteriostatic concentration, medium, experiment

**Funders Acknowledgements:**

Human Frontier Science Program

Grant ID: LT000537/2015

CNRS Momentum Program

Fyssen Foundation Research Grant

Gore Family Foundation start-up grant

SEVAB PhD School (Toulouse III University)

## Abstract

Protocol is mainly inspired from: WormBook Methods

[http://www.wormbook.org/chapters/www\\_strainmaintain/strainmaintain.html](http://www.wormbook.org/chapters/www_strainmaintain/strainmaintain.html).

This protocol is for making medium for foraging experiments with the worm *Caenorhabditis elegans*. The medium has the same basis as the standard Nematode Growth Medium (NGM), but contains no peptone (as no long-term worm growth is required for our experiments), and contains antibiotics that prevent our bacterial strains from growing (bacteriostatic concentrations).


It describes both medium preparation and pouring.

## Materials

Agar powder:  Agar **Merck MilliporeSigma (Sigma-Aldrich) Catalog #A1296-10KG**

Cholesterol:  Cholesterol powder **Merck MilliporeSigma (Sigma-Aldrich) Catalog #C3045-5G**

Absolute ethanol:

 Ethanol absolute  $\geq 99.8\%$  AnalaR NORMAPUR® ACS Reag. Ph. Eur. Analytical Reagent **VWR International (Avantor) Catalog #20821.310-1L**

Sodium chloride (NaCl):  Sodium chloride **Merck MilliporeSigma (Sigma-Aldrich) Catalog #S5886-1KG**


Peptone:  Peptone (Bacto) **Becton Dickinson (BD) Catalog #211677-500G**

Chloramphenicol:  Chloramphenicol **Merck MilliporeSigma (Sigma-Aldrich) Catalog #CO378-25G**

Novobiocin:  Novobiocin **Merck MilliporeSigma (Sigma-Aldrich) Catalog #N1628-5G**






### Tissue Culture Dishes:

100×20mm:  Falcon® 100 mm TC-treated Cell Culture Dish **Corning Catalog #353003**





60×15mm:  Petri Dishes 60×15mm with vent **greiner bio-one Catalog #628102**

35×10mm:  Plastic petri dish 35 × 10 mm style **Falcon Catalog #353001**

## Troubleshooting

- 1 In order to perform this protocol, you will need the following solutions:
  - 5 mg/mL cholesterol in absolute ethanol (stored in freezer)
  - pH=6 phosphate buffer (as prepared in: <https://www.protocols.io/view/preparation-of-0-5l-of-phosphate-buffer-ph-6-0-n2bvj8r7bgk5/v1>)
  - 1 M magnesium sulfate (as prepared in: <https://www.protocols.io/view/preparation-of-1m-magnesium-sulfate-solution-mgso4-ca4rsgv6>)
  - 1 M calcium chloride (as prepared in: <https://www.protocols.io/view/preparation-of-1m-calcium-chloride-solution-cacl2-b8parvie>)
  - 10 mg/mL chloramphenicol in absolute ethanol (stored in freezer)
  - 50 mg/mL novobiocin in milliQ water (filter-sterilized and then stored in freezer, in 1.8 mL aliquots to avoid defreezing-refreezing the whole stock)
- 2 Add  16 g  $\pm 0.2$  of agar powder to a 1L clean bottle.
- 3 Add  2.32 g  $\pm 0.02$  of sodium chloride (NaCl).
- 4 Add  780 mL of milliQ water.
- 5 Add a clean magnet (for stirring) and autoclave.
- 6 Put the bottle on a stirrer, and wait until it cools down to around 60°C. At this temperature, the bottle should feel very hot to the touch, but one can hold it for a while without feeling too uncomfortable. The agar should still be liquid.
- 7 While the mix is being stirred, do the following steps in that order, dispensing the reagents right above the surface (to avoid bubbles) with a sterile pipette:
- 8 Add  20 mL of phosphate buffer as prepared in <https://www.protocols.io/view/preparation-of-0-5l-of-phosphate-buffer-ph-6-0-n2bvj8r7bgk5/v1>. This should be added first to prevent any unnecessary precipitation of the next reagents. Wait for ~20 seconds after adding the buffer to give it time to mix well and stabilize the pH.
- 9 Add  0.8 mL of 1M magnesium sulfate (MgSO<sub>4</sub>) solution as prepared in <https://www.protocols.io/view/preparation-of-1m-magnesium-sulfate-solution-mgso4-ca4rsgv6>.



- 10 For this step only, you can put the pipette right below the surface of the agar, otherwise cholesterol tends to stick to the surface.  
Add  0.8 mL of cholesterol (5 mg/mL in absolute ethanol, stored in freezer).
- 11 Add  0.8 mL of 1M calcium chloride ( $\text{CaCl}_2$ ) solution as prepared in <https://www.protocols.io/view/preparation-of-1m-calcium-chloride-solution-cacl2-b8parvie>.  
This step should make the medium a bit cloudier, but remain translucent.
- 12 Add  0.8 mL of chloramphenicol (10 mg/mL in absolute ethanol, stored in freezer).
- 13 Add  1.6 mL of novobiocin (50 mg/mL in milliQ water, stored in freezer).
- 14 Pour in petri dishes before it solidifies, using a pipette and trying to have the same volume in all dishes. We usually pour the plates on the bench, but to minimize the chance of contamination one can pour them in a microbiology hood.  
These are our usual volumes:  
For 100×20mm (diameter x height) petri dishes, we pour 18 mL.  
For 60×15mm petri dishes, we pour 8 mL.  
For 35×10mm petri dishes, we pour 2 mL.
- 15 We store our experimental plates at room temperature for 2 to 8 days. For longer storage and better plate quality, the plates are sometimes stacked in boxes that are sealed with a plastic bag, in order to reduce evaporation.