

Jul 28, 2019

Version 1

Preparation of JF dye for retro-orbital injection in mice V.1

 In 1 collection

DOI

<https://dx.doi.org/10.17504/protocols.io.5u4g6yw>

Boaz Mohar¹

¹HHMI Janelia Research Campus



Boaz Mohar

HHMI Janelia Research Campus

Create & collaborate more with a free account

Edit and publish protocols, collaborate in communities, share insights through comments, and track progress with run records.

Create free account

OPEN  ACCESS



DOI: <https://dx.doi.org/10.17504/protocols.io.5u4g6yw>

Protocol Citation: Boaz Mohar 2019. Preparation of JF dye for retro-orbital injection in mice. **protocols.io**
<https://dx.doi.org/10.17504/protocols.io.5u4g6yw>

License: This is an open access protocol distributed under the terms of the **Creative Commons Attribution License**, which permits unrestricted use, distribution, and reproduction in any medium, provided the original author and source are credited

Protocol status: In development

We are still developing and optimizing this protocol



Created: July 26, 2019

Last Modified: July 28, 2019

Protocol Integer ID: 26236

Keywords: orbital injection in mice, preparation of jf dye, retro orbital injection, jf dye, orbital injection, dye, mice

Abstract

How to prepare **JF dyes** for retro orbital injection.

Materials

STEP MATERIALS

⊗ Dimethyl sulfoxide (DMSO) **Merck MilliporeSigma (Sigma-Aldrich) Catalog #D2650**

⊗ Pluronic™ F-127 (20% Solution in DMSO) **Merck MilliporeSigma (Sigma-Aldrich) Catalog #P3000MP**

⊗ PBS **Invitrogen - Thermo Fisher**

⊗ 29G x 12.7mm 0.3ml syringe **Ulticare Catalog #09230**

Protocol materials

⊗ Dimethyl sulfoxide (DMSO) **Merck MilliporeSigma (Sigma-Aldrich) Catalog #D2650**

⊗ Pluronic™ F-127 (20% Solution in DMSO) **Merck MilliporeSigma (Sigma-Aldrich) Catalog #P3000MP**

⊗ PBS **Invitrogen - Thermo Fisher**

⊗ 29G x 12.7mm 0.3ml syringe **Ulticare Catalog #09230**

Dye preparation

1 Take an aliquote with [M] 100 nmol of dye from  -20 °C storage.

2 Add  20 µL DMSO and do ~10 up and down with the pipeter 



Dimethyl sulfoxide (DMSO) **Merck MilliporeSigma (Sigma-Aldrich) Catalog #D2650**

3 Add 20ul Pluronic F-127 20% in DMSO ~10 up and down with the pipeter 



Pluronic™ F-127 (20% Solution in DMSO) **Merck MilliporeSigma (Sigma-Aldrich) Catalog #P3000MP**

4 Add 200ul 1x PBS ~10 up and down with the pipeter 



PBS **Invitrogen - Thermo Fisher**

5 Drew 200ul without bubbles to a 29G syringe



29G x 12.7mm 0.3ml syringe **Ulticare Catalog #09230**

6 Follow the Retro-orbital injection protocol



Protocol

NAME

Retro-orbital injection of virus or dye in mice

CREATED BY

Boaz Mohar

Preview

- 6.1 Prepare either virus dilution (5.0×10^{11} – 1.0×10^{12}) with PBS prepare or a **dye aliquot**. Injection is up to 200ul per mouse.
- 6.2 Take one mouse from the cage into an induction chamber and start isoflurane at 3% with flow rate of ~1L/min
- 6.3 After the mouse has slow breathing (~1/s) move the mouse to a nose cone delivery of isoflurane at 1.5% with around the same flow rate.
- 6.4 Expose the eye ball with two fingers (above and below) and make sure the you are not pressing on the trachea and the mouse's breathing is regular. 
- 6.5 Use a 29G syringe at a 30-degree angle beveled downwards until you hit the bone, retract a bit and slowly inject. See cited paper for more details. 

Citation

Tal Yardeni, Michael Eckhaus, H. Douglas Morris, Marjan Huizing, Shelley Hoogstraten-Miller

(Invalid date). Retro-orbital injections in mice. LabAnimal.

[10.1038/labani0511-155](https://doi.org/10.1038/labani0511-155)

LINK



29G x 12.7mm 0.3ml syringe **Ulticare Catalog #09230**



- 6.6 Remove the needle slowly and press the eyelid shut for a few seconds.
- 6.7 Validating the injection was successful is possible with JF525 by looking at the pee of the animal after it wakes up.
Leaving the animal in a clean cup with kim-wips will collect the pee and it would seem pink.

