



Mar 19, 2025

Preparation of Negative Stain Transmission Electron Microscopy Grids

DOI

<https://dx.doi.org/10.17504/protocols.io.4r3l29x1jv1y/v1>

Patricia Yuste-Checa¹, Chenchen Mi¹, F Ulrich Hartl¹

¹Department of Cellular Biochemistry, Max Planck Institute of Biochemistry, Martinsried, Germany



Patricia Yuste-Checa

MPI Biochemistry

Create & collaborate more with a free account

Edit and publish protocols, collaborate in communities, share insights through comments, and track progress with run records.

Create free account

OPEN  ACCESS



DOI: <https://dx.doi.org/10.17504/protocols.io.4r3l29x1jv1y/v1>

Protocol Citation: Patricia Yuste-Checa, Chenchen Mi, F Ulrich Hartl 2025. Preparation of Negative Stain Transmission Electron Microscopy Grids. **protocols.io** <https://dx.doi.org/10.17504/protocols.io.4r3l29x1jv1y/v1>

License: This is an open access protocol distributed under the terms of the **Creative Commons Attribution License**, which permits unrestricted use, distribution, and reproduction in any medium, provided the original author and source are credited

Protocol status: Working

We use this protocol and it's working

Created: March 10, 2025



Last Modified: March 19, 2025

Protocol Integer ID: 124243

Keywords: ASAPCRN, preparation of negative stain transmission electron microscopy grid, negative stain protein samples for transmission electron microscopy, negative stain transmission electron microscopy grid, negative stain protein sample, transmission electron microscopy, imaging, protein sample

Funders Acknowledgements:

Aligning Science Across Parkinson's

Grant ID: ASAP-000282

Abstract

This protocol details how to prepare negative stain protein samples for transmission electron microscopy (TEM) imaging.



Preparation of Negative Stain

3m 30s

1 Using an EM-grade forceps, place carbon film supported copper grids (Quantifoil) with the matte carbon side (other side, polished, shiny copper) facing up onto a filter paper inside a glass petri dish.

2 Glow-discharge the carbon grids using a plasma cleaner like PDC-3XG (Harrick) for  00:00:30 .

30s

3 Hold the grid with EM-grade forceps with the carbon side facing up above a filter paper.

4 **Blotting:** Pipette  4 μL of your sample onto the carbon grid and incubate for  00:01:00 .

1m



5 Wick away liquid with filter paper wedge.

6 **Stain:** Add  7 μL of 2% uranyl acetate solution (Electron Microscopy Sciences) on top of the carbon grid and incubate for  00:01:00 .

1m



Note

- Uranyl acetate is toxic and radioactive! Wear appropriate safety gear, prevent incorporation. Waste containing uranyl acetate must be discarded as radioactive waste.
- Before use, centrifuge the 2% uranyl acetate solution at  20000 x g,  00:10:00 in order to pellet possible precipitates.

7 Carefully dry the carbon grid using a filter paper.

8 Add  7 μL of 2% uranyl acetate solution (Electron Microscopy Sciences) on top of the carbon grid and incubate for  00:01:00 .

1m





- 9 Carefully dry completely the carbon grid using a filter paper wedge.
- 10 Air dry the carbon grid for a few minutes.
- 11 Store the carbon grid in a cassette until imaging by TEM.