

Aug 05, 2024

🌐 Preparation of primary hippocampal neurons

📖 [Science Advances](#)

DOI

<https://dx.doi.org/10.17504/protocols.io.j8nlkk241l5r/v1>

Arpine Sokratian¹, yuan.yuan², andrew.west west²


¹Duke Univeristy; ²Duke University

West lab protocols



Arpine Sokratian

Duke Univeristy

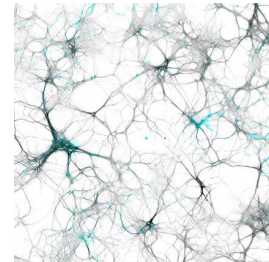

Create & collaborate more with a free account

Edit and publish protocols, collaborate in communities, share insights through comments, and track progress with run records.

Create free account

OPEN  ACCESS



DOI: <https://dx.doi.org/10.17504/protocols.io.j8nlkk241l5r/v1>

Protocol Citation: Arpine Sokratian, yuan.yuan , andrew.west west 2024. Preparation of primary hippocampal neurons. protocols.io <https://dx.doi.org/10.17504/protocols.io.j8nlkk241l5r/v1>

Manuscript citation:

DOI: [10.1126/sciadv.adi8716](https://doi.org/10.1126/sciadv.adi8716)

License: This is an open access protocol distributed under the terms of the **Creative Commons Attribution License**, which permits unrestricted use, distribution, and reproduction in any medium, provided the original author and source are credited

Protocol status: Working

We use this protocol and it's working

Created: May 30, 2022

Last Modified: August 05, 2024

Protocol Integer ID: 63449

Keywords: primary neuron, mouse hippocampus, ASAPCRN, preparation of primary hippocampal neuron, primary hippocampal neuron culture, primary hippocampal neuron, rodent embryos p1 from ntg, rodent embryos p1, transgenic mice, preparation, cell

Funders Acknowledgements:

ASAP

Grant ID: 020527

Abstract

This protocol details preparation of the primary hippocampal neuron culture. The protocol involves extraction of dissecting hippocampi from rodent embryos P1 from nTg or transgenic mice, enzymatic digestion to dissociate cells, and seeding onto poly-D-lysine-coated dishes.

Attachments



[451-951.docx](#)

23KB



Materials

Prepare solutions for plating

Borate buffer:

- [M] 50 millimolar (mM) ,  3.09 g/L ,  8.5
- Sterilize with 0.2 um filter.

Poly D lysine (PDL) hydrobromide (gibco, 0.1mg/ml, REF: A3890401. Life Technologies Corporation 3175 Staley Rd., Grand Island, NY 14072, USA):

	A	B
	40x stock	2 mg/mL
	Borate buffer pH 8.5	100 mg/50 mL

1. For 1x, dilute 40x with borate buffer.
2. Sterilize with 0.2 um filter.
3. Store in  -20 °C .

Prepare solutions for dissection

Note

The same filter unit can be applied for each step of dissection

HBSS (Hanks' Balanced Salt Solution 1X)  HBSS Gibco - Thermo Fisher Scientific Catalog #14025-092

	A	B
	HBSS	490 mL
	HEPES pH 7.4	6 mL
	pen/strep	3 mL
	100 mM pyruvic acid	6 mL
	distilled H2O	100 mL
	Total	600 mL

Filter (VWR Vacuum Filtration 250 ML 0.45 µm PES FILTER UNIT, Made in China Manufactured for VWR International, LLC 100 Matsonford Rd, Radnor, PA 19087)

Enzyme solution

	A	B
	HBSS	10 mL
	Papain suspension	200 µL
	L-cysteine	2 mg
	0.5 mM EDTA	22 µL

1. Rotate for  00:30:00 .
2. Filter (Millex-GP. Syringe-driven Filter Unit. 33mm, Pes Membrane, 0.22 µm, Sterilized. Merck Millipore Ltd. Tullagreen, Carrigtwohill, Co. Cork, IRELAND. Rev. 07/20).

Plating media (containing neurobasal media, supplemented with 5% FBS, 1x B27 supplement (ThermoFisher,17504044),  0.5 millimolar (mM) L-glutamine, and 100 unit per mL of Penicillin-Streptomycin)

	A	B
	Neurobasal	183 mL
	FBS	10 mL
	GlutaMAX	2 mL
	B27	4 mL
	Penicillin-Streptomycin	1 mL

1. Filter (VWR Vacuum Filtration 250ML 0.45 µm PES FILTER UNIT, Made in China Manufactured for VWR International, LLC 100 Matsonford Rd. Radnor, PA 19087) and keep in  4 °C .
2. Warm  Overnight /2 hours before in incubator in T75 flask (Thermo Fisher Scientific, Nunc™ EasYFlask™ 25 cm² Nuncion™ Delta Surface.Nunc A/S. Kamstrupvej 90. P.O. Box 280 DK-4000 Roskilde. Denmark) to calibrate pH.

Neuronal/culture media (containing neurobasal media supplemented with B27 and  0.5 millimolar (mM) L-glutamine)

	A	B
	Neurobasal	193 mL
	GlutaMAX	2 mL
	B27	4 mL

1. Filter (VWR Vacuum Filtration 250ML 0.45 µm PES FILTER UNIT, Made in China Manufactured for VWR International, LLC 100 Matsonford Rd. Radnor, PA 19087).
2. Warm  Overnight /2 hours before in incubator in T75 flask to calibrate pH.

Materials

-  Poly-D-Lysine **Thermo Fisher Scientific Catalog #A3890401**
-  HBSS **Gibco - Thermo Fisher Scientific Catalog #14025-092**
-  Papain **Worthington Biochemical Corporation Catalog #LS003126**
-  B-27 Supplement **Gibco - Thermo Fisher Scientific Catalog #17504044**
-  Neurobasal®; Medium, minus phenol red **Thermo Fisher Catalog #12348017**
-  Glutamax (100x) **Gibco - Thermo Fisher Scientific Catalog #35050-061**

Protocol materials

-  Neurobasal®; Medium, minus phenol red **Thermo Fisher Catalog #12348017**
-  Glutamax (100x) **Gibco - Thermo Fisher Scientific Catalog #35050-061**
-  Poly-D-Lysine **Thermo Fisher Scientific Catalog #A3890401**
-  HBSS **Gibco - Thermo Fisher Scientific Catalog #14025-092**
-  Papain **Worthington Biochemical Corporation Catalog #LS003126**
-  B-27 Supplement **Gibco - Thermo Fisher Scientific Catalog #17504044**
-  Nunc™ Cell-Culture Treated Multidishes 48 well **Thermo Scientific Catalog #12565322**
-  Water sterile-filtered **Merck MilliporeSigma (Sigma-Aldrich) Catalog #RNBK1827**



Plate preparation

1 Wash 48 well plates



Nunc™ Cell-Culture Treated Multidishes 48 well Thermo Scientific Catalog #12565322

with distilled H₂O



Water sterile-filtered Merck MilliporeSigma (Sigma-Aldrich) Catalog #RNBK1827

1.1 Wash 48 well plates with  0.5 mL distilled H₂O. (1/3)

1.2 Wash 48 well plates with  0.5 mL distilled H₂O. (2/3)

1.3 Wash 48 well plates with  0.5 mL distilled H₂O. (3/3)

2 Treat each well with  0.5 mL PDL at  37 °C for  01:00:00 .

1h

3 Wash with distilled H₂O.



3.1 Wash with  0.5 mL distilled H₂O. (1/3)

3.2 Wash with  0.5 mL distilled H₂O. (2/3)

3.3 Wash with  0.5 mL distilled H₂O. (3/3)

4 Place  0.5 mL plating media into each well at  37 °C .

Note

Note: This should be done right before dissection.

Dissection

- 5 Collect the hippocampi of 1-day postnatal CD1 mice pups.
- 6 Clean surgical instrument with 70% ethanol (Sharp fine scissors 14060-11 F.S.T, Sekmen Forceps, 11008-15 F.S.T).
- 7 Cut off head and using scissors, cut skin of the top of the head laterally.
- 8 Stabilizing head with tweezers, use another set of tweezers to pull skin to the side.
- 9 Cut through the bone.
- 10 Using a spoon, scoop brain from underneath and place into culture dish.
- 11 Separate the two hemispheres at the interhemispheric fissure from the brainstem.
- 12 Cut off the olfactory bulbs.
- 13 Carefully pull off the meninges but leave connected at the bottom.
- 14 Flip so that the bottom of the brain is facing up.
- 15 Completely pull off the meninges and associated tissues.
- 16 Hippocampus should be visible and make two cuts at either end.



- 17 Flip out hippocampus.
- 18 Trim hippocampus so that additional tissue is discarded.
- 19 Dissect hippocampus and using one pair of tweezers to push the tissue onto another, place hippocampus into the 15-mL tube with the  10 mL L-glutamin.

Cell culture

12h 45m

- 20 Wash with HBSS by pipetting to aspirate and NOT vacuuming.  
- 20.1 Wash with  10 mL HBSS by pipetting to aspirate and NOT vacuuming. (1/2)
- 20.2 Wash with  10 mL HBSS by pipetting to aspirate and NOT vacuuming. (2/2)
- 21 Leave  1 mL HBSS in tube.
- 22 After filtering the enzyme solution with rotator, add all  10 mL of solution to the tube. 
- 23 Incubate at  37 °C for around  00:45:00 and no longer than 1 hour.  45m
- 24 Remove media and leave  1 mL .
- 25 Pipette  10 mL plating media and  50 µL DNase (DNase stock is  50 µg/mL) 
into a 15-mL tube and invert 2-3 times.



- 26 Pipette  10 mL DNase solution to the  1 mL digestion solution with hippocampi and invert 2-3 times. 
- 27 Remove all media.
- 28 Wash with  10 mL plating media. 
- 29 Wash with HBSS by pipetting to aspirate and NOT vacuuming. 
- 29.1 Wash with  10 mL HBSS by pipetting to aspirate and NOT vacuuming. (1/2)
- 29.2 Wash with  10 mL HBSS by pipetting to aspirate and NOT vacuuming. (2/2)
- 30 Remove HBSS and leave  1 mL with HBSS with hippocampi.
- 31 Resuspend hippocampi with P1000 filter tip and pipet up and down 20 times until obvious chunks disappear. 
- 32 Add  5 mL plating media. 
- 33 Pass cells through strainer with 40 μm mesh size.
- 34 Count cells and plate 50,000 cells per well for a 48 well plate, and 25,000 cells per well for a 96 well plate.
- 34.1 Using cell counter.
1. Place glass slip on top of slide and pipette  10 μL cell solution under slip.
 2. Count number of cells in one 4 $\times$ 4 grid (live cells appear round with dark ring and transparent inside) $\rightarrow$ # $\times 10^4$ cells/mL.



- 34.2 Make sure not to add a small volume of cells to a large volume of plating media, aim for  250 μL of cell solution per well for a 48 well plate.
- 35 Leave plates in  37 °C incubator for  12:00:00 .
- 36 Remove the plating media and swap with culture media containing neurobasal media supplemented with B27 and  0.5 millimolar (mM) L-glutamine.
- 37 Culture the cells for an additional 7 days before use. At day-in-vitro (DIV) 10, add fibril strains to each well of neurons to a final estimated concentration of  0.62 nanomolar (nM) ( 1 $\mu\text{g}/\text{mL}$).

12h

