

Jun 03, 2019

Version 1

Prepare 2L of B-Broth for culturing bacteria V.1

DOI

<https://dx.doi.org/10.17504/protocols.io.3nmgmc6>

Cancer Research UK / Wellcome Gurdon Institute media kitchen¹

¹Cancer Research UK / Wellcome Gurdon Institute



Cristian Riccio

University of Cambridge

Create & collaborate more with a free account

Edit and publish protocols, collaborate in communities, share insights through comments, and track progress with run records.

Create free account

OPEN  ACCESS



DOI: <https://dx.doi.org/10.17504/protocols.io.3nmgmc6>

Protocol Citation: Cancer Research UK / Wellcome Gurdon Institute media kitchen 2019. Prepare 2L of B-Broth for culturing bacteria. [protocols.io https://dx.doi.org/10.17504/protocols.io.3nmgmc6](https://dx.doi.org/10.17504/protocols.io.3nmgmc6)

License: This is an open access protocol distributed under the terms of the [Creative Commons Attribution License](#), which permits unrestricted use, distribution, and reproduction in any medium, provided the original author and source are credited

Protocol status: Working

We use this protocol and it's working

Created: June 03, 2019

Last Modified: June 03, 2019

Protocol Integer ID: 23981

Keywords: culturing bacteria, broth, bacteria, 2l

Abstract

Prepare 2L of B-Broth for culturing bacteria

Materials

MATERIALS

 double distilled water (ddH₂O)

 Sodium chloride meets analytical specification of Ph.Eur Fisher Chemical **Fisher Scientific Catalog #S/3160/65**


 bacto tryptone

Safety warnings

 Make sure your know how to use the autoclave.



1

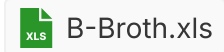

2

	Ingredients	Quantity	
	Bacto tryptone	20g	
	NaCl		10g
	Double distilled H2O	up to 2L	

3

1	Dissolve Bacto tryptone and NaCl in 1L double distilled H2O in a 2L measuring cylinder
2	Make up to final volume using double distilled H2O
3	Aliquot in 100 ml bottles, label and autoclave