

Jun 03, 2019

Version 1

Prepare NGM plates for nematode, with peptone, without fungizone V.1

DOI

<https://dx.doi.org/10.17504/protocols.io.3ntgmen>

Cancer Research UK / Wellcome Gurdon Institute media kitchen¹

¹Cancer Research UK / Wellcome Gurdon Institute



Cristian Riccio

University of Cambridge

Create & collaborate more with a free account

Edit and publish protocols, collaborate in communities, share insights through comments, and track progress with run records.

Create free account

OPEN  ACCESS



DOI: <https://dx.doi.org/10.17504/protocols.io.3ntgmen>

Protocol Citation: Cancer Research UK / Wellcome Gurdon Institute media kitchen 2019. Prepare NGM plates for nematode, with peptone, without fungizone. **protocols.io** <https://dx.doi.org/10.17504/protocols.io.3ntgmen>

License: This is an open access protocol distributed under the terms of the **Creative Commons Attribution License**, which permits unrestricted use, distribution, and reproduction in any medium, provided the original author and source are credited

Protocol status: Working

We use this protocol and it's working

Created: June 03, 2019



Last Modified: June 03, 2019

Protocol Integer ID: 23987

Keywords: fungizone prepare ngm plates for nematode, fungizone prepare ngm plate, ngm plates for nematode, nematode, fungizone, ngm plate, peptone

Abstract

Prepare NGM plates for nematodes, with peptone, without fungizone

Materials

MATERIALS



Agar



Difco Bacto peptone **Fisher Scientific**



double distilled water (ddH₂O)



Sodium chloride meets analytical specification of Ph.Eur Fisher Chemical **Fisher Scientific Catalog #S/3160/65**



SYCHEM autoclave **Syschem**

Safety warnings

 Make sure you know how to use the autoclave before starting this protocol.



1

 NGM_agar_1L.xls

2

	Ingredients		Quantity
	NaCl		3g
	bacto peptone		2.5g
	Agar		17g
	Double distilled H2O	972ml	

3

	Measure 972ml double distilled H2O and put in a 1L duran bottle with a magnetic flea
	Add 3g NaCl, 2.5g bacto peptone and stir.
	Stir until all solutes are dissolved then add the agar.
	Leave magnetic flea in bottle.
	Label, date and autoclave