

Jul 13, 2022

## Primary microglial culture

DOI

[dx.doi.org/10.17504/protocols.io.b4ubqwsn](https://dx.doi.org/10.17504/protocols.io.b4ubqwsn)

Xiqun Chen<sup>1</sup>, Qing Ye<sup>2</sup>

<sup>1</sup>Department of Neurology Massachusetts General Hospital Harvard Medical School Charlestown MA 02129 USA, Aligning Science Across Parkinson's (ASAP) Collaborative Research Network Chevy Chase MD 20815 USA;

<sup>2</sup>Department of Neurology Massachusetts General Hospital Harvard Medical School Charlestown MA 02129 USA, Department of Neurology Longhua Hospital Shanghai University of Traditional Chinese Medicine Shanghai 200032 China



Pranay Srivastava

### Create & collaborate more with a free account

Edit and publish protocols, collaborate in communities, share insights through comments, and track progress with run records.

Create free account

OPEN  ACCESS



DOI: <https://dx.doi.org/10.17504/protocols.io.b4ubqwsn>

**Protocol Citation:** Xiqun Chen, Qing Ye 2022. Primary microglial culture. **protocols.io**  
<https://dx.doi.org/10.17504/protocols.io.b4ubqwsn>

**License:** This is an open access protocol distributed under the terms of the **Creative Commons Attribution License**, which permits unrestricted use, distribution, and reproduction in any medium, provided the original author and source are credited



**Protocol status:** Working

**We use this protocol and it's working**

**Created:** February 09, 2022

**Last Modified:** May 31, 2024

**Protocol Integer ID:** 57955

**Keywords:** ASAPCRN, primary microglial culture mixed glial cell, primary microglial culture, mixed glial cell culture, microglial cell, mixed glial cell, mechanical shaking of the mixed glial cell culture, glial cell

## Abstract

Mixed glial cells were obtained from C57BL/6 mice embryonic day 17. Cells were cultured in high-glucose DMEM/F12 supplemented with 10% FBS in humidified air containing 5% CO<sub>2</sub> at 37°C. The culture medium was replaced with fresh medium 24 h after the initial preparation and every 3 days thereafter. After 1 week, microglial cells were obtained by mechanical shaking of the mixed glial cell cultures for 1 h. Cells were routinely monitored for purity by ionized calcium-binding adaptor molecule 1 (Iba1) staining and the population of Iba1<sup>+</sup> cells was >95%.

## Protocol materials

⊗ DMEM, high glucose Thermo Fisher Scientific Catalog #11965092

⊗ DMEM, high glucose Thermo Fisher Scientific Catalog #11965092

⊗ Fetal Bovine Serum (FBS) ATCC Catalog #30-2020

⊗ DMEM, high glucose Thermo Fisher Scientific Catalog #11965092

⊗ DMEM, high glucose Thermo Fisher Scientific Catalog #11965092

## Troubleshooting



- 1 Primary microglial culture - Use C57BL/6J mice at embryonic day 17
- 2 Anesthetized pregnant mice (1% sodium pentobarbital, 80mg/kg), dissect their embryos and collect the cortex.  
(Separate and remove the soft membrane and blood vessels, rinse the cerebral cortex in PBS, and use the ophthalmic scissor to cut pieces of the cortex)
- 3 Collect the cortices in PBS in a 50 ml tube on ice  
(The 50 ml tube contains 30 ml of PBS) On ice
- 4 Transfer the cortices to 15 ml tubes containing 1.5 ml trypsin-EDTA (0.25%) and incubate it at 37 °C for 00:15:00 Dissociate the cortices by triturating with a 10 mL serological pipette 10 – 15 times 15m
- 5 Centrifuge the dissociated cortices ( 400 x g , 00:05:00 ). Aspirate the media and resuspend the pellet in 5 ml 5m  
 DMEM, high glucose **Thermo Fisher Scientific Catalog #11965092**
- 6 Count the cells and plate them in a density of 50,000 cells/cm<sup>2</sup> into PLL coated T-75 flask. Makeup the volume to 15 ml with the  
 DMEM, high glucose **Thermo Fisher Scientific Catalog #11965092**
- 7 Change the culture medium the next day followed by the addition of a fresh culture medium every 168:00:00 till 14 days of culture 1w
- 8 Put the flask on a shaker for 06:00:00 at 37 °C (The microglia grow as a monolayer on the top) 8h 5m  
Collect the detached cells, centrifuge ( 400 x g, Room temperature , 00:05:00 )  
and resuspended in 4 mL of  
 DMEM, high glucose **Thermo Fisher Scientific Catalog #11965092**

Count the cells and plate at a density of 50,000 cells/cm<sup>2</sup> in a 24 well plate  
Incubate them at 37 °C in 5% CO<sub>2</sub>. After 02:00:00 microglia usually attaches to the bottom of the plate.


The purity of the cells was monitored by staining with ionized calcium-binding adaptor molecule 1 (Iba1) antibody






(Microglia obtained by this method have a purity of 90%-95% and can be identified by immunofluorescence staining.

## 9 Transduction with BRAF (Optional)

1d

The purified microglial cells were seeded at a density of 80% and transduced with BRAFV<sup>600E</sup>, or BRAF<sup>WT</sup>, or vector lentivirus with 8 µg/ml polybrene (Sigma–Aldrich, USA) for  24:00:00 .

10 After transduction, the cells were cultured in  DMEM, high glucose **Thermo Fisher Scientific Catalog #11965092** with or without  Fetal Bovine Serum (FBS) **ATCC Catalog #30-2020** for  120:00:00 , and the medium, as well as the cells, were collected for subsequent experiments.

5d