

Mar 03, 2025

Primary Neuron Cultures

DOI

<https://dx.doi.org/10.17504/protocols.io.3byl4wpyovo5/v1>

Emily Tabaie¹

¹University of California, Riverside



Emily Tabaie

University of California, Irvine

Create & collaborate more with a free account

Edit and publish protocols, collaborate in communities, share insights through comments, and track progress with run records.

Create free account

OPEN  ACCESS



DOI: <https://dx.doi.org/10.17504/protocols.io.3byl4wpyovo5/v1>

Protocol Citation: Emily Tabaie 2025. Primary Neuron Cultures. protocols.io
<https://dx.doi.org/10.17504/protocols.io.3byl4wpyovo5/v1>

License: This is an open access protocol distributed under the terms of the **Creative Commons Attribution License**, which permits unrestricted use, distribution, and reproduction in any medium, provided the original author and source are credited

Protocol status: Working

We use this protocol and it's working

Created: March 03, 2025



Last Modified: March 03, 2025

Protocol Integer ID: 123717

Keywords: preparation of primary cortical neuron culture, primary cortical neuron culture, primary neuron cultures this protocol, primary neuron culture, embryo, mice

Abstract

This protocol is for the isolation and preparation of primary cortical neuron cultures from the embryos of C57BL/6 mice.

Primary Neuron Cultures

- 1 Set up a breeder pair 18 days prior to harvest
- 2 When the female is close to giving birth coat a 6 well plate days before the dissection occurs and make the media
- 3 Prior to coating, sterilize coverslips by autoclaving
- 4 Using sterile forceps, transfer one pre-sterilized glass coverslip per well of a 6 well plate
- 5 Dispense 1-2 mL of Poly-L-Lysine (PLL) solution into each well containing a coverslip
- 6 Incubate for a minimum of 2 hours at 37°C or overnight at 4°C
- 7 Wash each well with 1 mL Tissue Grade Water two times
- 8 Add 2 mL of complete neurobasal media to each well of a 6 well plate
- 9 Prepare Sacrifice pregnant female by isoflurane/cervical dislocation
- 10 Surgically open peritoneum and remove embryonic sacks
- 11 Prepare 1950 µL of Enzyme Mix 1 for up to 400 mg of tissue (5 brains)
- 12 Pre-heat the mixture at 37°C for 10-15 min prior to use



- 13 Carefully open each sack and separate the embryo in a petri dish filled with cold HBSS without Ca^{2+} or Mg^{2+}
- 14 Remove the head and the skullcap to excise the brain and get rid of any red blood cells
- 15 Separate the forebrain from the cerebellum and place the forebrain into a 15 mL conical tube containing 14 mL of HBSS (without Ca^{2+} or Mg^{2+})
- 16 Centrifuge at 300 x g for 2 min at room temperature and aspirate the supernatant
- 17 Add 1950 μL of enzyme mix 1
- 18 Incubate in a closed tube for 15 min at 37°C under slow, continuous rocking/rotation
- 19 Prepare 30 μL of enzyme mix 2 per eppendorf tube and add to the 15 mL conical and invert the tubes gently to mix
- 20 Dissociate tissue mechanically by passing through an 18 gauge needle 10 times
- 21 Incubate in a closed tube for 10 min at 37°C under slow, continuous rocking/rotation
- 22 Dissociate tissue again by passing through an 20 gauge needle 10 times
- 23 Incubate in a closed tube for 10 min at 37°C under slow, continuous rocking/rotation
- 24 Apply the cell suspension to a 40 m strainer placed on a 50 mL tube
- 25 Wash the filter with 10mL HBSS (with Ca^{2+} or Mg^{2+}) (Sigma 55037C)



- 26 Centrifuge cell suspension at 300 x g for 10 min at room temperature and aspirate supernatant
- 27 Resuspend pellet in 2 mL of warm Neurobasal Media and count cells
- 28 Plate 500,000 cells per well in a 6 well plate
- 29 Change ONLY 50% of the media every three days with Neurobasal media without glutamic acid