



Feb 25, 2025

Processing Serum, Plasma, Plasma EDTA from Whole Blood

DOI

<https://dx.doi.org/10.17504/protocols.io.eq2ly64wrgx9/v1>

Sara Lucas Del Pozo¹, David Chau¹, Jane Macnaughtan¹, Giuseppe Uras¹, Roxana Mezabrovschi¹

¹UCL and ASAP



Sofia Koletsi

University College London, University of London

Create & collaborate more with a free account

Edit and publish protocols, collaborate in communities, share insights through comments, and track progress with run records.

Create free account

OPEN  ACCESS



DOI: <https://dx.doi.org/10.17504/protocols.io.eq2ly64wrgx9/v1>

Collection Citation: Sara Lucas Del Pozo, David Chau, Jane Macnaughtan, Giuseppe Uras, Roxana Mezabrovschi 2025. Processing Serum, Plasma, Plasma EDTA from Whole Blood. **protocols.io**

<https://dx.doi.org/10.17504/protocols.io.eq2ly64wrgx9/v1>

License: This is an open access collection distributed under the terms of the [Creative Commons Attribution License](#), which permits unrestricted use, distribution, and reproduction in any medium, provided the original author and source are credited

Protocol status: Working

We use this collection and it's working



Created: February 05, 2025

Last Modified: February 25, 2025

Collection Integer ID: 119995

Keywords: ASAPCRN, plasma edta from whole blood, edta from whole blood sample, plasma edta, processing of serum, processing serum, whole blood sample, serum, edta, whole blood, processing, plasma, collection of protocol, blood, protocol

Funders Acknowledgements:

Aligning Science Across Parkinson's

Grant ID: ASAP-000420

Abstract

This collection of protocols describes the processing of serum, plasma, plasma EDTA from whole blood sample.

Guidelines

The protocol needs prior approval by the users' Institutional Review Board (IRB), Institutional Animal Care and Use Committee (IACUC) or equivalent ethics committee(s) as applicable.

Safety warnings

 The protocol needs prior approval by the users' Institutional Review Board (IRB), Institutional Animal Care and Use Committee (IACUC) or equivalent ethics committee(s) as applicable.

Before start

The protocol needs prior approval by the users' Institutional Review Board (IRB), Institutional Animal Care and Use Committee (IACUC) or equivalent ethics committee(s) as applicable.

Files

 SEARCH

Protocol

NAME

Processing Serum from Whole Blood

VERSION 1

CREATED BY



Sofia Koletsi
University College London, University of London

OPEN →

Protocol

NAME

Processing plasma from Whole Blood

VERSION 1

CREATED BY



Sofia Koletsi
University College London, University of London

OPEN →

Protocol

NAME

Dry blot spot cards - Processing plasma EDTA from Whole Blood

VERSION 1

CREATED BY



Sofia Koletsi
University College London, University of London

OPEN →