

Sep 17, 2024

Version 2

Processing wastewater samples for bacterial & viral targets enrichment V.2

DOI

<https://dx.doi.org/10.17504/protocols.io.4r3l2qb7pl1y/v2>

Victor Mabasa¹

¹NICD



Victor Mabasa

National Institute for Communicable Diseases

Create & collaborate more with a free account

Edit and publish protocols, collaborate in communities, share insights through comments, and track progress with run records.

Create free account

OPEN  ACCESS



DOI: <https://dx.doi.org/10.17504/protocols.io.4r3l2qb7pl1y/v2>

Protocol Citation: Victor Mabasa 2024. Processing wastewater samples for bacterial & viral targets enrichment . protocols.io <https://dx.doi.org/10.17504/protocols.io.4r3l2qb7pl1y/v2>Version created by [Victor Mabasa](#)

License: This is an open access protocol distributed under the terms of the [Creative Commons Attribution License](#), which permits unrestricted use, distribution, and reproduction in any medium, provided the original author and source are credited

Protocol status: Working

We use this protocol and it's working

Created: September 16, 2024



Last Modified: September 17, 2024

Protocol Integer ID: 107675

Keywords: wastewater genomics syndicate, processing wastewater sample, wastewater sample processing protocol, wastewater sample, supernatant components of wastewater, viral targets enrichment, wastewater supernatant, wastewater, nucleic acid extraction, pathogen, different pathogen, such as ultrafiltration, ultrafiltration, bacteria, virus, extraction, national institute of communicable disease, enrichment, processing

Funders Acknowledgements:

Bill & Melinda Gates Foundation

Abstract

This wastewater sample processing protocol by the Wastewater Genomics Syndicate at the National Institute of Communicable Diseases (NICD) is designed to enrich bacteria and viruses from the settled solids and supernatant components of wastewater, as different pathogens partition differently within this matrix. The end products are waterless solids and clarified wastewater supernatants, ready for downstream applications such as ultrafiltration, bead enrichment and nucleic acid extraction.

Materials

Equipment

–80°C freezer to store settled solids samples

4°C refrigerator storage of samples

Biological safety cabinet (Class II)

Refrigerated centrifuge (4°C) with adapters to accommodate 500 mL centrifuge bottles

Consumables

2 mL cryo-tubes

250 mL Conical centrifuge bottles

Labels

Waste container

Plain wooden sticks

Solutions/Reagents

70% ethanol

1% Virkon



Sample Collection & Handling

23h

- 1 Using correct personal protective equipment (PPE), collect a 1 000 mL grab wastewater sample from each location in a sterile bottle and transfer them to the laboratory in a cold chain (keep in a lab fridge at 4°C). It is important to proceed to the next step within the next 23 hours after sample collection to prevent major degradation of microorganisms.

23h



Preparing for centrifugation

10m

- 2 All work should be performed in a Class II biological safety cabinet while wearing appropriate PPE. Clean your workspace with a 1% Virkon solution and a 70% ethanol solution. Double-line your waste bucket with biohazard waste plastic, and ensure all your filter tips are refilled. Finally, label 250 mL conical centrifuge bottles, one for each wastewater sample.

10m

Centrifugation

50m

- 3 Transfer 250 mL of the wastewater samples to the corresponding conical tubes
- 4 Centrifuge at $4\,650 \times g$ (deceleration speed set at 5), at 4°C for 10 minutes.
- 5 Carefully transfer the supernatant from the 250 mL conical bottle to a clean, marked 1 000 mL bottle without disturbing the settled solids.
- 6 Repeat steps 3 to 5 until the original sample is finished.
- 7 Keep the clarified wastewater sample at 4°C for different downstream applications (check our page for different nucleic acid extraction protocols from clarified wastewater).

5m

15m

5m

20m

5m

Transfer settled solids

20m

- 8 Use clean, plain wooden sticks to pick out settled solids and pack them into marked 2 mL cryo-tubes. Centrifuge briefly to remove water.
- 9 Store the settled solids in the -80°C freezer until needed for downstream applications. (check our page for different nucleic acid extraction protocols from settled solids).

15m

5m

