

Oct 01, 2023

Protein Digestion and Mass Spectrometry Analysis Protocol

DOI

<https://dx.doi.org/10.17504/protocols.io.4r3l2261ql1y/v1>

Anthony Iavarone¹, Annan Cook^{1,2}

¹University of California, Berkeley; ²ASAP CRN



Annan SI Cook

University of California, Berkeley

Create & collaborate more with a free account

Edit and publish protocols, collaborate in communities, share insights through comments, and track progress with run records.

Create free account

OPEN  ACCESS



DOI: <https://dx.doi.org/10.17504/protocols.io.4r3l2261ql1y/v1>

Protocol Citation: Anthony Iavarone, Annan Cook 2023. Protein Digestion and Mass Spectrometry Analysis Protocol. protocols.io <https://dx.doi.org/10.17504/protocols.io.4r3l2261ql1y/v1>

License: This is an open access protocol distributed under the terms of the **Creative Commons Attribution License**, which permits unrestricted use, distribution, and reproduction in any medium, provided the original author and source are credited

Protocol status: Working

We use this protocol and it's working

Created: September 25, 2023



Last Modified: May 31, 2024

Protocol Integer ID: 88519

Keywords: ASAPCRN, protocol for the trypsin digestion, digested peptide, trypsin digestion, mass spectrometry analysis protocol this protocol detail, mass spectrometry analysis protocol, protein digestion, mass spectrometry, protein, pi3kc3, protocol detail

Funders Acknowledgements:

Aligning Science Across Parkinson's

Grant ID: ASAP-000350

Abstract

This protocol details a protocol for the trypsin digestion of PI3KC3-C1 and subsequent analysis of digested peptides using LC-MS/MS.

Attachments



[852-2203.pdf](#)

39KB

Materials

Materials

- PI3KC3-C1 (mCherry-ATG14|VPS15-TSF) protein sample
- Urea
- tris(hydroxymethyl)aminomethane (TRIS) pH 8.0
- tris(2-carboxyethyl)phosphine (TCEP)
- Iodoacetamide (IAA)
- Trypsin (Gold Trypsin, 0.5 mg/mL Promega V5280)
- CaCl_2
- $\text{NH}_4\text{CH}_3\text{CO}_2$
- Acetonitrile
- Formic acid (Optima LC-MS grade, 99.9% minimum)
- Water (purified using the Milli-Q Gradient system)
- Zorbax 300SB-C8 Micro Bore Rapid Resolution column (150 mm length, 1.0 mm inner diameter, 3.5 μm particle size, Agilent)
- LC-MS system (e.g., Agilent 1200 series LC and Thermo Fisher Scientific LTQ-Orbitrap-XL mass spectrometer)
- Proteome Discoverer software (version 1.3, Thermo Fisher Scientific)



Protein Denaturation and Reduction

20m

- 1 Mix PI3KC3-C1 (mCherry-ATG14|VPS15-TSF) protein with [M] 8 Mass Percent urea, [M] 50 millimolar (mM) TRIS pH 8.0 , and [M] 10 millimolar (mM) TCEP. 
- 2 Add Iodoacetamide (IAA) to achieve a final concentration of [M] 15 millimolar (mM) . 
- 3 Incubate the mixture at ° Room temperature for ⌚ 00:20:00 to alkylate cysteine residues. 

Trypsin Digestion Setup

- 4 Prepare a trypsin digestion solution containing Gold Trypsin (0.5 mg/mL), [M] 50 millimolar (mM) TRIS pH 8 , and [M] 100 millimolar (mM) CaCl_2 .

Trypsin digestion solution

	A	B
	Gold Trypsin (0.5 mg/mL)	8 uL
	50 mM TRIS pH 8	160 uL
	100 mM CaCl_2	2 uL

- 5 Mix the trypsin digestion solution with the alkylated protein sample. 

Digestion Incubation

20m

- 6 Incubate the digestion mixture at ° 37 °C ⌚ Overnight with shaking at ↺ 200 rpm .  

Confirmation by SDS-PAGE

- 7 Analyze the digested proteins by SDS-PAGE to confirm successful digestion. The absence of bands corresponding to the PI3KC3-C1 components indicates successful digestion. 

Liquid Chromatography (LC) Setup

- 8 Prepare LC mobile phase solvents: Solvent A and Solvent B.

Solvent A

	A
	99.9% water
	0.1% formic acid

Solvent B

	A
	99.9% acetonitrile
	0.1% formic acid

- 9 Set up the LC system with a Zorbax 300SB-C8 Micro Bore Rapid Resolution column (150 mm length, 1.0 mm inner diameter, 3.5 μ m particle size).
- 10 Maintain the column compartment at 50 °C .

LC-MS Analysis

2h

- 11 Load a 10 μ L sample onto the column.

- 12 Implement the following gradient elution program:

2h

- Isocratic flow at 1% (volume/volume) B for 00:02:00 .
- Linear gradient to 35% B over 01:30:00 .
- Linear gradient to 95% B over 00:01:00 .
- Isocratic flow at 95% B for 00:06:00 .
- Linear gradient to 1% B over 00:01:00 .
- Isocratic flow at 1% B for 00:20:00 .



- 13 Set the flow rate to  120 μL /min.

Mass Spectrometry Analysis

- 14 Acquire full-scan, high-resolution mass spectra in positive ion mode over the m/z range of 340 to 1800 using the Orbitrap mass analyzer with a mass resolution of 60,000 (at m/z = 400, FWHM).

Data-Dependent MS/MS Analysis

- 15 Select the ten most intense ions exceeding an intensity threshold of 10,000 raw ion counts from each full-scan mass spectrum for tandem mass spectrometry (MS/MS) analysis using collision-induced dissociation (CID).
- 16 Acquire MS/MS spectra using the linear ion trap.

Data Analysis

- 17 Use Proteome Discoverer software (version 1.3, SEQUEST algorithm) to search raw data files against the amino acid sequences of mCherry-labeled Class III phosphatidylinositol 3-kinase complex I (mCherry-PI3KC3-C1) proteins.
- 18 Consider tryptic peptides with up to three missed cleavages and specify post-translational modifications.
- 19 Validate peptide assignments by manually inspecting MS/MS spectra.