



Jul 10, 2024

Protein extraction and Western blotting

DOI

<https://dx.doi.org/10.17504/protocols.io.rm7vzb44xvx1/v1>

Shiyi Wang¹

¹Duke University



Shiyi Wang

Duke University

Create & collaborate more with a free account

Edit and publish protocols, collaborate in communities, share insights through comments, and track progress with run records.

Create free account

OPEN  ACCESS



DOI: <https://dx.doi.org/10.17504/protocols.io.rm7vzb44xvx1/v1>

Document Citation: Shiyi Wang 2024. Protein extraction and Western blotting. **protocols.io**
<https://dx.doi.org/10.17504/protocols.io.rm7vzb44xvx1/v1>

License: This is an open access document distributed under the terms of the **Creative Commons Attribution License**, which permits unrestricted use, distribution, and reproduction in any medium, provided the original author and source are credited

Created: May 23, 2023

Last Modified: July 10, 2024



Document Integer ID: 82326

Keywords: ASAPCRN, blotting protein extraction, protein extraction, extraction

Funders Acknowledgements:

Aligning Science Across Parkinson's (ASAP) initiative

Grant ID: ASAP-020607

Disclaimer

DISCLAIMER – FOR INFORMATIONAL PURPOSES ONLY; USE AT YOUR OWN RISK

The protocol content here is for informational purposes only and does not constitute legal, medical, clinical, or safety advice, or otherwise; content added to [protocols.io](https://www.protocols.io) is not peer reviewed and may not have undergone a formal approval of any kind. Information presented in this protocol should not substitute for independent professional judgment, advice, diagnosis, or treatment. Any action you take or refrain from taking using or relying upon the information presented here is strictly at your own risk. You agree that neither the Company nor any of the authors, contributors, administrators, or anyone else associated with [protocols.io](https://www.protocols.io), can be held responsible for your use of the information contained in or linked to this protocol or any of our Sites/Apps and Services.

Abstract

Protein extraction and Western blotting

1. Protein was extracted from cultured rat astrocytes using membrane solubilization buffer (25 mM Tris pH 7.4, 150 mM NaCl, 1 mM CaCl_2 , 1 mM MgCl_2 , 0.5% NP-40, and protease inhibitors).
2. Cultured astrocytes were washed twice with ice-cold TBS containing 1 mM CaCl_2 and 1 mM MgCl_2 and incubated on ice in membrane solubilization buffer for 20 minutes with occasional agitation.
3. Cell lysates were collected, vortexed briefly, and centrifuged at 4°C at high speed for 10 minutes to pellet non-solubilized material. The supernatant was collected and stored at -80°C.
4. Pierce BSA Protein Assay Kit (Thermo Fisher) was used to determine protein concentration, and lysates were mixed with 4× Pierce™ Reducing Sample Buffer (Thermo Scientific) and incubated at 45°C for 45 minutes to denature proteins.
5. 7-10 µg (cultured astrocyte lysates) of protein was loaded into Bolt™ 4-12% Bis-Tris Plus gels (Thermo Scientific) and run at 150 V for 1 hour.
6. Proteins were transferred at 100 V to PVDF membrane (Millipore) for 1.5 hours, blocked in 5% BSA in TBST (137 mM NaCl, 2.68 mM KCl, 24.7 mM Tris-Base, 0.1% (w/v) Tween 20) for 1 hour and incubated in primary antibodies overnight at 4°C.
7. Primary antibodies used were: anti-LRRK2 (Rabbit, 1:500; ab133474, abcam), GAPDH (mouse, 1:5000; ab8245, abcam), β-actin (mouse, 1:5000; A5441, Millipore Sigma), phospho-ERM (Rabbit, 1:500; #3141, Cell Signaling), ERM (Rabbit, 1:500; #3142, Cell Signaling), phospho-Rab10 (Rabbit, 1:500; ab230261, abcam), Rab10 (Rabbit, 1:500; ab237703, abcam).
8. The next day, membranes were washed with TBST, incubated in HRP secondary antibodies (Thermo Fisher Scientific) for 2 hours, washed in TBST, and imaged on a Biorad Gel Doc imaging system.
9. Protein levels were quantified using FIJI.