

May 10, 2024

Protein Extraction for Amyloid Beta Fractionation

DOI

<https://dx.doi.org/10.17504/protocols.io.j8nlk8ky1l5r/v1>

Lisa Blackmer-Raynolds¹, Tim Sampson¹

¹Emory University



Lisa Blackmer-Raynolds

Emory University

Create & collaborate more with a free account

Edit and publish protocols, collaborate in communities, share insights through comments, and track progress with run records.

Create free account

OPEN  ACCESS



DOI: <https://dx.doi.org/10.17504/protocols.io.j8nlk8ky1l5r/v1>

Protocol Citation: Lisa Blackmer-Raynolds, Tim Sampson 2024. Protein Extraction for Amyloid Beta Fractionation. **protocols.io** <https://dx.doi.org/10.17504/protocols.io.j8nlk8ky1l5r/v1>

License: This is an open access protocol distributed under the terms of the **Creative Commons Attribution License**, which permits unrestricted use, distribution, and reproduction in any medium, provided the original author and source are credited

Protocol status: Working

We use this protocol and it's working

Created: May 09, 2024



Last Modified: May 10, 2024

Protocol Integer ID: 99473

Keywords: amyloid beta, fractionation, solubility, protein, protein extraction for amyloid beta fractionation, amyloid beta fractionation, evaluating amyloid beta pathology, insoluble amyloid beta, protein extraction, amyloid beta, associated amyloid beta, different amyloid beta structure, protein fraction, amyloid beta pathology as different amyloid beta structure, protein from murine brain tissue, soluble fraction, amyloid plaque, formic acid soluble fraction, extraction, protein, fraction, solubility, murine brain tissue, other two fraction, brain tissue, hippocampal sample, tris soluble

Funders Acknowledgements:

NIH

Grant ID: F31AG076332-02

ASAP

Grant ID: ASAP-020527

Abstract

This protocol can be used to isolate protein from murine brain tissue (volumes listed are for hippocampal samples) by solubility. It creates three protein fractions: tris soluble, triton soluble, and formic acid soluble. This protocol can be helpful when evaluating amyloid beta pathology as different amyloid beta structures can be found within each fraction. For example, insoluble amyloid beta associated with amyloid plaques will be found within the formic acid soluble fraction, whereas more soluble non-plaque associated amyloid beta will be found within the other two fractions.



Preparation

- 1 **5x Homogenization Buffer (store 4°C):**
125 ml of 1.0 M Tris
30 ml of 0.5 M MgCl₂
25 ml of 0.1 M EDTA (pH to 7.2 and store at 4°C for up to 6 months)
20ml of MiliQ Water
- 2 **1x Homogenization Buffer (dilute day of):**
Dilute 5x Homogenization buffer to 1x in MiliQ Water and add 1 tablet protease inhibitor (cat#11697498001 Roche) per 10 mL
- 3 **1x Homogenization Buffer with 1% Triton X (dilute day of):**
Add 1% Triton X-100 to 1x Homogenization buffer made above
- 4 **70% Formic Acid (store RT) **Caution Acidic****
7mL formic acid
3mL water
- 5 **1.5x Homogenization Buffer pH 14 (store 4°C) **Caution Corrosive****
Dilute 5x Homogenization buffer to 1.5x in MiliQ Water and pH to 14

Tissue Processing

- 6 Weigh frozen tissue sample and record for later normalization
- 7 Homogenize the tissue on ice using quick bursts of a sonicator in 400µl 1x Homogenization buffer
- 8 Centrifuge maximum speed at 4°C for 1hr
- 9 Collect supernatant as tris soluble fraction and store -80°C for later use
- 10 Homogenize the pellet on ice using quick bursts of a sonicator in 200µl 1x Homogenization Buffer with 1% Triton X-100
- 11 Centrifuge maximum speed at 4°C for 1hr



- 12 Collect supernatant as triton soluble fraction and store -80°C for later use
- 13 Homogenize the pellet on ice using quick bursts of a sonicator in 50µl 70% formic acid
Be careful with acid
- 14 Centrifuge maximum speed at 4°C for 1hr
- 15 Collect supernatant as formic acid soluble fraction and neutralize with 1.5x
Homogenization buffer pH14 on ice *Be careful with acid and base**. Store -80°C for
later use