



Jul 11, 2019

Protein interaction analysis of KaiC3 with various Kai homologs via yeast two-hybrid experiments (Growth Assay)

DOI

dx.doi.org/10.17504/protocols.io.wcnfave

Anika Wiegard¹, Christin Köbler², Katsuaki Oyama³, Anja K. Dörrich⁴, Chihiro Azai^{5,3}, Kazuki Terauchi^{5,3}, Annegret Wilde², Ilka Maria IM Axmann⁶

¹Heinrich-Heine Universität Düsseldorf;

²Institute of Biology III, Faculty of Biology, University of Freiburg, 79104 Freiburg, Germany;

³Graduate School of Life Sciences, Ritsumeikan University, Kusatsu, Shiga 525-8577, Japan;

⁴Institute for Microbiology and Molecular Biology, Justus-Liebig University, 35392 Giessen, Germany;

⁵College of Life Sciences, Ritsumeikan University, Kusatsu, Shiga 525-8577, Japan;

⁶Institute for Synthetic Microbiology, Cluster of Excellence on Plant Sciences (CEPLAS), Heinrich Heine University Duesseldorf, 40225 Duesseldorf, Germany

CyanoWorld



Christin Köbler

Create & collaborate more with a free account

Edit and publish protocols, collaborate in communities, share insights through comments, and track progress with run records.

Create free account

OPEN  ACCESS





DOI: <https://dx.doi.org/10.17504/protocols.io.wcnfave>

Protocol Citation: Anika Wiegard, Christin Köbler, Katsuaki Oyama, Anja K. Dörrich, Chihiro Azai, Kazuki Terauchi, Annegret Wilde, Ilka Maria IM Axmann 2019. Protein interaction analysis of KaiC3 with various Kai homologs via yeast two-hybrid experiments (Growth Assay). **protocols.io** <https://dx.doi.org/10.17504/protocols.io.wcnfave>

License: This is an open access protocol distributed under the terms of the **[Creative Commons Attribution License](#)**, which permits unrestricted use, distribution, and reproduction in any medium, provided the original author and source are credited

Protocol status: Working

We use this protocol and it's working

Created: December 11, 2018

Last Modified: July 11, 2019

Protocol Integer ID: 18542

Keywords: protein interaction analysis of kaic3, various kai homologs via yeast, protein interaction via yeast, various kai homolog, protein interaction analysis, protein interaction, kaic3, yeast cell, protein, yeast, hybrid experiment, activity of the histidine, growth assay

Abstract

This protocol can be used to investigate protein-protein interaction via yeast two-hybrid experiments. It describes the yeast-two hybrid method relying on the activity of the histidine-synthetase. If the proteins interact, the yeast cells are able to grow on selective medium without histidine.



Materials

List of Materials

- Yeast cells (AH109)
- Frozen-EZ Yeast Transformation II Kit (Zymo Research)
- Yeast Nitrogen Base without amino acids (Formedium, CYN0401)
- Drop-out mixture (-Leu -Trp; MP Biomedicals, 114520012)
- Drop-out mixture (-Leu -Trp -his; MP Biomedicals, 114530112)
- Adenine-hemisulfate (Sigma-Aldrich, A9126)
- 3-amino-1,2,4-triazole (3-AT, Sigma-Aldrich, A8056)
- Bacto Agar (BD Diagnostics, 214010)
- D-Glucose (Roth, X997.2)

Complete supplement medium (CSM)

	Components for CSM Agar	
	Yeast Nitrogen Base with ammonium /without amino acids	6.7 g/L
	D-Glucose	20 g/L
	Bacto Agar	20g/L
	Drop-Out-Mix (amino acid mixture)	0.60-0.64 g/L
	Adenine Hemisulfate	50 mg/L
	dd H ₂ O	

- Autoclave 15 min, 121°C or filter sterilize before using

Troubleshooting

Preparation of buffers and media

- 1 Prepare
 - CSM -Leu -Trp agar (complete supplement mixture lacking leucine and tryptophan)
 - CSM -Leu -Trp -His agar (complete supplement mixture lacking leucine, tryptophan and histidine)
 - 3-AT solution, 1M

Transformation of yeast cells

- 2 Perform the transformation of yeast cells according to manufacturer's guidelines using the Frozen-EZ Yeast Transformation II Kit (Zymo Research) and select transformed cells on complete supplement mixture lacking leucine and tryptophan (CSM -Leu -Trp) at 30 °C for 3–4 days.

Restreaking of colonies

- 3
 - Prepare plates: CSM -Leu -Trp -His + 3-AT (interaction plate)
 - Determine the appropriate concentration of 3-AT by adding different concentrations of 3-AT to the interaction plate
 - Possible concentrations of 3-AT: 0 mM, 3 mM, 6.25 mM, 12.5 mM, 25 mM and 50mM
 - Let the agar cool down before adding 3-AT, it is heat sensitive
 - Mark and restreak 1-3 formed colonies from step 2 onto the interaction plates
suggestion: only use a small amount of cells for restreaking to reduce false positives
 - Wrap parafilm around the plates
 - Incubate plates at 30°C for 3-6 days
 - Use the lowest concentration of 3-AT that allows only small (<1-mm) colonies to grow after 3-4 days for further interaction plates

Growth assay

- 4
 - Prepare plates: CSM -Leu -Trp (control) and CSM -Leu -Trp -His + 3-AT (interaction plate), use determined concentration of 3-AT
 - Restreak 1 colony from step 2 onto the control plate and afterwards onto the interaction plate (reduces amount of cells and therefore false positives)
 - Wrap parafilm around the plates
 - Incubate plates at 30°C for 3-4 days
 - Result: Interaction of two proteins is positive if there is growth on both, the control and interaction plate, but without any growth of the controls on the interaction plate.