



Mar 19, 2025

Protein Sample Preparation from Acrylamide Gel for Mass Spectrometry

DOI

<https://dx.doi.org/10.17504/protocols.io.36wgg6r33lk5/v1>

Patricia Yuste-Checa¹, Barbara Steigenberger², Andreas Bracher¹, F Ulrich Hartl¹

¹Department of Cellular Biochemistry, Max Planck Institute of Biochemistry, Martinsried, Germany;

²Mass Spectrometry Core Facility, Max Planck Institute of Biochemistry, Martinsried, Germany



Patricia Yuste-Checa

MPI Biochemistry

Create & collaborate more with a free account

Edit and publish protocols, collaborate in communities, share insights through comments, and track progress with run records.

Create free account

OPEN  ACCESS



DOI: <https://dx.doi.org/10.17504/protocols.io.36wgg6r33lk5/v1>

Protocol Citation: Patricia Yuste-Checa, Barbara Steigenberger, Andreas Bracher, F Ulrich Hartl 2025. Protein Sample Preparation from Acrylamide Gel for Mass Spectrometry. **protocols.io**

<https://dx.doi.org/10.17504/protocols.io.36wgg6r33lk5/v1>

License: This is an open access protocol distributed under the terms of the [Creative Commons Attribution License](#), which permits unrestricted use, distribution, and reproduction in any medium, provided the original author and source are credited

Protocol status: Working

We use this protocol and it's working

Created: March 11, 2025

Last Modified: March 19, 2025

Protocol Integer ID: 124566

Keywords: ASAPCRN, protein sample preparation from acrylamide gel, protein samples from polyacrylamide gel, mass spectrometry proteomic, protein sample preparation, protein sample, mass spectrometry, mass spectrometry this protocol detail, acrylamide gel, polyacrylamide gel, sample, proteomic, sample preparation

Funders Acknowledgements:

Aligning Science Across Parkinson's

Grant ID: ASAP-000282

Abstract

This protocol details how to prepare protein samples from polyacrylamide gels for mass spectrometry proteomics.

Materials

Digestion Buffer:

	A	B
	Tris-HCl	25 mM
	Acetonitrile	10%
	Trypsin	10 ng μl^{-1}



Protein Sample Preparation from Acrylamide Gel

- 1 Slice the protein band of interest from a Coomassie blue-stained polyacrylamide gel using a scalpel.

Note

Avoid contamination by human keratin by using clean apparel and reagents. Wear gloves.

- 2 Wash thoroughly the gel pieces with  150 μL of de-staining buffer ( 25 millimolar (mM) ammonium bicarbonate, 50% ethanol). 
- 3 Dehydrate the gel pieces with  150 μL of pure ethanol.
- 4 Remove the ethanol and dry the gel pieces using a SpeedVac.
- 5 Add  50 μL of digestion buffer (25 mM Tris-HCl, 10% acetonitrile, 10 ng μL^{-1} trypsin). 

Digestion Buffer:

	A	B
	Tris-HCl	25 mM
	Acetonitrile	10%
	Trypsin	10 ng μL^{-1}

- 6 Cool  On ice for  00:20:00 .

20m

- 7 Add  50 μL of  25 millimolar (mM) ammonium bicarbonate buffer. 



- 8 Incubate Overnight at 37 °C . 8h
- 9 Collect the peptides released to the supernatant.
- 10 Extract additional peptides by incubation at 25 °C for 00:20:00 in 100 µL of extraction buffer (3% trifluoroacetic acid (TFA), 30% acetonitrile). 20m
- 11 Centrifuge at 4000 rpm, 00:02:00 . 2m
- 12 Collect the supernatant. Repeat steps 10-12 twice.
- 13 Dehydrate the gel pieces with 100 µL of pure acetonitrile at 25 °C during 00:10:00 . 10m
- 14 Centrifuge at 4000 rpm, 00:02:00 . 2m
- 15 Collect the supernatant and combine with previous supernatants.
- 16 Remove the acetonitrile using a SpeedVac.
- 17 Add 70 µL of 2 Mass Percent Tris-HCl pH 8.0, 10 millimolar (mM) TCEP, and 40 millimolar (mM) chloroacetamide (CAA).
- 18 Incubate 00:30:00 at 37 °C . 30m
- 19 Acidify the peptides to 1% TFA.



- 20 For the desalting of the peptides, prepare a SDB-XC stage tip by pushing 3 plugs of 0.8 mm diameter from C18/SDB-XC and push them into a 300 µl pipette tip.
- 21 Wash with  20 µL of methanol to complete dryness by spinning. 
- 22 Wash with  20 µL of buffer B (0.1% Formic acid and 80% acetonitrile in water). Spin the tips but make sure that the tip is not dry. A small volume of buffer B should be visible on the top. 
- 23 Equilibrate the tips 2 times with  200 µL of buffer A (0.1% formic acid in water).
- 24 Load the sample onto the stage tips and spin at  1500 x g . 
- 25 Wash the tip with  100 µL of buffer A and spin to dryness. 
- 26 Elute the peptides from the tip using  60 µL of buffer B.
- 27 After elution, speed-vac the samples at  30 °C for  00:40:00 and store at  -20 °C until loading on the LC-MS system. 