

Oct 21, 2023

Proteolytic Peptide Desalting with C18 Hydrophilic–Lipophilic Balance (HLB) Cartridges

DOI

<https://dx.doi.org/10.17504/protocols.io.3byl4q382vo5/v1>

J P Rose¹, M Watson¹, B Schilling¹, Joanna Bons¹

¹Buck Institute for Research on Aging



M A Watson

Buck Institute

Create & collaborate more with a free account

Edit and publish protocols, collaborate in communities, share insights through comments, and track progress with run records.

Create free account

OPEN  ACCESS



DOI: <https://dx.doi.org/10.17504/protocols.io.3byl4q382vo5/v1>

Protocol Citation: J P Rose, M Watson, B Schilling, Joanna Bons 2023. Proteolytic Peptide Desalting with C18 Hydrophilic–Lipophilic Balance (HLB) Cartridges. **protocols.io** <https://dx.doi.org/10.17504/protocols.io.3byl4q382vo5/v1>

License: This is an open access protocol distributed under the terms of the **Creative Commons Attribution License**, which permits unrestricted use, distribution, and reproduction in any medium, provided the original author and source are credited

Protocol status: Working

We use this protocol and it's working



Created: October 21, 2023

Last Modified: October 21, 2023

Protocol Integer ID: 89689

Keywords: Desalting, HLB, Peptides, Proteomics, Mass Spectrometry, cartridges desalting of proteolytic peptide elution, proteolytic peptide desalting with c18 hydrophilic, proteolytic peptide elution, proteolytic peptide desalting, downstream proteomic profiling, using c18 hydrophilic, lipophilic balance, c18 hydrophilic, cartridges in preparation, cartridges desalting

Abstract

Desalting of proteolytic peptide elution using C₁₈ Hydrophilic–Lipophilic Balance (HLB) Cartridges in preparation for downstream proteomic profiling.

Materials

- Oasis HLB 1cc cartridges, 10 mg sorbent (Waters, Cat. #WAT058882)
- Vacuum manifold
- Centrifuge
- Centrifugal vacuum concentrator
- 1.5-mL microcentrifuge tubes
- Condition buffer/Elution buffer: 50% acetonitrile (ACN), 0.2% formic acid (FA) in water
- Equilibration buffer/Wash buffer: 0.2% formic acid (FA) in water
- HPLC-grade water

- 1 Centrifuge the samples at 1,850 x *g* for 5 minutes at room temperature to pellet insoluble material.
- 2 Desalt the samples using Oasis HLB solid-phase extraction cartridges placed on top of a vacuum manifold as follows:
 - 2.1 Condition each cartridge two times with 800 µL of the condition buffer (50% ACN, 0.2% FA).
 - 2.2 Equilibrate each cartridge three times with 800 µL of the equilibration buffer (0.2% FA).
 - 2.3 Load the peptide samples.
 - 2.4 Wash each cartridge three times with 800 µL of the wash buffer (0.2% FA).
 - 2.5 Elute peptides with 800 µL of the elution buffer (50% ACN, 0.2% FA) followed by a further 400 µL of the same elution buffer.
- 3 Vacuum dry the eluted peptide solution in a centrifugal vacuum concentrator.
- 4 Reconstitute the dried peptides in 0.2% FA for a final protein concentration of 1 µg/µL and thoroughly mix the solution.
- 5 Centrifuge at 12,000 x *g* at room temperature for min, and store at -20°C.