

Mar 29, 2024

# Proteomic Analysis of Human Ovarian Cortex and Medulla Secretome Using Liquid Chromatography-Tandem Mass Spectrometry (LC-MS/MS) Acquisition by Data-Independent Acquisition (DIA) on an Orbitrap Eclipse Tribrid Mass Spectrometer

DOI

[dx.doi.org/10.17504/protocols.io.n92ld83bnv5b/v1](https://dx.doi.org/10.17504/protocols.io.n92ld83bnv5b/v1)

Joanna Bons<sup>1</sup>, J P Rose<sup>1</sup>, M A Watson<sup>1</sup>, B Schilling<sup>1</sup>

<sup>1</sup>Buck Institute for Research on Aging

Metabolomics Protocols ...

Cellular Senescence Net...



**M A Watson**

Buck Institute

## Create & collaborate more with a free account

Edit and publish protocols, collaborate in communities, share insights through comments, and track progress with run records.

Create free account

OPEN  ACCESS



DOI: <https://dx.doi.org/10.17504/protocols.io.n92ld83bnv5b/v1>

**Protocol Citation:** Joanna Bons, J P Rose, M A Watson, B Schilling 2024. Proteomic Analysis of Human Ovarian Cortex and Medulla Secretome Using Liquid Chromatography-Tandem Mass Spectrometry (LC-MS/MS) Acquisition by Data-Independent Acquisition (DIA) on an Orbitrap Eclipse Tribrid Mass Spectrometer. **protocols.io**

<https://dx.doi.org/10.17504/protocols.io.n92ld83bnv5b/v1>

**Manuscript citation:**

[dx.doi.org/10.17504/protocols.io.e6nvw1px7lmk/v](https://dx.doi.org/10.17504/protocols.io.e6nvw1px7lmk/v)

**License:** This is an open access protocol distributed under the terms of the **[Creative Commons Attribution License](#)**, which permits unrestricted use, distribution, and reproduction in any medium, provided the original author and source are credited

**Protocol status:** Working

**We use this protocol and it's working**

**Created:** March 25, 2024

**Last Modified:** April 08, 2024

**Protocol Integer ID:** 97359

**Keywords:** Human, Ovary, Secretome, SASP, Senescence, Doxorubicin, LC-MS/MS, DIA, Orbitrap Eclipse, Mass Spectrometry, Conditioned Media, proteomic analysis of human ovarian cortex, human ovarian cortex, secretome proteomics profiling, orbitrap eclipse tribrid mass spectrometer ovary, hours for secretome proteomics profiling, proteomic analysis, proteolytic peptide measurement, tandem mass spectrometry, ml doxorubicin, medulla secretome, medulla explant, using liquid chromatography

**Funders Acknowledgements:**

**NIH Grant**

Grant ID: 5U54AG075932-02

**NIH Grant**

Grant ID: 5U54AG075932-03

## Disclaimer

The **protocols.io** team notes that research involving animals and humans must be conducted according to internationally-accepted standards and should always have prior approval from an Institutional Ethics Committee or Board.



## Abstract

Ovaries from human donors were cut into 3–5 mm sections. These sections were further processed into 500  $\mu\text{m}$  slices containing cortex and medulla. The slices were then processed into pieces (1 mm x 1 mm x 500  $\mu\text{m}$ ), and cortex and medulla pieces were cultured separately as explants in static cultures.

Explants were cultured and treated with either DMSO vehicle control or 0.1  $\mu\text{g}/\text{mL}$  doxorubicin to induce senescence. After 10 days, cortex and medulla explants were thoroughly washed with serum-free basal media and transferred to a clean plate with pre-equilibrated serum-free basal media and inserts. The conditioned media were collected after 24 hours for secretome proteomics profiling.

The concentrated conditioned media was subjected to tryptic digestion using S-trap Spin columns. The reconstituted peptide elution was desalted with C18 hydrophilic-lipophilic balance (HLB) cartridges. The final reconstituted peptides were diluted with 2% ACN and 0.1% FA. Proteolytic peptide measurement was completed using liquid chromatography-tandem mass spectrometry (LC-MS/MS) acquisition by Data-Independent Acquisition (DIA) on an Orbitrap Eclipse Tribrid mass spectrometer for peptide/protein identification and quantification.

## Troubleshooting



## Conditioned Media Concentration with Amicon Ultra Centrifugal Filters

1 [dx.doi.org/10.17504/protocols.io.e6nvw1px7lmk/v1](https://dx.doi.org/10.17504/protocols.io.e6nvw1px7lmk/v1)

## Protein Digestion with S-trap Spin Columns using Conditioned Concentrated Media

2 [dx.doi.org/10.17504/protocols.io.x54v928eml3e/v1](https://dx.doi.org/10.17504/protocols.io.x54v928eml3e/v1)

## Proteolytic Peptide Desalting with C18 HLB Cartridges

3 [dx.doi.org/10.17504/protocols.io.eq2lywdzpvx9/v1](https://dx.doi.org/10.17504/protocols.io.eq2lywdzpvx9/v1)

## LC-MS/MS Acquisition by DIA on an Orbitrap Eclipse Tribrid Mass Spectrometer

4 [dx.doi.org/10.17504/protocols.io.36wgq3m15lk5/v1](https://dx.doi.org/10.17504/protocols.io.36wgq3m15lk5/v1)

## DIA Data Processing using Spectronaut/directDIA (Biognosys): Secretome Analysis

5 [dx.doi.org/10.17504/protocols.io.q26g71rz3gwz/v1](https://dx.doi.org/10.17504/protocols.io.q26g71rz3gwz/v1)