

Apr 20, 2024

Version 2

Protocol for Mouse IP Anesthesia (Ketamine-Xylazine) V.2

DOI

<https://dx.doi.org/10.17504/protocols.io.14egn9qg5d/v2>



Haowei Li¹, Markus Holzl¹, Mohsen Khosravi-Maharlooeei¹, Nichole Danzl¹, Austin Chen¹, Megan Sykes¹

¹Columbia Center for Translational Immunology, Columbia University, New York

Human Islet Research N...



Sandy Beshir

City of Hope

Create & collaborate more with a free account

Edit and publish protocols, collaborate in communities, share insights through comments, and track progress with run records.

Create free account

OPEN  ACCESS



DOI: <https://dx.doi.org/10.17504/protocols.io.14egn9qg5d/v2>

Protocol Citation: Haowei Li, Markus Holzl, Mohsen Khosravi-Maharlooeei, Nichole Danzl, Austin Chen, Megan Sykes 2024. Protocol for Mouse IP Anesthesia (Ketamine-Xylazine). **protocols.io**

<https://dx.doi.org/10.17504/protocols.io.14egn9qg5d/v2> Version created by **Sandy Beshir**

License: This is an open access protocol distributed under the terms of the **[Creative Commons Attribution License](#)**, which permits unrestricted use, distribution, and reproduction in any medium, provided the original author and source are credited

Protocol status: Working

We use this protocol and it's working

Created: April 20, 2024

Last Modified: April 20, 2024

Protocol Integer ID: 98525

Keywords: protocol for mouse ip anesthesia, mouse ip anesthesia, anesthesia cocktail, ketamine, xylazine, kidney capsule, protocol, protocol detail, mice, thymectomy, mouse

Funders Acknowledgements:

NIH

Grant ID: U01 DK123559

Abstract

This protocol details the constitution of our anesthesia cocktail that we use for sedating our mice when performing surgeries. The surgeries we typically perform are thymectomies and implantation of material under the kidney capsule (through open abdomen or minimally invasive approaches).

Note

The changes of this protocol was made by Austin Chen on Markus Holzl original protocol.

Note

Corresponding Authors

Mohsen Khosravi-Maharlooei

Email: mkm2182@cumc.columbia.edu

Austin Chen

Email: ac4274@cumc.columbia.edu

Tel: 425-283-6900



Materials

Materials:

Anesthesia is a mixture of ketamine and xylazine in 1x PBS.

- **Ketamine:** Ketathesia (100 mg/mL; cat: 056344; Henry Schein)
- **Xylazine:** Xylazine sterile solution (20 mg/mL; Columbia ICM)
- **1x PBS:** Phosphate Buffered Solution (1X) (cat: **BW17516F24**; Lonza)
- **0.22 um filter:** Syringe Filter Unit, 0.22 µm, polyethersulfone (**Millex-GP**; cat: **SLGPR33RS**)

Preparation:

Ketamine stock: 100 mg/mL

Xylazine stock: 20 mg/mL

Desired ketamine concentration: 0.08 mg/gram of mouse

Desired xylazine concentration: 0.012 mg/gram of mouse

- We will make anesthesia solution to dose  10 µL anesthesia/gram of mouse for IP injections

For 10 mL of anesthesia:

1.  0.8 mL Ketamine (100 mg/mL) +  0.6 mL Xylazine (20 mg/mL) +  8.6 mL 1x PBS
2. Filter solution with 0.22 um Millex-GP filter

Calculations:

- Ketamine: 0.8 mL * 100 mg/mL in 10 mL → 8 mg/mL Ketamine → 0.08 mg/10 uL
- Xylazine 0.6 mL * 20 mg/mL in 10 mL → 1.2 mg/mL Xylazine → 0.012 mg/10 uL



For Intraperitoneal (IP) Approach:

- 1 Use 10 uL/gram of anesthesia

Comments:

1. It takes time until the mice fall asleep

2. Very often, mice do not fall asleep but move around like zombies. This requires another small dose of IP ketamine-xylazine injection or a small dose of IV injection

- a. When doing IP injection again, we have injected IP with a volume of about 1/3 of original volume

- b. For IV approach after IP, it takes very little anesthesia for them to fall asleep. We usually use around 10-20 uL. Inject this slowly into the mouse. The mouse should fall asleep rapidly. It is very easy to overdose with IV injection so be careful.