

Jul 03, 2024

Version 1

Protocols for Recinto et al. "A rewiring of the earliest immune events leading to T-cell mediated disease following intestinal microbial infection in a PINK1 KO mouse model of Parkinson's disease" V.1

DOI

<https://dx.doi.org/10.17504/protocols.io.kxygxy77ol8j/v1>

Sherilyn Junelle Recinto¹, adam.macdonald¹, Moein Yaqubi¹, Alexandra Kazanova²

¹Department of Neurology and Neurosurgery, Montreal Neurological Institute-Hospital, McGill University, Montréal, QC, Canada;

²Department of Microbiology and Immunology, McGill University, Montréal, QC, Canada



Lilia Rodriguez

Université de Montréal

Create & collaborate more with a free account

Edit and publish protocols, collaborate in communities, share insights through comments, and track progress with run records.

Create free account

OPEN  ACCESS



DOI: <https://dx.doi.org/10.17504/protocols.io.kxygxy77ol8j/v1>

Collection Citation: Sherilyn Junelle Recinto, adam.macdonald , Moein Yaqubi, Alexandra Kazanova 2024. Protocols for Recinto et al. "A rewiring of the earliest immune events leading to T-cell mediated disease following intestinal microbial infection in a PINK1 KO mouse model of Parkinson's disease". protocols.io <https://dx.doi.org/10.17504/protocols.io.kxygxy77ol8j/v1>

License: This is an open access collection distributed under the terms of the **Creative Commons Attribution License**, which permits unrestricted use, distribution, and reproduction in any medium, provided the original author and source are credited

Protocol status: Working

We use this collection and it's working

Created: June 20, 2024

Last Modified: July 30, 2024

Collection Integer ID: 102180

Keywords: ASAPCRN, immune cell type in pink1 ko, initiating immune event, dysregulated immune cell type, rewiring of the earliest immune event, pink1 ko mouse model of parkinson, earliest immune event, targeted pathogen, immune events at the site, induced pd mouse modeling system, pink1 ko mice with gram, enhanced proinflammatory profile, pink1 ko mice, infection at the earliest stage, cell interaction pathway, pink1 ko mouse model, pd mouse modeling system, intestinal microbial infection, mediated disease, antigen presentation, mice, following intestinal microbial infection, proinflammatory profile, infection, myeloid cell, capacity for antigen presentation, cell, disease, parkinson

Funders Acknowledgements:

Aligning Science Across Parkinson's

Grant ID: ASAP-000525

Disclaimer

The protocols.io team notes that research involving animals and humans must be conducted according to internationally-accepted standards and should always have prior approval from an Institutional Ethics Committee or Board.

Abstract

Our group has developed GI-targeted pathogen-induced PD mouse modeling systems (in PINK1 KO mice with gram negative bacterial infections) and found that T cells are a major player in driving PD-like motor symptoms at late stages following infection. Herein, we now map the initiating immune events at the site of infection at the earliest stages with the goal of shedding light on the earliest mechanisms triggering T cell-mediated pathological processes relevant to PD. Using unbiased single cell sequencing, we demonstrate that myeloid cells are the earliest dysregulated immune cell type in PINK1 KO infected mice (at 1-week post-infection) followed by a dysregulated T cell response shortly after (at 2 weeks post-infection). We find that these myeloid cells have an enhanced proinflammatory profile, are more mature, and develop enhanced capacity for antigen presentation. Using unbiased prediction analysis, our data suggests that cytotoxic T cells and myeloid cells are particularly poised for interacting with each other, and we identify possible direct cell-cell interaction pathways that might be implicated.



Disclaimer

The [protocols.io](#) team notes that research involving animals and humans must be conducted according to internationally-accepted standards and should always have prior approval from an Institutional Ethics Committee or Board.

Files

 SEARCH



SR2024_Immunophenotyping_flowcytometry.docx

DOWNLOAD ↓



SR2024_gutLP_prep.docx

DOWNLOAD ↓



SR2023_LipocalinELISA.docx

DOWNLOAD ↓



SR2022_10XChromium single cell 3'RNAseq preparation.docx

DOWNLOAD ↓



SR2024_RNAextraction+qPCR.docx

DOWNLOAD ↓



SR2024_Immunophenotyping_flowcytometry.docx

DOWNLOAD ↓



Procedure_Citrobacter rodentium infection model SOP.docx

DOWNLOAD ↓