Nov 07, 2024

🌐 Purificación de GFP en reacciones cell-free para actividad docente // GFP purification from cell-free reactions for educational purpose

DOI

<https://dx.doi.org/10.17504/protocols.io.6qpvr81nplmk/v1>

Severine Cazaux¹, Valentina Ferrando¹, Paola Larrauri¹, Quinn Matthews², Mohammad Simchi², Severino Jefferson Ribeiro Da Silva², Keith Pardee², Fernan Federici³

¹Pontificia Universidad Católica; ²University of Toronto; ³PUC ReClone

Laboratorio de Tecnolog...

Severine Cazaux

Pontificia Universidad Católica

Create & collaborate more with a free account

Edit and publish protocols, collaborate in communities, share insights through comments, and track progress with run records.

Create free account

OPEN  ACCESS

DOI: <https://dx.doi.org/10.17504/protocols.io.6qpvr81nplmk/v1>

Protocol Citation: Severine Cazaux, Valentina Ferrando, Paola Larrauri, Quinn Matthews, Mohammad Simchi, Severino Jefferson Ribeiro Da Silva, Keith Pardee, Fernan Federici 2024. Purificación de GFP en reacciones cell-free para actividad docente // GFP purification from cell-free reactions for educational purpose. **protocols.io**

<https://dx.doi.org/10.17504/protocols.io.6qpvr81nplmk/v1>

Manuscript citation:

This section will be updated when the manuscript is published (currently in writing stage).

License: This is an open access protocol distributed under the terms of the **Creative Commons Attribution License**, which permits unrestricted use, distribution, and reproduction in any medium, provided the original author and source are credited

Protocol status: Working

We use this protocol and it's working

Created: October 22, 2024

Last Modified: November 07, 2024

Protocol Integer ID: 110482

Keywords: gfp purification from cell, gfp purification, basics of protein purification, protein purification, proteína con materiales que, purification, protein of interest, traditional lab infrastructure, protein, free reactions for educational purpose este, una proteína, centrifugation step, gfp

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This protocol describes an educational activity to explain the basics of protein purification with materials that can be used without traditional lab infrastructure. In few words, the protein of interest (in this case, 6xHis-sfGFP) is produced in cell-free lysates overnight. The day after, it is purified using NEBExpress Spin Columns and all the centrifugation steps (for binding, washing and elution) are performed using a 3D-printed hand-powered centrifuge. The activity lasts two hours.

Materials

Hardware:

- 3D-printed hand centrifuge
- 500 lb fishing line ([link](#))
- Neodymium magnets with ¼ inch diameter and 1/8 inch thickness ([link](#)).
- Blue LED Transilluminator

Wetware:

- 500uL cell-free reactions of the GFP to be purified

Solutions for purification (described in NEBExpress Ni Spin column kit NEB #S1427):

- Binding Buffer: 20 mM sodium phosphate, 300 mM NaCl, pH 7.4
- Wash Buffer: 20 mM sodium phosphate, 300 mM NaCl, 5 mM Imidazole, pH 7.4
- Elution Buffer: 20 mM sodium phosphate, 300 mM NaCl, 500 mM Imidazole, pH 7.4
- Spin columns from NEB loaded with Ni-NTA resin

Disposables:

- 2mL tubes for recollection of the fractions
- Pipette tips 1000uL and 200uL
- Parafilm + Scissors

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#####

Add 20nM of plasmid DNA sfGFP to a 500uL cell-free reaction in a 5mL centrifuge tube:
167 uL of cell-free lysate
133 uL of CFPS supplementation solutions
20nM of DNA
Top up to 500uL with nuclease-free water

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#####


Incubate overnight at 30°C and 80 rpm (it can handle up to 37°C and 220rpm but with higher risk of protein aggregation).

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####

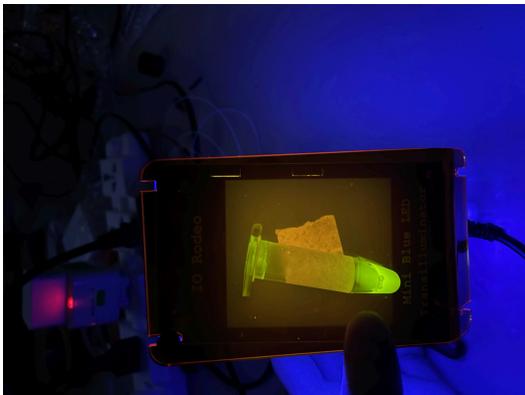Check the fluorescence of the cell-free reaction in the blue led transilluminator

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
####

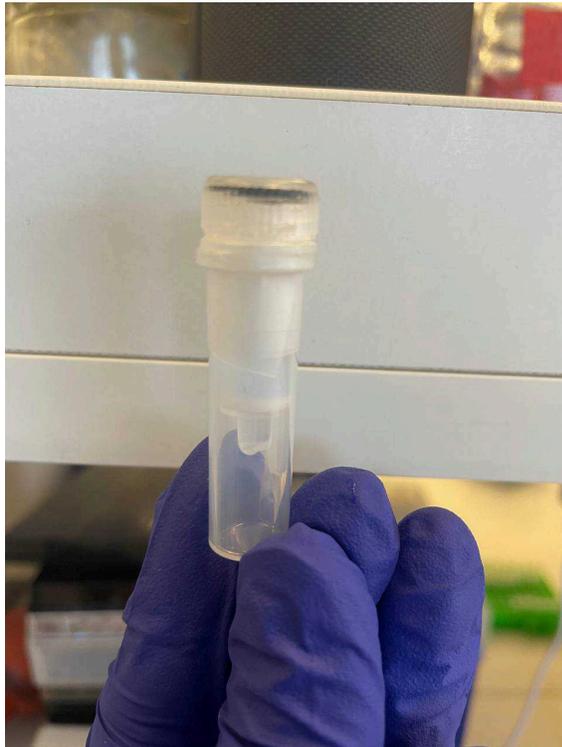Get the Ni Spin columns ready to use. To do that, remove the bottom tab, loosen the cap of the column and place it into a 2mL collection tube. At this point, the column is filled with storage buffer.

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####

Seal the joint between the collection tube and the column with parafilm to avoid the liquid to pour out during centrifugation.

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
####

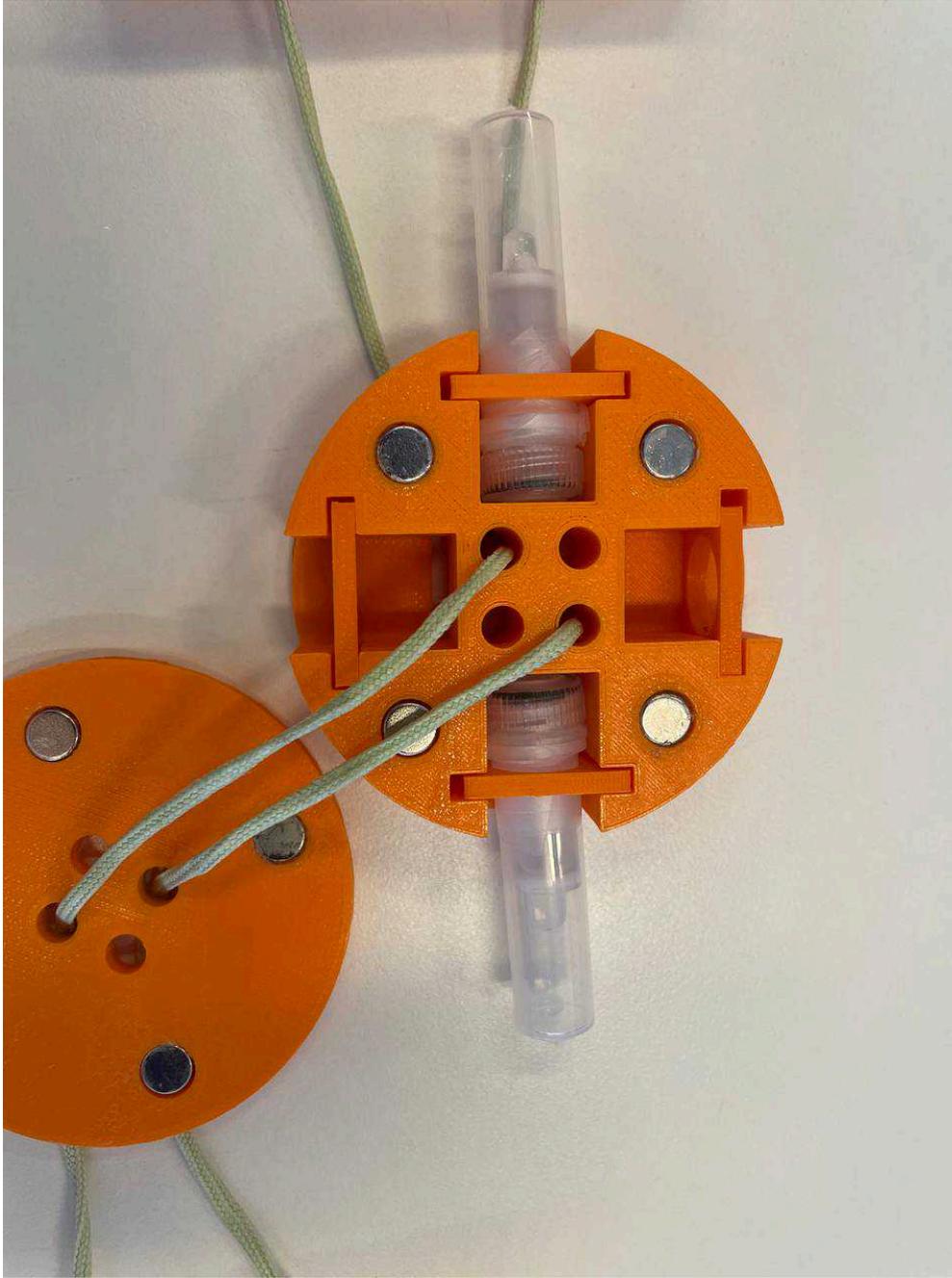The centrifuge must always be used equilibrated, which means that opposing sides should have the same weight. Therefore, make sure to work with two spin columns even if the second one has no protein (you can fill it with water to match the content of the other column).

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####

Place the two columns in opposing sides of the centrifuge, with the collection tube pointing toward the outside and the cap toward the center of the centrifuge.

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####

Centrifuge. To do that, coil the fishing line as much as possible then stretch it to start centrifuging. Check the progress every 10s by stopping the centrifuge in a super-coiled configuration so that it is easier to start again. Once all the liquid has gone through the resin and reached the bottom of the collection tube, stop centrifuging.

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####

Still in a coiled configuration, open the centrifuge and remove the columns.

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####

Remove the parafilm and the cap. Place the columns in a new collection tube.

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####

Add 250 uL of Binding Solution. Place the cap back with new parafilm between the collection tube and the column.

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####

Add 500uL of the cell-free reaction containing the GFP. Place the cap back with new parafilm between the collection tube and the column.

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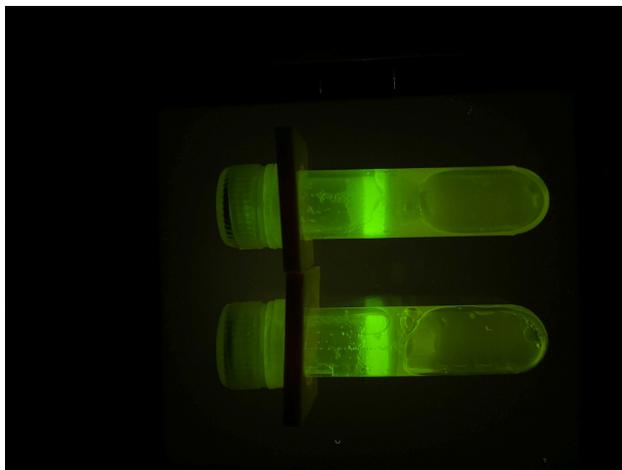
####

Look at the column in the transilluminator. The resin is now loaded with GFP, which means the 6x-His tag has bound to the Ni²⁺ in the resin.

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####

Note: Steps 18-21 are optional for educational purposes. If time lacks these steps can perfectly be skipped.

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####

Add 250uL of Washing Solution. Place the cap back with new parafilm between the collection tube and the column.

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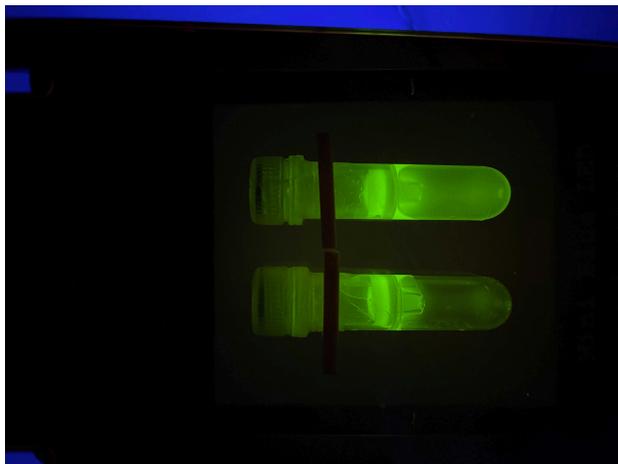
####

Look at the column in the transilluminator. The protein should still be bound to the resin with no or minimal loss in the flow-through.

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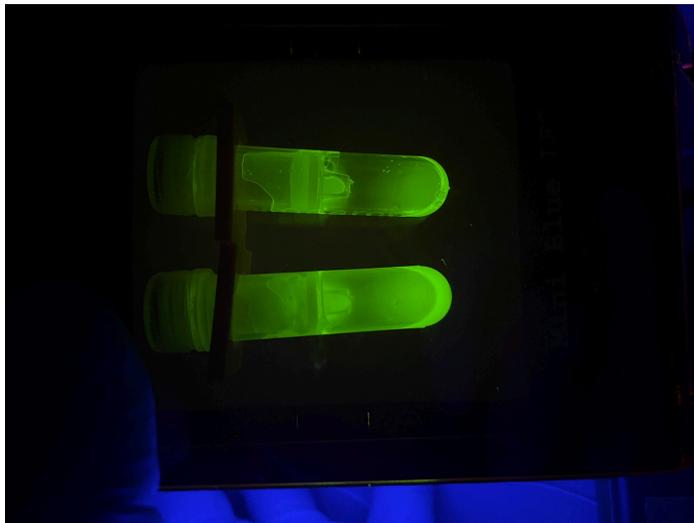
####

Look at the column in the transilluminator. The high concentration of imidazole during the elution step should have detached the protein from the resin, which should now be located at the bottom of the collection tube. The resin should look transparent. If relevant amounts of fluorescence are observed in the resin, a second step of elution can be performed.

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
####

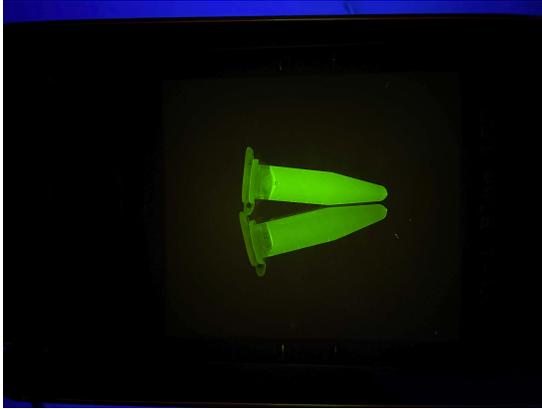Optional: You can transfer the flow-through to a new centrifuge tube for higher clarity. You have purified GFP!

Instrucciones para el ensamblaje de la centrífuga // Instructions for building the 3D-printed, hand-powered centrifuge

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####

Print the .stl files attached using PLA filament:

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####

Glue the magnets in the holes in both the tap and the body of centrifuge (8 magnets x centrifuge). Make sure to use a strong glue as the magnetic force can be more powerful than the glue.

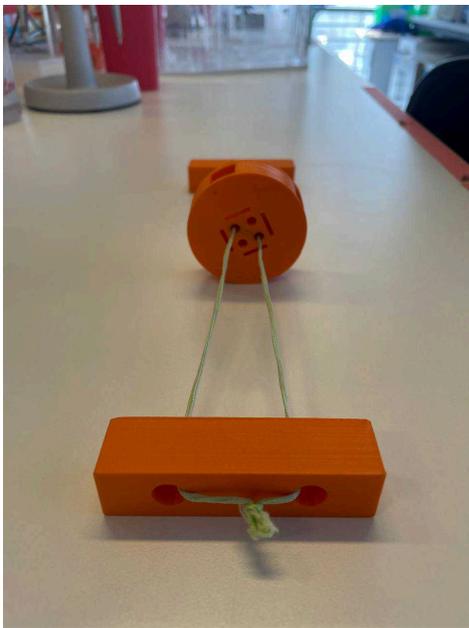
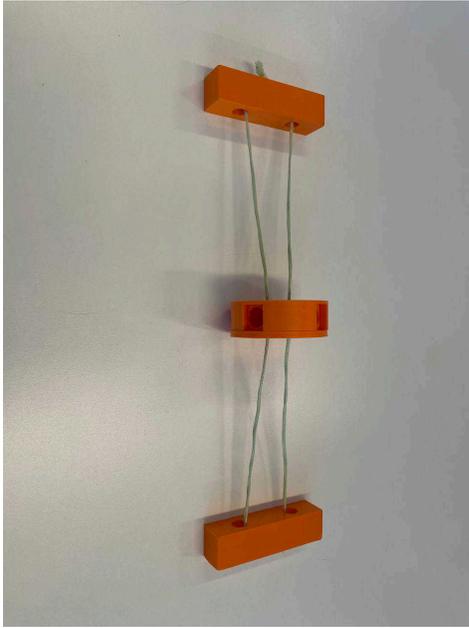
Mind the polarity of the magnets! Any mistake can prevent the correct assembly of the centrifuge.

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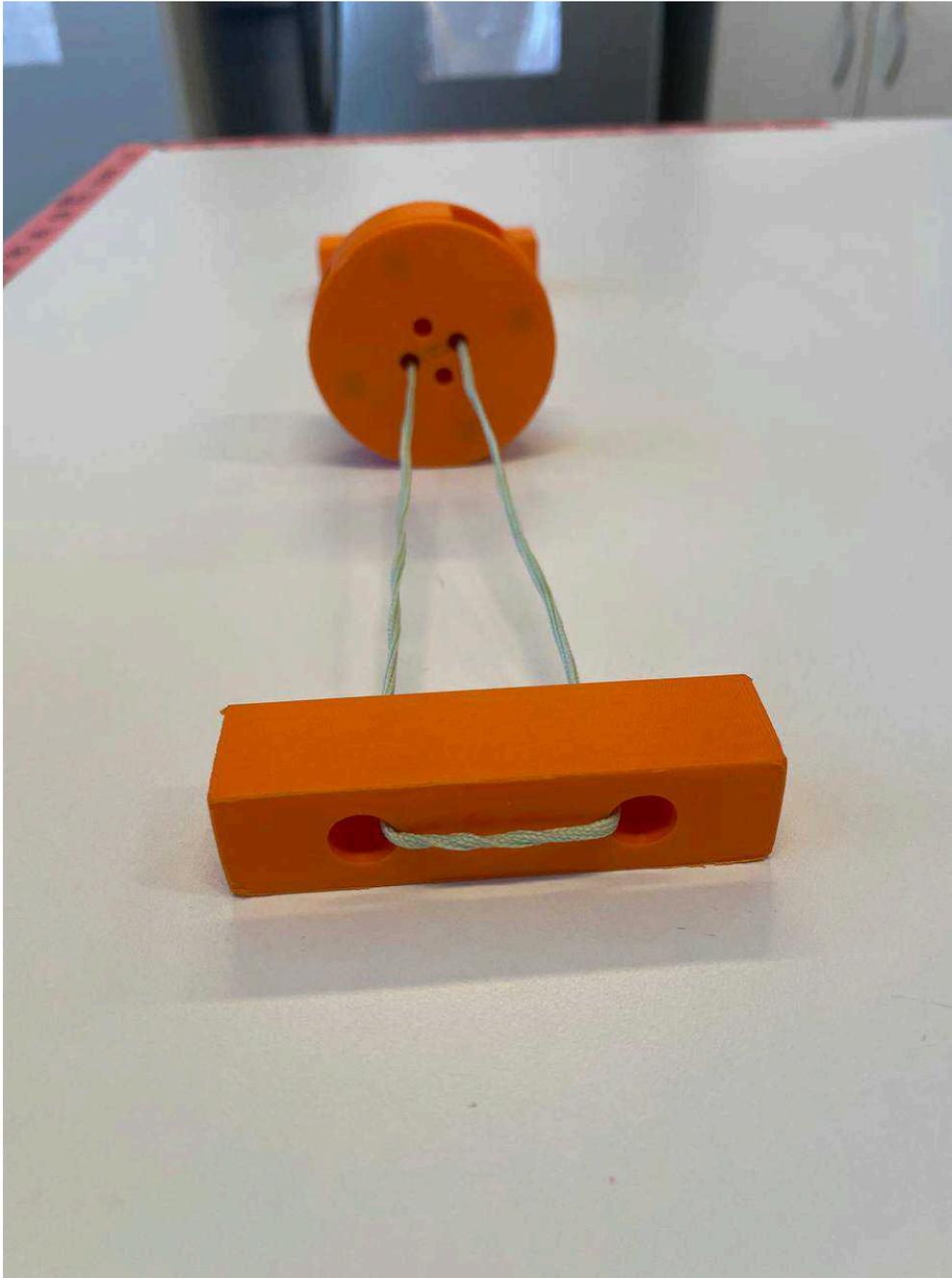
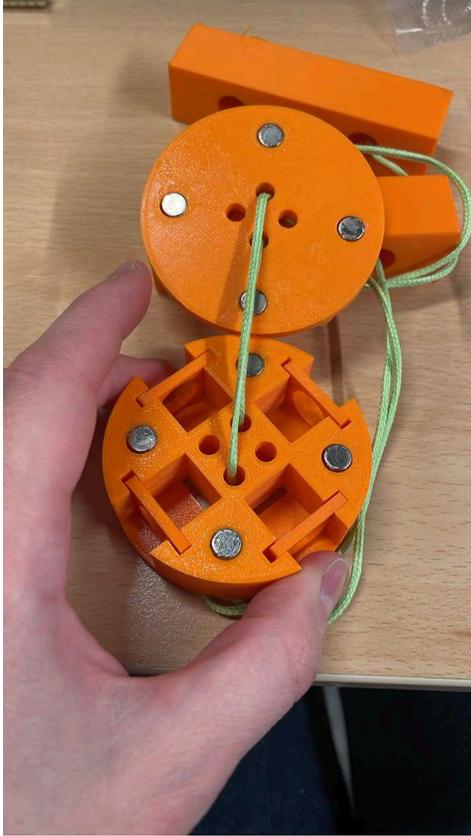
####

Insert the fishing line through the holes of the centrifuge as follows. Use a solid knot: when operating the centrifuge, the knot could unfold/break! (It has happened to us - twice during the same activity!).

Protocol references

1. Byagathvalli G, Pomerantz A, Sinha S, Standeven J, Bhamla MS (2019) A 3D-printed hand-powered centrifuge for molecular biology. PLOS Biology 17(5): e3000251.
2. IO Rodeo LED Transilluminator
https://iorodeo.github.io/led_transilluminator_docs/?ref=blog.iorodeo.com
2. NEBExpress Ni Spin Columns (NEB #S1427):
<https://doi.org/10.1371/journal.pbio.3000251> <https://www.neb.com/en/protocols/2019/08/28/nebexpress-ni-spin-column-reaction-protocol-neb-s1427>

Acknowledgements

As part of this protocol we want to acknowledge specific contributions of the following authors:

Mohammad Simchi modified the 3D-fuge by replacing the nut-and-bolt closure with a magnetic closure.

Quinn Matthews adapted the 3D-printed centrifuge (3D-fuge) for cell-free protein purification and designed the initial purification protocol.

Severino Jefferson Ribeiro da Silva provided the 3D-fuge to Federici's lab through IDRC collaborative project.