



Aug 29, 2024

Purification of FUNDC1-GFP

 [Nature Cell Biology](#)

DOI

<https://dx.doi.org/10.17504/protocols.io.4r3l2qnx3l1y/v1>

Elias Adriaenssens¹

¹Sascha Martens lab, University of Vienna, Max Perutz Labs - Vienna



Elias Adriaenssens

Sascha Martens lab, University of Vienna, Max Perutz Labs - ...

Create & collaborate more with a free account

Edit and publish protocols, collaborate in communities, share insights through comments, and track progress with run records.

Create free account

OPEN  ACCESS



DOI: <https://dx.doi.org/10.17504/protocols.io.4r3l2qnx3l1y/v1>

External link: <https://doi.org/10.1038/s41556-025-01712-y>

Protocol Citation: Elias Adriaenssens 2024. Purification of FUNDC1-GFP. **protocols.io**

<https://dx.doi.org/10.17504/protocols.io.4r3l2qnx3l1y/v1>

Manuscript citation:

Adriaenssens, E., Schaar, S., Cook, A.S.I. *et al.* Reconstitution of BNIP3/NIX-mitophagy initiation reveals hierarchical flexibility of the autophagy machinery. *Nat Cell Biol* **27**, 1272–1287 (2025). <https://doi.org/10.1038/s41556-025-01712-y>

License: This is an open access protocol distributed under the terms of the **Creative Commons Attribution License**, which permits unrestricted use, distribution, and reproduction in any medium, provided the original author and source are credited

Protocol status: Working

We use this protocol and it's working

Created: May 24, 2024

Last Modified: August 29, 2024

Protocol Integer ID: 103091

Keywords: ASAPCRN, purification of fundc1, gfp this protocol, purification, fundc1, gfp, protocol detail, protocol

Funders Acknowledgements:

Aligning Science Across Parkinson's (ASAP)

Grant ID: ASAP-000350

Marie Skłodowska-Curie MSCA Postdoctoral fellowship

Grant ID: 101062916

Abstract

This protocol details the purification of FUNDC1-GFP.

Materials

Lysis buffer:

	Tris-HCl	50 mM
	pH	7.4
	NaCl	300 mM
	Triton X-100	1%
	glycerol	5%
	MgCl ₂	2 mM
	DTT	1 mM
	β-mercaptoethanol	2mM
	cOmplete EDTA-free protease inhibitors (Roche)	
	CIP protease inhibitor (Sigma)	
	DNase (Sigma)	

Wash buffer:

	Tris-HCl	50 mM
	pH	7.4
	NaCl	300 mM
	DTT	1 mM

High salt buffer:

	Tris-HCl	50 mM
	pH	7.4

	NaCl	700 mM
	DTT	1 mM

SEC buffer:

	Tris-HCl	25 mM
	pH	7.4
	NaCl	300 mM
	DTT	1 mM

- FUNDC1-GFP (available from Addgene) or FUNDC1(Y18A/L21A)-GFP (Δ LIR) (available from Addgene)
- pET-DUET1 vector (available on Addgene)  pETDuet-1 TIM9,10 **addgene Catalog #170280**
- FUNDC1 Y18A/L21A (Δ LIR)(available on Addgene)
- Rosetta pLysS cells (Novagen Cat# 70956-4)

 Rosetta™(DE3)pLysS Competent Cells - Novagen **Merck Catalog #70956-4**
- SORVAL RC6+ centrifuge with an F21S-8×50Y rotor (Thermo Scientific)
- 10 kDa cut-off Amicon filter (Merck Millipore)

 Amicon® Ultra Centrifugal Filter, 10 kDa MWCO **Merck MilliporeSigma (Sigma-Aldrich) Catalog #UFC801008**



Purification - FUNDC1-GFP

20h 46m

- 1 To purify GFP-tagged
 - FUNDC1-GFP (available from Addgene) or FUNDC1(Y18A/L21A)-GFP (Δ LIR) (available from Addgene),

fuse the cytosol-exposed domain of FUNDC1 (1-50aa) to a C-terminal GFP-tag through cloning into a pET-DUET1 vector (available on Addgene).
- 2 Introduce the point mutants by in vitro mutagenesis to generate
 - FUNDC1 Y18A/L21A (Δ LIR)(available on Addgene).
- 3 After the transformation of the pET-DUET1 vector encoding FUNDC1-GFP wild-type or mutants in *E. coli* Rosetta pLysS cells (Novagen Cat# 70956-4), grow the cells in 2x Tryptone Yeast extract (TY) medium at  37 °C until an OD₆₀₀ of 0.4 and then continue at  18 °C .
- 4 Once the cells reaches an OD₆₀₀ of 0.8, induce protein expression with  100 micromolar (μ M) isopropyl β -D-1-thiogalactopyranoside (IPTG) for  16:00:00 at  18 °C .
- 5 Collect the cells by centrifugation and resuspend in lysis buffer.

Lysis buffer:

A	B
Tris-HCl pH 7.4	50 mM
NaCl	300 mM
Triton X-100	1%
Glycerol	5%
MgCl ₂	2 mM
DTT	1 mM
β -mercaptoethanol	2mM
cComplete EDTA-free protease inhibitors (Roche)	
CIP protease inhibitor (Sigma)	



A	B
DNase (Sigma)	

6 Sonicate the cell lysates twice for 30 s and clear by centrifugation at  18000 rpm, 4°C, 00:45:00 in a SORVAL RC6+ centrifuge with an F21S-8×50Y rotor (Thermo Scientific).

45m



6.1 Sonicate the cell lysates for  00:00:30 (1/2).

30s

6.2 Sonicate the cell lysates for  00:00:30 (2/2).

30s

7 Collect the supernatant and incubate with pre-equilibrated Glutathione Sepharose 4B beads (GE Healthcare) for  02:00:00 at  4 °C with gentle shaking to bind FUNDC1-GFP.

2h



8 Centrifuge the samples to pellet the beads and remove the unbound lysate.



9 Wash the beads twice with wash buffer, once with high salt wash buffer, and two more times with wash buffer.



Wash buffer:

A	B
Tris-HCl pH 7.4	50 mM
NaCl	300 mM
DTT	1 mM

High salt buffer:

A	B
Tris-HCl pH 7.4	50 mM
NaCl	700 mM



A	B
DTT	1 mM

10 Cleave off the GST-tag  Overnight by eluting the GFP-tagged cargo receptor from the GSH beads by the addition of TEV protease in wash buffer at  4 °C



11 To collect the supernatant, collect the beads by centrifugation.



12 Wash the beads twice with  4 mL of wash buffer, and collect the supernatant.



13 Pool the supernatant fractions, filtered through a 0.45 µm syringe filter, concentrated with 10 kDa cut-off Amicon filter (Merck Millipore), and load onto a pre-equilibrated Superdex 200 Increase 10/300 GL column (Cytiva).

14 Elute the proteins with SEC buffer.

SEC buffer:

A	B
Tris-HCl pH 7.4	25 mM
NaCl	300 mM
DTT	1 mM

15 Analyze fractions by SDS-PAGE and Coomassie staining. Pool the fractions containing purified FUNDC1-GFP.

16 After concentrating the purified protein, aliquote the protein and snap-frozen in liquid nitrogen.



Note

Store the proteins at  -80 °C .

