




Apr 01, 2022

Purification of lysosomes using dextran-conjugated SPIONs

 [Journal of Cell Biology](#)

DOI

dx.doi.org/10.17504/protocols.io.bp2l61dr1vqe/v1

William Hancock-Cerutti^{1,2,3,4,5}, Arun Kumar Tharkeshwar¹, Shawn Ferguson^{1,5}, Pietro De Camilli^{1,2,3,5}

¹Departments of Neuroscience and of Cell Biology, Yale University School of Medicine, New Haven, Connecticut 06510, USA;

²Howard Hughes Medical Institute;

³Program in Cellular Neuroscience, Neurodegeneration and Repair, Yale University School of Medicine, New Haven, Connecticut 06510, USA;

⁴Interdisciplinary Neuroscience Program and MD-PhD Program, Yale University School of Medicine, New Haven, Connecticut 06510, USA;

⁵Aligning Science Across Parkinson's (ASAP) Collaborative Research Network, Chevy Chase, MD, 20815



William Hancock-Cerutti

Create & collaborate more with a free account

Edit and publish protocols, collaborate in communities, share insights through comments, and track progress with run records.

Create free account

OPEN  ACCESS



DOI: <https://dx.doi.org/10.17504/protocols.io.bp2l61dr1vqe/v1>



External link: <https://doi.org/10.1083/jcb.202106046>

Protocol Citation: William Hancock-Cerutti, Arun Kumar Tharkeshwar, Shawn Ferguson, Pietro De Camilli 2022. Purification of lysosomes using dextran-conjugated SPIONs. **protocols.io** <https://dx.doi.org/10.17504/protocols.io.bp2l61dr1vqe/v1>

Manuscript citation:

William Hancock-Cerutti, Zheng Wu, Peng Xu, Narayana Yadavalli, Marianna Leonzino, Arun Kumar Tharkeshwar, Shawn M. Ferguson, Gerald S. Shadel, Pietro De Camilli (2022) ER-lysosome lipid transfer protein VPS13C/PARK23 prevents aberrant mtDNA-dependent STING signaling. *Journal of Cell Biology* doi: [10.1083/jcb.202106046](https://doi.org/10.1083/jcb.202106046)

License: This is an open access protocol distributed under the terms of the **Creative Commons Attribution License**, which permits unrestricted use, distribution, and reproduction in any medium, provided the original author and source are credited

Protocol status: Working

We use this protocol and it's working

Created: April 01, 2022

Last Modified: May 31, 2024

Protocol Integer ID: 60196

Keywords: ASAPCRN, purification of lysosome, lysosome fraction from cultured cell, conjugated superparamagnetic iron nanoparticle, superparamagnetic iron nanoparticle, lysosome, lysosome fraction, spions this method, conjugated spion, using dextran, cultured cell, purification, spion, usage of dextran, dextran

Abstract

This method describes the usage of dextran-conjugated superparamagnetic iron nanoparticles (SPIONs) to purify a lysosome fraction from cultured cells.

Attachments



HC-Protocol-
Lysosome...

16KB

Troubleshooting

