

Sep 24, 2023

Purification of NDP52 (untagged)

DOI

<https://dx.doi.org/10.17504/protocols.io.36wgg35xklk5/v1>

Elias Adriaenssens¹

¹Sascha Martens lab, University of Vienna, Max Perutz Labs - Vienna



Elias Adriaenssens

Sascha Martens lab, University of Vienna, Max Perutz Labs - ...

Create & collaborate more with a free account

Edit and publish protocols, collaborate in communities, share insights through comments, and track progress with run records.

Create free account

OPEN  ACCESS



DOI: <https://dx.doi.org/10.17504/protocols.io.36wgg35xklk5/v1>

Protocol Citation: Elias Adriaenssens 2023. Purification of NDP52 (untagged). protocols.io
<https://dx.doi.org/10.17504/protocols.io.36wgg35xklk5/v1>

License: This is an open access protocol distributed under the terms of the **Creative Commons Attribution License**, which permits unrestricted use, distribution, and reproduction in any medium, provided the original author and source are credited

Protocol status: Working

We use this protocol and it's working

Created: July 07, 2023

Last Modified: June 01, 2024

Protocol Integer ID: 84642

Keywords: Purification of NDP52 (untagged), ASAPCRN, purification of ndp52, purification, ndp52, protocol

Funders Acknowledgements:

Aligning Science Across Parkinson's (ASAP)

Grant ID: ASAP-000350

Marie Skłodowska-Curie MSCA Postdoctoral fellowship

Grant ID: 101062916

Abstract

This protocol describes purification of NDP52 (untagged).

Attachments



[775-1961.pdf](#)

56KB

Materials

Materials

- Human NDP52 cDNA in pGST2 vector with an N-terminal GST tag (RRID:Addgene #187828)
- IPTG
- SORVAL RC6+ centrifuge with an F21S8×50Y rotor (Thermo Scientific)
- Glutathione Sepharose 4B beads (GE Healthcare)
- 30 kDa cut-off Amicon filter (Merck Millipore)
- Superdex 200 Increase 10/300 GL column (Cytiva)

Lysis buffer

	A	B
	Tris-HCl pH 7.4	50 mM
	NaCl	300 mM
	DTT	1 mM
	MgCl ₂	2 mM
	β-mercaptoethanol	2 mM
	cOmplete EDTA-free protease inhibitors (Roche)	
	DNase (Sigma)	

Wash buffer

	A	B
	Tris-HCl pH 7.4	50 mM
	NaCl	300 mM
	DTT	1 mM

High salt wash buffer

	A	B
	Tris-HCl pH 7.4	50 mM



	A	B
	NaCl	700 mM
	DTT	1 mM

SEC buffer

	A	B
	Tris-HCl pH 7.4	25 mM
	NaCl	150 mM
	DTT	1 mM



Purification of NDP52 (untagged)

16h

- 1 The human NDP52 cDNA into a pGST2 vector with an N-terminal GST tag followed by a TEV cleavage site is available from Addgene (RRID: Addgene #187828).
- 2 After the transformation of the pGST2 vector encoding GST-TEV-NDP52 in *E. coli* Rosetta pLySS cells, grow cells in 2xTY medium at 37 °C until an OD₆₀₀ of 0.4 and then continued at 18 °C .
- 3 Once the cells reached an OD₆₀₀ of 0.8, induce protein expression with 50 micromolar (μM) IPTG for 16:00:00 at 18 °C . 16h
- 4 Collect cells by centrifugation and resuspend in lysis buffer.
- 5 Sonicate cell lysates.
- 5.1 Sonicate cell lysates for 00:00:30 . (1/2) 30s
- 5.2 Sonicate cell lysates for 00:00:30 . (1/2) 30s
- 6 Clear lysates by centrifugation at 18000 rpm, 4°C, 00:45:00 in a SORVAL RC6+ centrifuge with an F21S8×50Y rotor (Thermo Scientific). 45m
- 7 Collect the supernatant and incubate with preequilibrated Glutathione Sepharose 4B beads (GE Healthcare) for 02:00:00 at 4 °C with gentle shaking to bind GST-NDP52. 2h
- 8 Centrifuge the samples to pellet the beads and remove the unbound lysate.
- 9 Wash the beads.
- 9.1 Wash the beads twice with wash buffer.



9.2 Wash the beads with high salt wash buffer.

9.3 Again, wash the beads twice with wash buffer.

10 Incubate beads  Overnight with TEV protease at  4 °C .



11 After the GST tag was cleaved off, filter the protein through a  0.45 µm syringe filter, concentrated using a 30 kDa cut-off Amicon filter (Merck Millipore), and load onto a pre-equilibrated Superdex 200 Increase 10/300 GL column (Cytiva).

12 Elute proteins with SEC buffer.

13 Analyze fractions by SDS-PAGE and Coomassie staining.



14 Pool fractions containing purified NDP52.

15 After concentrating the purified protein, aliquot the protein and snap-freeze in liquid nitrogen. Store the proteins at  -80 °C .

