

Nov 14, 2023

Version 1

QMask Hemispheric Separation V.1

 In 1 collection

DOI

dx.doi.org/10.17504/protocols.io.n2bvj3qdbl5/v1

Michael X. Henderson¹

¹Van Andel Institute

Michael X. Henderson: ORCID: 0000-0001-9710-0726



Michael Henderson

Van Andel Research Institute

Create & collaborate more with a free account

Edit and publish protocols, collaborate in communities, share insights through comments, and track progress with run records.

Create free account

OPEN  ACCESS



DOI: <https://dx.doi.org/10.17504/protocols.io.n2bvj3qdbl5/v1>

Protocol Citation: Michael X. Henderson 2023. QMask Hemispheric Separation. **protocols.io**
<https://dx.doi.org/10.17504/protocols.io.n2bvj3qdbl5/v1>

License: This is an open access protocol distributed under the terms of the **Creative Commons Attribution License**, which permits unrestricted use, distribution, and reproduction in any medium, provided the original author and source are credited

Protocol status: Working

We use this protocol and it's working

Created: August 16, 2023

Last Modified: May 31, 2024

Protocol Integer ID: 90663

Keywords: ASAPCRN, qmask hemispheric separation this protocol, qmask hemispheric separation, hemispheric separation, qmask

Funders Acknowledgements:

Aligning Science Across Parkinson's

Grant ID: ASAP-020616

National Institute on Aging

Grant ID: R01-AG077573

Abstract

This protocol describes QMask hemispheric separation.

Attachments




[812-2118.pdf](#)

4.6MB

Troubleshooting

QMask Hemispheric Separation

- 1 Open the Qmask tool and click "pick XML" and open the **QuickN XML** file generated in QuickNII.
- 2 Enter the appropriate coordinates from the **QMask Coordinates file**.
- 3 Click "Destination" and navigate to the QVN\Mask folder. Click "Go".

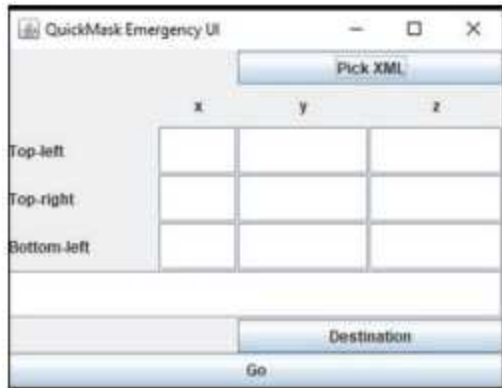

The mask output should be a black and White PNG that looks something like this:



- 4 Rename the mask files using the appropriate naming convention. Example: NG18I_s061_mask.