

Sep 10, 2024

R2C2 protocol using 10x Single-Cell RNA-seq Assay cDNA

DOI

<https://dx.doi.org/10.17504/protocols.io.81wgbzonygpk/v1>

Xian Adiconis¹, Joshua Levin¹

¹Broad Institute of MIT and Harvard



xian Adiconis

Broad Institute

Create & collaborate more with a free account

Edit and publish protocols, collaborate in communities, share insights through comments, and track progress with run records.

Create free account

OPEN  ACCESS



DOI: <https://dx.doi.org/10.17504/protocols.io.81wgbzonygpk/v1>

Protocol Citation: Xian Adiconis, Joshua Levin 2024. R2C2 protocol using 10x Single-Cell RNA-seq Assay cDNA. **protocols.io** <https://dx.doi.org/10.17504/protocols.io.81wgbzonygpk/v1>

Manuscript citation:

Simmons SK, Adiconis X, Haywood N, Parker J, Lin Z, Liao Z, Tuncali I, Al'Khafaji AM, Shin A, Jagadeesh K, Gosik K, Gatzem M, Smith JT, El Kods DN, Kuras Y, Baecher-Allan C, Serrano GE, Beach TG, Garimella K, Rozenblatt-Rosen O, Regev A, Dong X, Scherzer CR, Levin JZ. Experimental and Computational Methods for Allelic Imbalance Analysis from Single-Nucleus RNA-seq Data. bioRxiv [Preprint]. 2024 Aug 16:2024.08.13.607784. doi: 10.1101/2024.08.13.607784. PMID: 39185246; PMCID: PMC11343128.

License: This is an open access protocol distributed under the terms of the [Creative Commons Attribution License](#), which permits unrestricted use, distribution, and reproduction in any medium, provided the original author and source are credited

Protocol status: Working

We use this protocol and it's working

Created: August 31, 2024

Last Modified: September 10, 2024

Protocol Integer ID: 106781

Keywords: R2C2, single cell RNA-seq, ONT, long-read sequencing, cdna this protocol, r2c2 protocol, using cdna, rna, rolling circle to concatemeric consensus, single cell rna, seq assay, cell rna, 10x splint fragment, cdna, r2c2, sequencing, nanopore, sequencing platform, oxford nanopore technology, concatemeric consensus, circle amplification, rolling circle amplification, major modification from the original protocol, seq assay cdna this protocol, seq assay cdna

Funders Acknowledgements:

Aligning Science Across Parkinson's

Grant ID: ASAP-000301

Abstract

This protocol is based on the published Rolling Circle to Concatemeric Consensus (R2C2) method (1). The aim is to make elongated cDNA via rolling circle amplification using cDNA generated from the 10x Chromium 3' (or 5') single-cell RNA-seq assay to increase the accuracy of downstream sequencing on an Oxford Nanopore Technologies sequencing platform. The major modification from the original protocol is to add an additional cDNA re-amplification step (see section "10x cDNA re-amplification") to enrich for fragments with the proper sequence (poly(T) at 3' and TSO at 5') at both ends, which is essential in circularization with the 10x Splint fragment.

Image Attribution

Xian Adiconis

Materials

Oligos

	A	B	C
	Oligo Name	Vendor	Sequence
	10X_UMI_Splint_Forward	IDT	5'- AGATCGGAAGAGCGTCGTGTAGTGAGGCTGATGAGTTCCATA NNNNNTATATNNNNNATCACTACTTAGTTTTTTGATAGCTTC AAGCCAGAGTTGTCTTTTTCTCTTTGCTGGCAGTAAAAG -3'
	10X_UMI_Splint_Reverse	IDT	5'- CTCTGCGTTGATACCACTGCTTAAAGGGATATTTTCGATCGC NNNNNATATANNNNNTTAGTGCAATTTGATCCTTTTACTCCTC CTAAAGAACAACCTGACCCAGCAAAAGGTACACAATACTTTT ACTGCCAGCAAAGAG -3'
	PCR_primer1	IDT	5'- CTACACGACGCTCTTCCGATCT -3'
	PCR_primer2	IDT	5'- AAGCAGTGGTATCAACGCAGAGT -3'

Reagents

	A	B	C
	Reagents	Vendor	Part Number
	KAPA HiFi HotStart ReadyMix	KAPA Bioscience	KK2602
	NEBNext® High-Fidelity 2X PCR Master Mix	New England Biolabs	M0541S
	Select-a-Size DNA Clean & Concentrator Kit	Zymo Research	D4080
	ProNex® Size-Selective Purification System, 10ml	Promega	NG2001
	NEBuilder® HiFi DNA Assembly Master Mix	New England Biolabs	E2621S
	NEB Buffer 2	New England Biolabs	B7002S
	Exonuclease I (20 U/μl)	New England Biolabs	M0293S
	Exonuclease III (100 U/μl)	New England Biolabs	M0206S
	Lambda Exonuclease (5 U/μl)	New England Biolabs	M0262S
	Agencourt AMPure® XP SPRI beads, 60 ml	Beckman Coulter Genomics	A63881
	phi29 DNA Polymerase (10 U/μl)	New England Biolabs	M0269S

	A	B	C
	Deoxynucleotide (dNTP) Solution Mix (10mM)	New England Biolabs	N0447L
	Exo-Resistant Random Primer (10 μ M)	Thermo Fisher Scientific	SO181
	T7 Endonuclease I (10 U/ μ l)	New England Biolabs	M0302S

Safety warnings

 For hazard information and safety warnings, please refer to the MSDSs (Material Safety Data Sheets).



Splint generation

2h

- 1 Prepare the following on ice

10m

A	B
	1X (μl)
KAPA HiFi HotStart ReadyMix	25
H2O	23
10x_UMI_Splint_forward (100 μM)	1
10x_UMI_Splint_reverse (100 μM)	1
Total	50

- 2 Run the following program

11m

Thermocycler Conditions: (Lid@ 105 °C , 50 μl)

95 °C , 00:03:00

followed by 1 cycle of:

98 °C , 00:01:00

62 °C , 00:01:00

72 °C , 00:06:00

then

4 °C , 00:00:00 hold

- 3 Cleanup with Select-a-Size columns (Zymo) and a size cut-off of ~125 bp. Follow the Select-a-Size cleanup protocol for single-size selection, but in the buffer preparation add 85 μL of 100% ethanol to 500 μL of Select-a-Size DNA binding buffer. Elute in 15 μL DNA elution buffer (from the kit). QC with 1/5x dilution for Quant-it (Thermo Fisher Scientific) and 1/50x dilution for BioAnalyzer DNA HS assay (Agilent).

1h 30m

Label this product as "10x Splint", store at -20 °C for later use.

10x cDNA re-amplification

2h

4 Enrichment PCR

10m

Prepare the follow reaction mix on ice

A	B
	1X (μ l)
NEBNext® High-Fidelity 2X PCR Master Mix	50
12 μ M PCR Primer 1	2
12 μ M PCR Primer 2	2
20 ng 10x cDNA in H2O	46
Total	100

5 Run the following program

9m 10s

Thermocycler Conditions: (Lid@ 105 °C , 100 μ l)

98 °C , 00:00:45

followed by 10 cycle of:

98 °C , 00:00:10

62 °C , 00:00:15

72 °C , 00:03:00

then

72 °C , 00:05:00

4 °C , 00:00:00 hold

6 cDNA purification

1h 30m

- 6.1 Add 95 μ L of resuspended, room-temperature ProNex beads to the PCR mix.
Pipette mix 10 times. Perform a quick spin to collect all liquid from the sides of the tube.

- 6.2 Incubate sample on bench top for 00:05:00 at Room temperature .

5m

- 6.3 Place the tube on a magnetic stand to separate the beads from the supernatant. Use a P200 pipettor to remove the supernatant.
- 6.4 Wash 2 times with  200 μL of freshly prepared 80% ethanol. After removal of the second wash of  200 μL of ethanol, spin the tube strip briefly, return to magnetic stand and remove residual ethanol with a P20 pipettor. Do not let the beads dry out.
- 6.5
 - Remove the tube from the magnetic stand. Immediately add  15 μL of buffer EB (10 mM Tris-HCl, pH 8.5) and pipette mix to resuspend. Perform a quick spin to collect all liquid from the sides of the tube. Place at  Room temperature for  00:05:00 to elute the DNA from the beads.
- 6.6 Place the tube on a magnetic stand to separate the beads from the supernatant. Transfer the eluted cDNA samples to a new tube.
- 6.7 QC with 1/5x dilution for Quant-it and BioAnalyzer DNA HS assay. Store the remaining product at  -20 $^{\circ}\text{C}$ for later use.

5m

cDNA circularization

8h

7 Circularization

7h 20m

Prepare the following mix

A	B
	1X (μl)
10x Splint (200 ng)	2
Re-amplified cDNA (200 ng)	8
NEBuilder® HiFi DNA Assembly Master Mix	10
Total	20

Incubate @  50 $^{\circ}\text{C}$ for  01:00:00

Then add the following mix

A	B
	1X (μl)



A	B
NEB Buffer 2	5
H2O	16
Exonuclease I (20 U/μl)	3
Exonuclease III (100 U/μl)	3
Lambda Exonuclease (5 U/μl)	3
Total	30

Incubate @ 37 °C for 06:00:00 , then @ 80 °C for 00:20:00

Cleanup the reaction with 40 μL SPRI beads (0.8x)

8 Circularization product purification

30m

8.1 Add 40 μL (0.8x) of resuspended, room-temperature AMPure XP SPRI beads to the reaction mix. Pipette mix 10 times. Perform a quick spin to collect all liquid from the sides of the tube.

8.2 Incubate sample on bench top for 00:05:00 at Room temperature .

5m

8.3 Place the tube on a magnetic stand to separate the beads from the supernatant. Use a P200 pipettor to remove the supernatant.

8.4 Wash 2 times with 200 μL of freshly prepared 80% ethanol. After removal of the second wash of 200 μL of ethanol, spin the tube strip briefly, return to magnetic stand and remove residual ethanol with a P20 pipettor. Do not let the beads dry out.

8.5 Remove the tube from the magnetic stand. Immediately add 30 μL of H₂O and pipette mix to resuspend. Perform a quick spin to collect all liquid from the sides of the tube. Place at Room temperature for 00:05:00 to elute the cDNA from the beads.

5m

8.6 Place the tube on a magnetic stand to separate the beads from the supernatant. Transfer the eluted cDNA samples to a new tube.

Rolling circle amplification

16h

- 9 Prepare the following mix, then split into 3 reaction wells of  50 µL each.

A	B
	1X (µl)
phi29 Buffer (10x)	15
phi29 DNA Polymerase (10 U/µl)	3
dNTP (10 mM)	7.5
Exo-Resistant Random Primer (10 µM)	7.5
Circularized cDNA	30
H2O	87
Total	150

Incubate in a thermocycler @  30 °C for  Overnight

- 9.1 From this point on, be really careful with the DNA since it will shear pretty easily. Do not pipet up and down and try not to vortex. 

10 SPRI clean and T7 endonuclease treatment

2h

Pool all reactions and adjust volume to  300 µL with H₂O. Add  150 µL

 150 µL SPRI beads (0.5x) to your DNA and leave on a rotator for 5 minutes to mix. If not homogeneous, flick the tube to mix.

Notes: After adding the AMPure XP SPRI beads and mixing by inverting the tubes, there's uneven distribution of the beads, mostly the sign of longer cDNA strand.

Place on a magnet stand until a complete pellet forms, wash twice with  500 µL 70% ethanol. After the ethanol removal, add the following mix to re-suspend the beads.


A	B
	1X (µl)
NEB Buffer 2	10
T7 Endonuclease I (10 U/µl)	5

A	B
H ₂ O	90
Total	105


Incubate on a thermal shaker at 37°C at 1000 RPM for 02:00:00 . Agitate the tube occasionally to help the solution homogenize.

Place on a magnet stand to pellet the beads. Recover the supernatant that contains the DNA to a new tube. Purify the DNA with 0.5x volume of AMPure XP SPRI beads, wash twice with 500 μL 70% ethanol. Elute in 20 μL H₂O.

QC with Quant-it for concentration, run BioAnalyzer DNA 12000 assay (Agilent) or 1% agarose gel to visualize the size distribution. The majority of the fragments should be > 10kb.

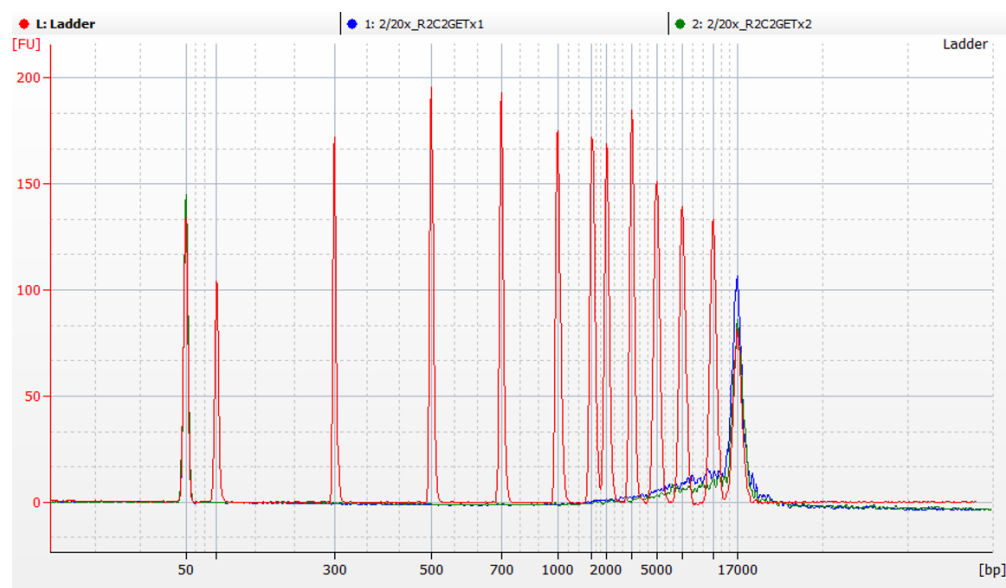

Rolling circle amplified DNA trace with BioAnalyzer DNA 12000 assay

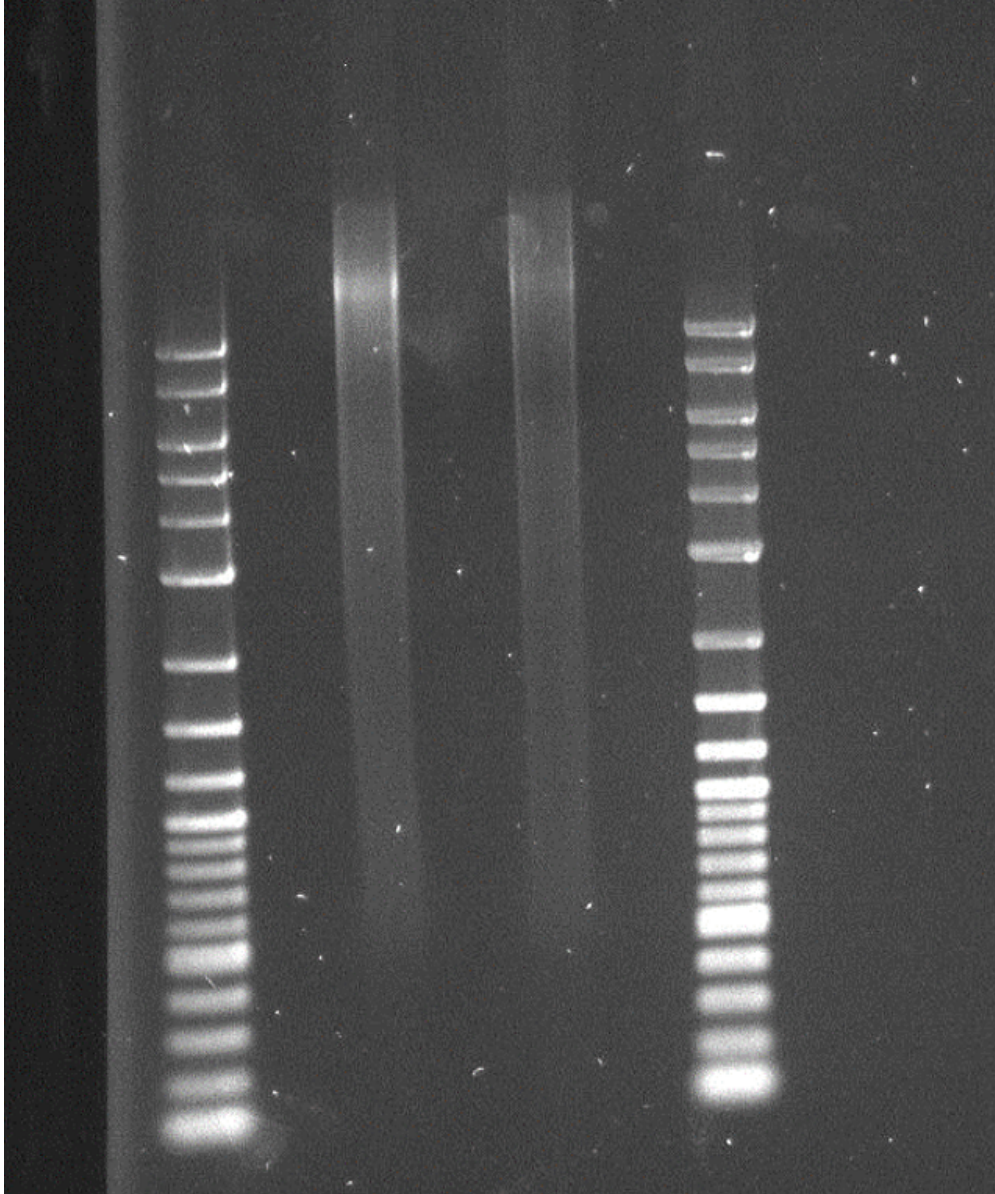

Rolling circle amplified DNA on a 1% agarose gel; the top band of the ladder is 10kb.

The elongated cDNA is now ready for downstream Oxford Nanopore Technologies (ONT) library prep.

Protocol references

1. Volden R, Vollmers C. Single-cell isoform analysis in human immune cells. *Genome Biol.* 2022;23(1):47.