


Mar 10, 2024 Version 2

Rapid transposase based total DNA library preparation and sequencing using MinION V.2

DOI

dx.doi.org/10.17504/protocols.io.j8nlko796v5r/v2

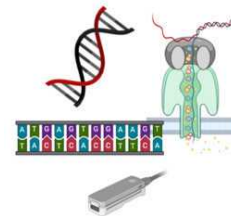

Lavi Singh¹, Ashley Jones¹, Benjamin Schwessinger¹

¹Australian National University



Benjamin Schwessinger

Australian National University

Create & collaborate more with a free account

Edit and publish protocols, collaborate in communities, share insights through comments, and track progress with run records.

[Create free account](#)

OPEN  ACCESS



DOI: <https://dx.doi.org/10.17504/protocols.io.j8nlko796v5r/v2>

Protocol Citation: Lavi Singh, Ashley Jones, Benjamin Schwessinger 2024. Rapid transposase based total DNA library preparation and sequencing using MinION. **protocols.io** <https://dx.doi.org/10.17504/protocols.io.j8nlko796v5r/v2> Version created by **Benjamin Schwessinger**

License: This is an open access protocol distributed under the terms of the **Creative Commons Attribution License**, which permits unrestricted use, distribution, and reproduction in any medium, provided the original author and source are credited



Protocol status: Working

We use this protocol and it's working

Created: March 10, 2024

Last Modified: March 10, 2024

Protocol Integer ID: 96431

Keywords: different wheat pathogen, wheat pathogen, uninfected wheat plant, plant pathogen diagnostic, week old wheat plant, major pathogen, old wheat plant, total dna library preparation, sequencing adapter, pathogen, based total dna library preparation, sequencing, native dna sample, genomic, sequencing run, infective agent for each treatment group, dna sample, total native dna, specific dna barcode, rapid transposase, dna samples to same concentration

Abstract

This protocols is part of the ANU Biosecurity mini-research project #1 "Plant Pathogen Diagnostics: Visuals, subcultures, and genomics".

You will be provided four pots of 3-4 week old wheat plants that have been infected with different wheat pathogens. Each pot has been infected with one major pathogen. You will not know which pot has been infected with which pathogen. However, you will be provided a compendium of 10-15 wheat pathogens that will guide you to identify the infective agent for each treatment group. The fifth treatment group will be uninfected wheat plants which will be clearly identified. You can use treatment group #5 as negative control for your experiments.

In total, each group will obtain five pots each:


	A	B
	Treatment group 1	Unknown infective agent
	Treatment group 2	Unknown infective agent
	Treatment group 3	Unknown infective agent
	Treatment group 4	Unknown infective agent
	Treatment group 5	Uninfected control

This specific protocol describes the molecular biology and a step-by-step guide for rapid barcoding of total native DNA extracted from the five treatment groups followed by Nanopore sequencing. We will perform sequencing of all five native DNA samples from three research groups to enable us to have three independent replicates per sample. Replications make our observation and analyses more robust. Total native DNA contains all the DNA extracted from your leaf material plus any contamination that was part of the extraction process. We did not include a negative extraction control this time round.

We will select the samples of three research groups based on their quality control values such as DNA concentration and DNA size measured by gel electrophoresis. Most of the class need not be present for these library preparations [Step 1-3] but will join in for loading the final library and starting sequencing runs. We will also explain more about the theory during the lab.

This protocol is applicable for week 6.

Conceptual overview:

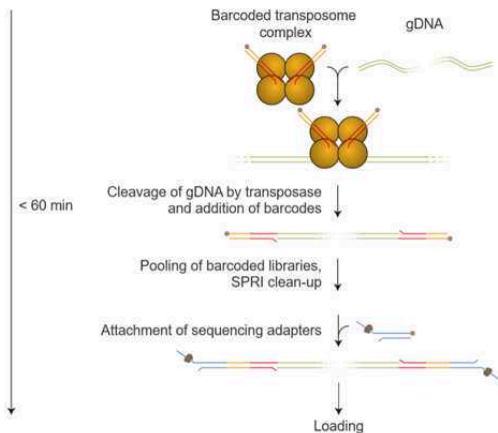

Adapted from Nanopore Technologies

- Bring all DNA samples to same concentration and same volume.
- Barcode individual samples with transposase loaded with specific DNA barcodes.
- Pool all samples and clean up with magnetic beads.
- Add sequencing adapters with click chemistry applied to the barcoded and pooled DNA samples.
- Prepare for loading onto the flowcell.
- Prime flowcell to make it ready for loading.
- Load library.
- Start of sequencing run.
- Basecall during or after the sequencing run.

You can cite this protocol in the methods section of your report as for all other protocols. No need to write it all up again :).

Attachments



Rapid sequencing

DNA...

96KB



rapid-sequencing-gdn...

1.5MB

Image Attribution

The logo image was generated with biorender.



Guidelines

This protocol describes how to carry out native barcoding of total native DNA using the Rapid Barcoding Kit 96 (SQK-RBK114.96). There are 96 unique barcodes available, allowing the user to pool up to 96 different samples in one sequencing experiment. We will use it to sequence 15 samples as once.

Rapid Barcoding Kit features. This kit is recommended for users who:

- wish to multiplex up to 96 samples to reduce sequencing cost per sample
- need a PCR-free method of multiplexing
- want to rapidly prepare libraries without aiming for ultralong reads

As always you need to bring a lab notebook, a printed version of this protocol, and a pen to record your adventures in the lab.

You must have read, understood, and follow the health and safety instructions provided in the "Overview Mini-Research Project #1 BIOL3106/6106" provided on Wattle (ANU learning portal).

You must have signed and returned one copy of the "Student Safety Declaration Form For Practical Class Work" before starting any laboratory work.

You must have read and understood the Hazard Sheets (Risk assessment) of all chemicals listed bellow in the "Safety Warnings" section. These Hazard Sheets are provided on Wattle as part of the "Overview Mini-Research Project #1 BIOL3106/6106" document.



Materials

Materials

- Pure DNA

Consumables

- R10.4.1 MinION flowcell **Flow Cell (R10.4.1) (nanoporetech.com)**
- SQK-RBK114.96 **Rapid Barcoding Kit 96 V14 (nanoporetech.com)**
- magnetic beads (AMPure or equivalent) <https://www.beckman.com/reagents/genomic/cleanup-and-size-selection/pcr>
- 1.5 ml Eppendorf DNA LoBind tubes
- 0.2 ml thin-walled PCR tubes
- Nuclease-free water (e.g. ThermoFisher)
- Freshly prepared 80% ethanol in nuclease-free water
- Qubit Assay Tubes (ThermoFisher)
- Qubit dsDNA BR Assay Kit (ThermoFisher)

Equipment

- Hula mixer (gentle rotator mixer)
- Microfuge Magnetic rack
- Vortex mixer
- Thermal cycler/PCR machine
- Multichannel pipette
- P1000 pipette and tips
- P200 pipette and tips
- P100 pipette and tips
- P20 pipette and tips
- P10 pipette and tips
- P2 pipette and tips
- Ice bucket with ice Timer
- Qubit fluorometer (or equivalent for QC check)
- Computer with the latest version of MinKNOW installed.
- MinION Nanopore sequencer

Troubleshooting



Safety warnings

! This protocol does not require any hazardous substances.

You need to wear safety equipment at all times including lab coats, gloves, and safety goggles when handling chemicals and biological agents. While the major biological agents used for the wheat infection are pathogens commonly found in Australia, you must treat them as they were infective agents of general concern. Treat them with care. Do not remove them from the laboratory. Do not spread them via clothing. Use a dedicated notebook and pen to make notes during the mini-research project. Do not put anything into your mouth while in the laboratory. Wash your hands each time you leave the laboratory.

Preparation of input DNA

- 1 This first series of steps [1.x] will be performed by the three groups we selected in week 5. We will use their total native DNA samples for sequencing. We selected these samples bases on DNA concentrations and DNA integrity. We assessed the former using Qubit and the latter agarose gel electrophoresis
 - 1.1 Retrieve the DNA concentrations of your five samples from your Qubit measurements in week 4.

To note these are the original DNA samples and their concentrations BEFORE dilution to 10ng/ul.
 - 1.2 Label a PCR strip tube of eight with your group number and treatment group sample numbers.
 - 1.3 For each sample prepare 500 ng as input for the library preparation in 10ul. This means calculate the volume in uL you need for each sample to reach 500 ng given your concentration measured in week 4. Pipette this volume into the labelled tube and top it up to 10 ul with nuclease free water.

Remember:

$n = c \cdot V$ with n being the amount in ng, c the concentration in ng/ul, and V the volume in ul.

Consult previous protocols if you struggle with these calculations and check with a demonstrator.
 - 1.4 Mix the samples by short pipetting up and down. Collect the samples at the bottom of the tubes using a benchtop mini-centrifuge.
- 2 Once all three groups are done we will have 15 samples in total that we can process for library preparation. These include three replicates for each treatment group.

Rapid barcoding library preparation

- 3 For these steps we can rotate between groups and people. Only one person will perform the (sub) step at any one point in time.
 - 3.1 Program the thermal cycler: 30°C for 2 minutes, then 80°C for 2 minutes. Ramp it up and keep it at 30°C so we can start the program immediately.



3.2 Note down which sample will receive which barcode.

3.3 To each sample add one rapid barcode.

Pipette 1.5 ul of the barcode into the 10 ul of 500 ng of DNA.

Immediately mix all the components by pipetting up and down.

Collect all liquid at the bottom of the tube via a short centrifuge spin.

We will add the barcodes with a mutli-channel pipette and the demonstrator will help you with it or perform the task for you.

3.4 Incubate the tubes at 30°C for 2 minutes and then at 80°C for 2 minutes using the PCR machine we already set up. Once done, briefly put the tubes on ice to cool.

3.5 Pool all barcoded samples in a clean 1.5 ml Eppendorf DNA LoBind tube, noting the total volume of $15 \times 11.5\text{uL} = 172.5\text{ul}$

3.6 Resuspend the magentic beads for DNA binding by vortexing.

3.7 Add an equal volume of magnetic beads to the pooled barcoded DNA samples.

172.5ul DNA + 172.5ul magnetic beads.

3.8 Incubate while rotating the tube with your hand for 10 minutes at room temperature

3.9 Spin down the sample and pellet on a magnet. Keep the tube on the magnet, and pipette off the supernatant.

A demonstrator will help you with this step.

3.10 Keep the tube on the magnet and wash the beads with 1 ml of freshly prepared 80% ethanol without disturbing the pellet. Remove the ethanol using a pipette and discard.

3.11 Repeat the previous step 3.10.

- 3.12 Briefly spin down and place the tube back on the magnet. Pipette off any residual ethanol. Allow to dry for 1 - 2 mins, but do not dry the pellet to the point of cracking
- 3.13 Remove the tube from the magnetic rack and resuspend the pellet in 15 µl Elution Buffer (EB).
- 3.14 Incubate for 10 minutes while warming the tube with your fingers.
- 3.15 Pellet the beads on a magnet until the eluate is clear and colourless, for at least 1 minute.
- 3.16 Remove and retain the full volume of eluate into a clean 200ul PCR tube.
- 3.17 Quantify 1ul of library using Qubit BR as described in previous protocols. Note down the concentration
- 3.18 Transfer 11 µl of the sample into a clean 200ul PCR tube.
- 3.19 Add 0.5uL of Rapid Adapter (RA) and 1.2uL of Adapter Buffer (ADB) to the DNA.

A demonstrator will help you with this step.
- 3.20 Mix gently by flicking the tube, and spin down
- 3.21 Incubate the reaction for 5 minutes at room temperature.
- 3.22 Add 37.6uL of Sequencing Buffer (SB) and 25.5uL of Library Beads (LIB) to the sample.

Note – the library beads settle on the bottom of the tube very quickly. Make sure to mix well before pipetting.
- 3.23 We now have our library ready to load and will wait for the rest of the class to come in.

Flowcell preparation for sequencing aka priming



- 4 These steps will be performed by one group member while the rest of the class is watching.

If you are not familiar with MinION flowcells check out this explanation video here [Community - Flow cell introduction \(nanoporetech.com\)](#).


This is a useful video for flowcell priming and loading [Community - Priming and loading your flow cell \(nanoporetech.com\)](#).

- 4.1 Perform a flowcell check as shown here [Community - MinKNOW: Flow cell check \(nanoporetech.com\)](#).
- 4.2 To prepare the flow cell priming mix with BSA, combine the following reagents in a fresh 1.5 ml Eppendorf DNA LoBind tube. Mix by inverting the tube and pipette mix at room temperature.

Reagents	Volume [ul]
Flow Cell Flush (FCF)	1170
Bovine Serum Albumin (BSA) at 50 mg/ml	5
Flow Cell Tether (FCT)	30
Total volume	1205


- 4.3 Open the MinION device lid and slide the flow cell under the clip. Press down firmly on the flow cell to ensure correct thermal and electrical contact. This step might have been already performed if you did the flow cell check.

1a
Insert the flow cell into the device under the clip and press down firmly.

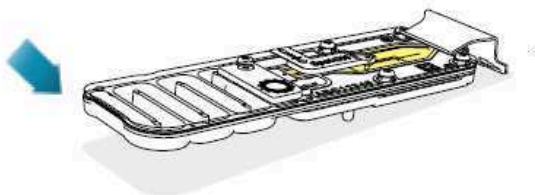


Key:
● Storage buffer
● Priming mix
● DNA library


1b
Insert the flow cell into the device under the clip and press down firmly.

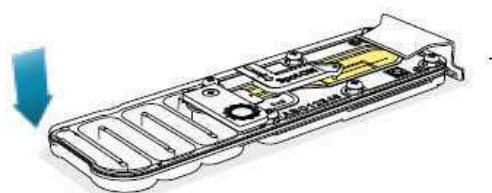

- 4.4 Slide the flow cell priming port cover clockwise to open the priming port. The priming port cover has the Nanopore logo.

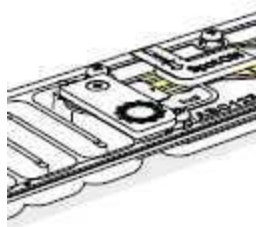

Closed priming port cover.

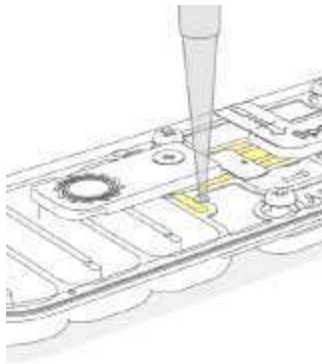

Open priming port cover with pipette entering the priming port.

4.5 IMPORTANT


Take care when drawing back buffer from the flow cell. Do not remove more than 20-30 μl , and make sure that the array of pores are covered by buffer at all times. Introducing air bubbles into the array can irreversibly damage pores.

4.6 After opening the priming port, check for a small air bubble under the cover. Draw back a small volume to remove any bubbles:

1. Set a P1000 pipette to 200 μl
2. Insert the tip into the priming port
3. Turn the wheel until the dial shows 220-230 μl , to draw back 20-30 μl , or until you can see a small volume of buffer entering the pipette tip.


Note: Visually check that there is continuous buffer from the priming port across the sensor array.

3 Insert a P1000 pipette with an empty tip into the **Priming port**. Turn the pipette wheel to draw back 20–30 μl or until you can see a small volume of buffer entering the pipette tip.

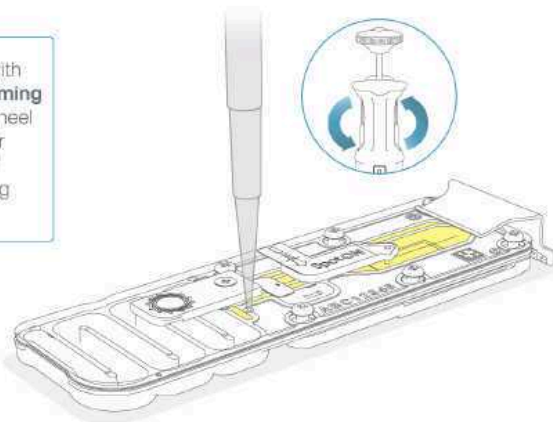

- 4.7** Load 800 μl of the priming mix into the flow cell via the priming port, avoiding the introduction of air bubbles. This is best done by turning the wheel slowly and ejecting the priming buffer slowly into the priming port. Close the priming port cover during incubation. Wait for five minutes.

4 Slowly load 800 μl of the priming mix into the **Priming port**. Ensure there are no air bubbles in the pipette tip.

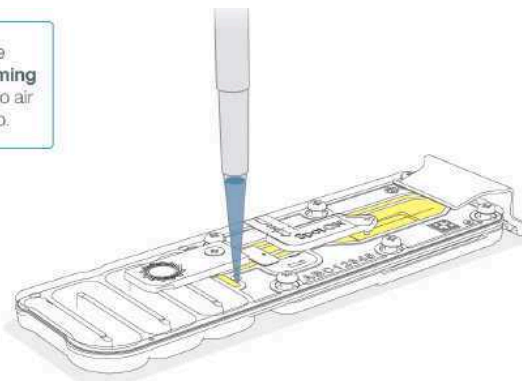

Wait 5 minutes before proceeding to the next step.

- 4.8** The flowcell is now ready for loading and sequencing.

Loading of DNA library onto MinION flowcell

- 5** Now we will load the DNA library from step 3.23 to start sequencing.


This is a useful video for flowcell priming and loading [Community - Priming and loading your flow cell \(nanoporetech.com\)](#).

5.1 IMPORTANT

The Library Beads (LIB) tube contains a suspension of beads. These beads settle very quickly. It is vital that they are mixed immediately before use.

5.2 Complete the flow cell priming:


1. priming port cover.
2. Gently lift the SpotON sample port cover to make the SpotON sample port accessible.
3. Load 200 μ l of the priming mix into the flow cell priming port (not the SpotON sample port), avoiding the introduction of air bubbles.

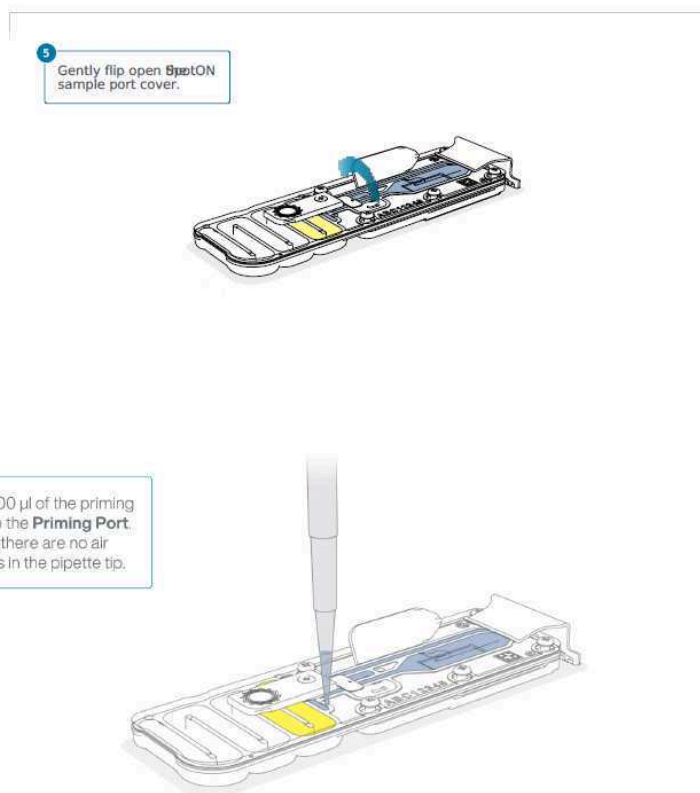

5.3 Mix the prepared library gently by pipetting up and down just prior to loading.


- 5.4 Add 75 μ l of the prepared library to the flow cell via the SpotON sample port in a dropwise fashion. Ensure each drop flows into the port before adding the next.

A demonstrator will assist you with this step.

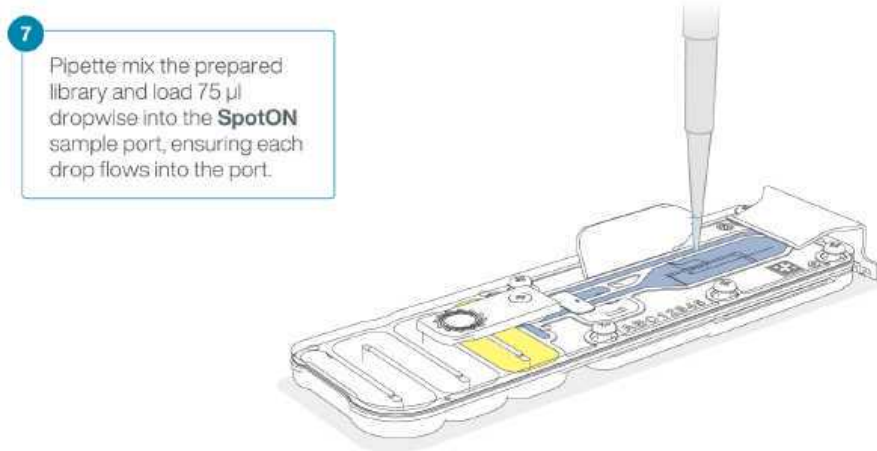


- 5.5 Gently replace the SpotON sample port cover, making sure the bung enters the SpotON port and close the priming port.

8 Gently replace the **SpotON** sample port cover.

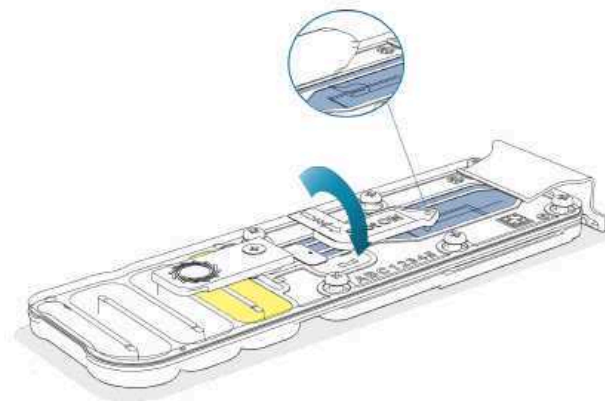

9 Gently close the **Priming port**.

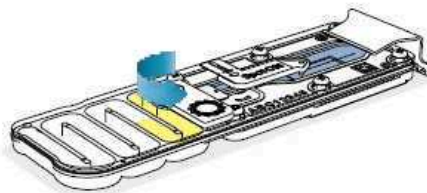

5.6 IMPORTANT

Install the light shield on your flow cell as soon as library has been loaded for optimal sequencing output.

We recommend leaving the light shield on the flow cell when library is loaded, including during any washing and reloading steps.

The shield can be removed when the library has been removed from the flow cell. Gently replace the SpotON sample port cover, making sure the bung enters the SpotON port and close the priming port.



5.7 Place the light shield onto the flow cell, as follows:

1. Carefully place the leading edge of the light shield against the clip.

Note: Do not force the light shield underneath the clip.



2. Gently lower the light shield onto the flow cell. The light shield should sit around the SpotON cover, covering the entire top section of the flow cell.

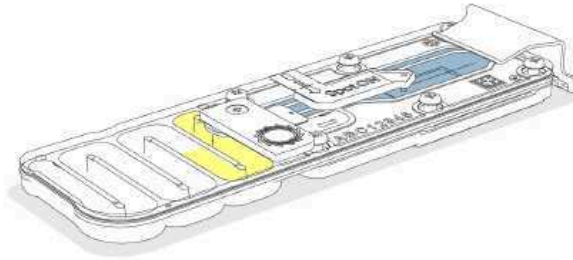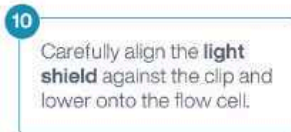

- 5.8 Now we are ready for DNA sequencing via the MinKNOW graphical user interface.

DNA sequencing and basecalling

- 6 The demonstrators will help setting up the sequencing runs. We will perform basecalling later on.

Here is a list of useful videos:

- Introduction to the MinKNOW graphic user interface [Community - MinKNOW: The graphical user interface \(GUI\) \(nanoporetech.com\)](#)
- How to set up a sequencing experiment [Community - MinKNOW: Setting up a new experiment \(nanoporetech.com\)](#)
- How to assess your run is going fine [Community - MinKNOW: Assessing your run \(nanoporetech.com\)](#)
- How MinKNOW selects the best channels [Community - MinKNOW: Mux scan and active channel selection \(nanoporetech.com\)](#)
- More about checking on if your sequencing run is going smoothly [Community - MinKNOW: Pore occupancy \(nanoporetech.com\)](#)

- 6.1 We will share sequencing results and initial analysis beginning of week 7.



Protocol references

This protocol is based on publicly available protocols from Oxford Nanopore Technology found here [Community - Rapid sequencing DNA V14 - barcoding \(SQK-RBK114.24 or SQK-RBK114.96\) Protocol - Rapid sequencing DNA V14 - barcoding \(SQK-RBK114.24 or SQK-RBK114.96\) \(nanoporetech.com\)](#).