

Oct 24, 2020

Restriction Digest

DOI

dx.doi.org/10.17504/protocols.io.bnwvmfe6

Zhujun Wei¹

¹2020 iGEM NEFU China

2020 iGEM NEFU China



Zhujun Wei

OPEN  ACCESS



DOI: dx.doi.org/10.17504/protocols.io.bnwvmfe6

Protocol Citation: Zhujun Wei 2020. Restriction Digest. **protocols.io** <https://dx.doi.org/10.17504/protocols.io.bnwvmfe6>

License: This is an open access protocol distributed under the terms of the **Creative Commons Attribution License**, which permits unrestricted use, distribution, and reproduction in any medium, provided the original author and source are credited

Protocol status: Working

We use this protocol and it's working

Created: October 24, 2020

Last Modified: October 24, 2020

Protocol Integer ID: 43701

Abstract


This protocol includes three different methods of digestion.

Materials

Restriction enzyme

Buffer 2.1/3.1/CutSmart

Safety warnings

 Please wear gloves for the experiment.



- 1 Add the following reagents to a tube.

Total	10ul	30ul	50ul
DNA	0.4ug	1.2ug	2ug
Enzyme	0.2ul	0.6ul	1ul
Buffer(10*)	1ul	3ul	5ul
ddwater Fill the rest with water			

- 2 Pipette up and down thoroughly.
- 3 Put at 37°C for about 2 hours (according to different enzymes).
- 4 Add the following reagents to a tube.

Total	10ul	30ul	50ul
DNA	0.4ug	1.2ug	2ug
Enzyme	0.2ul	0.6ul	1ul
Buffer(10*)	1ul	3ul	5ul
ddWater	Fill the rest with water	Fill the rest with water	Fill the rest with water

- 5 Pipette up and down thoroughly.
- 6 Put at 37°C for about 2 hours (according to different enzymes).
- 7 Put at 65°C for about 20 min.(according to different enzymes)
- 8 Purification of enzyme digestion products with Cycle Pure Kit D6492 (OMEGA bio-tek)



9 DNA product concentrations were measured using Nano

10 Add the following reagents to a tube:

Total	10ul	30ul	50ul
DNA product	0.4ug	1.2ug	2ug
CIP(NEB)	0.2ul	0.6ul	1ul
Cutsmart(10*)	1ul	3ul	5ul
ddWater	Fill the rest with water	Fill the rest with water	Fill the rest with water

11 Pipette up and down thoroughly.

12 Put at 37°C for about 2 hours (according to different enzymes).

13 Add the following reagents to a tube.

Total	10ul	30ul	50ul
DNA	0.4ug	1.2ug	2ug
Enzyme 1	0.2ul	0.6ul	1ul
Enzyme 2	0.2ul	0.6ul	1ul
Buffer(10*)	1ul	3ul	5ul
ddwater	Fill the rest with water	Fill the rest with water	Fill the rest with water

14 Pipette up and down thoroughly.

15 Put at 37°C for about 2 hours (according to different enzymes).



16 Add the following reagents to a tube.

Total	10ul	30ul	50ul
DNA product	0.4ug	1.2ug	2ug
DpnI(NEB)	0.2ul	0.6ul	1ul
Cutsmart(10*)	1ul	3ul	5ul
ddwater	Fill the rest with water	Fill the rest with water	Fill the rest with water

17 Pipette up and down thoroughly.

18 Put at 37°C for about 2 hours (according to different enzymes).