

Mar 30, 2019

RNAi by feeding in *Euplotes focardii* (povisional)

 In 1 collection

DOI

dx.doi.org/10.17504/protocols.io.zk8f4zw

Francesca Papi¹, Angela Piersanti¹

¹University of Camerino

Protist Research to Opti...



Angela Piersanti

University of Camerino

OPEN  ACCESS



DOI: dx.doi.org/10.17504/protocols.io.zk8f4zw

Protocol Citation: Francesca Papi, Angela Piersanti 2019. RNAi by feeding in *Euplotes focardii* (povisional). **protocols.io**
<https://dx.doi.org/10.17504/protocols.io.zk8f4zw>

License: This is an open access protocol distributed under the terms of the **Creative Commons Attribution License**, which permits unrestricted use, distribution, and reproduction in any medium, provided the original author and source are credited

Protocol status: In development

We are still developing and optimizing this protocol

Created: March 29, 2019

Last Modified: March 30, 2019

Protocol Integer ID: 21888

- 1 Grow RNase III deficient *E.coli* strain HT115 in LB with antibiotic selection o/n at 37°C.
- 2 Prepare a 1:100 dilution of the bacterial culture, and grow it at 37°C until it reaches an OD600 of 0.4.
- 3 Add 0.4 mM IPTG, and induce RNA transcription from the L4440 plasmid, in which it is included the gene of interest to be silenced, in the bacteria at 37°C, until they reach a OD600 of 1.
Bacteria with L4440 plasmid, empty vector (EV), were grown as well to be used as control.
- 4 Collect the bacteria by centrifugation at 7000 rpm for 8 minutes, wash them twice with ddH₂O, then resuspend in 1.5 ml of sea water (at 4°C) to an OD600 of 4 to feed 100 ml of *Euplotes focardii* culture with a cell density of ~3000 cells/ml.
- 5 Feed *Euplotes focardii* culture every 3 days, at least for 10 days and checked the phenotype.