

Apr 06, 2024

Version 2

Rotarod (40 RPM and 20 RPM with docking) V.2

DOI

<https://dx.doi.org/10.17504/protocols.io.3byl4qo5zvo5/v2>

daniel.dautan daniel^{1,2}, Per Svenningsson^{1,2}

¹Department of Clinical Neuroscience, Karolinska Institutet, 171 76 Stockholm, Sweden;

²Aligning Science Across Parkinson's (ASAP) Collaborative Research Network, Chevy Chase, MD 20815, USA



Jacquelyn Haytayan

Weill Cornell Medicine

Create & collaborate more with a free account

Edit and publish protocols, collaborate in communities, share insights through comments, and track progress with run records.

Create free account

OPEN  ACCESS



DOI: <https://dx.doi.org/10.17504/protocols.io.3byl4qo5zvo5/v2>

Protocol Citation: daniel.dautan daniel, Per Svenningsson 2024. Rotarod (40 RPM and 20 RPM with docking) . protocols.io <https://dx.doi.org/10.17504/protocols.io.3byl4qo5zvo5/v2> Version created by [Eileen Ruth Torres](#)

License: This is an open access protocol distributed under the terms of the **Creative Commons Attribution License**, which permits unrestricted use, distribution, and reproduction in any medium, provided the original author and source are credited

Protocol status: Working

We use this protocol and it's working



Created: February 23, 2024

Last Modified: September 23, 2024

Protocol Integer ID: 95631

Keywords: ASAPCRN, rpm with docking, rpm, motor function testing in mice, rotarod by the svenningsson lab, rotarod, motor function testing, docking, rpm constant speed after habituation, constant speed, mice, motor function

Funders Acknowledgements:

Aligning Science Across Parkinson's

Grant ID: 020608

Abstract

This protocol describes motor function testing in mice using the rotarod by the Svenningsson lab.

Two options for the protocol are included: 1) 40 RPM constant speed after habituation and 2) 20 RPM with docking.

Materials

Mouse Rotarod (Ugo Basile 47650)

50% ethanol



- 1 Transfer the cage into the experimental room at least 1 hr before testing.
- 2 To avoid stress, always put the cage into a ventilated cage rack.
- 3 Using a paper tube, transfer each mouse of the same cage onto the rotarod (Ugo Basile, 47650). Always maintain order, mice one into line one, mice 2 into line 2 etc...
- 4 Wait 2-3 minutes for the mice to habituate to the apparatus.
- 5 Lift the collecting platform and set up the protocol: either max speed of constant 40 RPM or 20 RPM docking.

STEP CASE

40 RPM 4 steps

- 6 Following habituation, start the rotarod program to reach a speed of 40 RPM within the first 5s of the protocol.
- 7 The time to fall is defined as the moment the mice fall in the receptacle or the moment the mice turn around the rotor more than 5 consecutive times. The maximum trial length is 2 min.
- 8 Repeat the experiment 3 times with at least 6 hr between each trial. If an animal does not fall or turn more than 5 times within 120s of the trial, stop the trial and record "120" as the individual time.
- 9 Analyze for each mouse, the time to fall for each session, the average and and max time.