

Oct 21, 2019

RPA DNA Amplification using Agdia AmplifyRP® Acceler8® Discovery Kit

DOI

dx.doi.org/10.17504/protocols.io.8ephtdn

Théo Nass¹, Luc Gabel¹

¹EPFL - EPF Lausanne

iGEM EPFL



Dana Mozaffari

OPEN  ACCESS



DOI: dx.doi.org/10.17504/protocols.io.8ephtdn

Protocol Citation: Théo Nass, Luc Gabel 2019. RPA DNA Amplification using Agdia AmplifyRP® Acceler8® Discovery Kit. [protocols.io](https://dx.doi.org/10.17504/protocols.io.8ephtdn) <https://dx.doi.org/10.17504/protocols.io.8ephtdn>

License: This is an open access protocol distributed under the terms of the [Creative Commons Attribution License](#), which permits unrestricted use, distribution, and reproduction in any medium, provided the original author and source are credited

Protocol status: Working

We use this protocol and it's working

Created: October 18, 2019

Last Modified: October 21, 2019

Protocol Integer ID: 28847

Abstract

This protocol describes the procedure to perform an RPA DNA Amplification using the Agdia AmplifyRP® Acceler8® Discovery Kit

Materials

Rehydration Buffer	5.9 μl/sample
Primer A	0.42 μl/sample
Primer B	0.42 μl/sample
Template (DNA)	1.0 μl/sample
MgOAc (280 mM)	0.50 μl/sample
dH ₂ O	1.76 μl/sample

Quantities for preparation of a 10 μl RPA reaction using Agdia AmplifyRP® Acceler8® Discovery Kit



- 1 Prepare the RPA pellet rehydration solution following the table in the Materials section. If possible do a Master Mix that will then be divided.
- 2 For each sample, transfer the entire volume (10 μ l) to a reaction pellet. Pipette up and down until the full pellet has been resuspended.
- 3 Cap the reaction tubes. Vortex and spin briefly two times.
- 4 Transfer the reaction tubes to a heat block. Incubate at 39°C for 20 min.
- 5 After the incubation, purify the sample if you want to run the RPA product on an agarose gel.

PCR purification kits work well, but since RPA primers can be very long (60+ bp), a high cutoff is recommended.