

Jul 10, 2023

RT-PCR amplification for the discrimination between wild type and mutation alleles in SARS-CoV-2 genomes from wastewater

DOI

dx.doi.org/10.17504/protocols.io.dm6gp3ww1vzp/v1

Ana C Reis¹, Daniela Pinto¹, Mónica V. Cunha¹

¹Centre for Ecology, Evolution and Environmental Changes (cE3c) & CHANGE - Global Change and Sustainability Institute, Faculdade de Ciências, Universidade de Lisboa, Lisboa, Portugal



Daniela Pinto

Centre for Ecology, Evolution and Environmental Changes (cE3c...)

Create & collaborate more with a free account

Edit and publish protocols, collaborate in communities, share insights through comments, and track progress with run records.

Create free account

OPEN  ACCESS



DOI: <https://dx.doi.org/10.17504/protocols.io.dm6gp3ww1vzp/v1>

Protocol Citation: Ana C Reis, Daniela Pinto, Mónica V. Cunha 2023. RT-PCR amplification for the discrimination between wild type and mutation alleles in SARS-CoV-2 genomes from wastewater. **protocols.io**

<https://dx.doi.org/10.17504/protocols.io.dm6gp3ww1vzp/v1>

License: This is an open access protocol distributed under the terms of the **Creative Commons Attribution License**, which permits unrestricted use, distribution, and reproduction in any medium, provided the original author and source are credited



Protocol status: Working

We use this protocol and it's working

Created: July 10, 2023

Last Modified: July 10, 2023

Protocol Integer ID: 84761

Keywords: mutation alleles in sar, reverse transcription of the sar, applied biosystems taqman snp genotyping assay, target sequence region in each mutation assay, isolated rna from sar, mutation assay, reverse transcription, pcr for the allelic discrimination, mutation sequence, pcr reaction, pcr amplification for the discrimination, wastewater this protocol, mutation allele, positive wastewater sample, mutation alleles in specific position, rna, wastewater sample, pcr amplification, genome, isolated rna, pcr, wastewater, known sar, reverse primer, sar, target sequence region

Funders Acknowledgements:

"Systematic Surveillance of SARS-CoV-2 and its variants in wastewaters, Portugal (Ref. 060701/2021/864489/SUB/ENV.C2)", financed by The European Commission through the Emergency Support Instrument
Grant ID: 060701/2021/864489/SUB/ENV.C2

Abstract

This protocol describes the procedure of RT-PCR for the allelic discrimination between wild type and mutation alleles in specific positions of the SARS-CoV-2 genome. Starting with isolated RNA from SARS-CoV-2 positive wastewater samples, this workflow combines Applied Biosystems TaqMan SNP Genotyping Assays with a one-step RT-PCR reaction to detect whether there are known SARS-CoV-2 mutations present in the wastewater samples.

Sequence-specific forward and reverse primers amplify the target sequence region in each mutation assay. The reverse primer also primes reverse transcription of the SARS-CoV-2 genomic RNA sequences. Each mutation assay includes two TaqMan minor groove binder probes with nonfluorescent quenchers: one VIC dye labelled probe to detect the reference sequence and one FAM dye labelled probe to detect the mutation sequence.

Materials

- TaqPath 1-Step RT-qPCR Master Mix CG, ThermoFisher Scientific, Catalog A15299
- AcroMterix SARS-CoV-2 Control, ThermoFisher Scientific, Catalog 954519
- Mutation Assay: S.G339D.GGT.GAT, ThermoFisher Scientific, Catalog CV47VRX
- Mutation Assay: S.Q493R.CAA.CGA, ThermoFisher Scientific, Catalog CVH49P2
- Mutation Assay: S.L452R.CTG.CGG, ThermoFisher Scientific, Catalog CVAAAAD
- Mutation Assay: S.T547K.ACA.AAA, ThermoFisher Scientific, Catalog CVYMJGA
- Mutation Assay: M.D3N.GAT.AAT, ThermoFisher Scientific, Catalog CVAAAAK
- Mutation Assay: ORF7b.L11F.TTG.TTT, ThermoFisher Scientific, Catalog CVCE3VH

Troubleshooting

Before start

Wipe bench surfaces with RNase Away and keep working environment clean. Maintain the RNA samples on ice, and prepare the RT-PCR mixture on ice.



RT-PCR amplification

- 1 Prepare the amplification mixture as follows:

Reagent	Concentration in the reaction
TaqPath 1-Step RT-qPCR Master Mix CG (4X)	1X
Mutation assay (40X)	2X
DNase/RNase free water	up to 15 µL

To the 15 µl of reaction mix, add 5 µl of wastewater RNA sample, or no-template control (DNase/RNase free water), or a wild-type AcroMetrix Coronavirus 2019 (COVID-19) RNA control, to complete 20 µl of reaction volume.

- 2 Thermocycling conditions

Step	Temperature (°C)	Time	No. Cycles
Reverse transcription	45	15'	
Taq polymerase activation	95	2'	
Denaturation	95	15''	45
Annealing	58	45''	
Post-read	60	30'	

Set the detection module for FAM and VIC labelled probes.

Data analysis

- 3 The allelic discrimination results can be graphed on a scatter plot contrasting reporter dye florescence (*i.e.*, allele X versus allele Y), and analyzed with the QuantStudio Design and Analysis Software, using the genotyping analysis module with real-time data.