

Aug 26, 2022

RT-QuIC alpha-synuclein

DOI

<https://dx.doi.org/10.17504/protocols.io.q26g7yzb3gwz/v1>

maria.xylaki¹

¹University Medical Center Goettingen



Mary Xylaki

UMG

Create & collaborate more with a free account

Edit and publish protocols, collaborate in communities, share insights through comments, and track progress with run records.

Create free account

OPEN  ACCESS



DOI: <https://dx.doi.org/10.17504/protocols.io.q26g7yzb3gwz/v1>

External link: <https://actaneurocomms.biomedcentral.com/articles/10.1186/s40478-021-01282-8>,
https://link.springer.com/protocol/10.1007/978-1-4939-9124-2_4

Protocol Citation: maria.xylaki 2022. RT-QuIC alpha-synuclein . protocols.io
<https://dx.doi.org/10.17504/protocols.io.q26g7yzb3gwz/v1>

Manuscript citation:

Russo MJ, Orru CD, Concha-Marambio L, Giaisi S, Groveman BR, Farris CM, et al. High diagnostic performance of independent alpha-synuclein seed amplification assays for detection of early Parkinson's disease. *Acta Neuropathol Commun. BioMed Central*; 2021;9:1–13. Concha-Marambio L, Shahnawaz M, Soto C. Detection of Misfolded α -Synuclein Aggregates in Cerebrospinal Fluid by the Protein Misfolding Cyclic Amplification Platform. 2019. p. 35–44.

License: This is an open access protocol distributed under the terms of the **Creative Commons Attribution License**, which permits unrestricted use, distribution, and reproduction in any medium, provided the original author and source are credited

Protocol status: Working

We use this protocol and it's working

Created: August 25, 2022

Last Modified: May 31, 2024

Protocol Integer ID: 69179

Keywords: ASAPCRN, synuclein forms in human cerebrospinal fluid, synuclein form, synuclein, detection of prionoid alpha, human cerebrospinal fluid, prionoid alpha, quic alpha, quic, diagnostic

Disclaimer

DISCLAIMER – FOR INFORMATIONAL PURPOSES ONLY; USE AT YOUR OWN RISK

The protocol content here is for informational purposes only and does not constitute legal, medical, clinical, or safety advice, or otherwise; content added to **protocols.io** is not peer reviewed and may not have undergone a formal approval of any kind. Information presented in this protocol should not substitute for independent professional judgment, advice, diagnosis, or treatment. Any action you take or refrain from taking using or relying upon the information presented here is strictly at your own risk. You agree that neither the Company nor any of the authors, contributors, administrators, or anyone else associated with **protocols.io**, can be held responsible for your use of the information contained in or linked to this protocol or any of our Sites/Apps and Services.

Abstract

This protocol is for the detection of prionoid alpha-synuclein forms in human cerebrospinal fluid using the Real Time Quacking-Induced Conversion method (RT-QulC). The protocol is adapted from Marco J. Russo, Christina D. Orru, Luis Concha-Marambio, Simone Giaisi et al., 2021 (doi:10.1186/s40478-021-01282-8) and Concha-Marambio et al., 2019 (doi:10.1007/978-1-4939-9124-2_4).

This assay is for research use only and not diagnostics.

Materials

- BMG Technologies Omega FLUOstar Reader or similar fluorometer
- regular lab equipment (balance, pHmeter, pipetts, tips, tubes)
- COSTAR 96-well ELISA plates (Corning, cat# 3916)
- MicroAmp Film (Applied Biosystems, cat# 4311971)
- Si_3N_4 beads 2.38 mm
- molecular biology grade nuclease free water
- 0.5 M PIPES pH 6.5
- 5 M NaCl
- Thioflavin T dissolven in water
- recombinant C-terminal His-tag alpha synuclein
- 1% BSA
- 5 N NaOH and 1N HCl for pH adjustment



Bead preparation

- 1 Beads are blocked with 1% BSA in 100mM PIPES for 1 hr and washed twice with PIPES.

Preparation of Reaction buffer

- 2 Reaction mixture is prepared calculating 200 µl per well with the following final concentrations:
 - 0.3 mg/mL recombinant alpha-synuclein
 - 0.5 M NaCl
 - 100 mM PIPES buffer
 - 5 µM ThT

Seeting up the assay

- 3 One bead is placed in each well of the assay plate. 160 µl reaction buffer and 40 µl CSF are carefully pipetted in each well. Samples are assessed in triplicates.
- 4 Plate is covered with the film and creases are removed manually.
- 5 Plate is placed in the plate reader and incubated for 240 hrs at 37 °C in cycles of 1 min shaking at 500rpm, 29 mins incubation and fluorescence measurements are taken after every incubation cycle at 440ex/490emm.

Data analysis

- 6 A sample is considered positive when it crosses a fluorescence threshold established at 3 standard deviations above baseline. 3/3 wells are positive and negative when 1/3 samples are positive while 2/3 is considered as incocclusive.
- 7 Relative fluorescence units measured at 490 are plotted versus the time in hours and should present a classical exponantial curve with lag phase and plateau. Kinetics parameters obtained from the curves include maximum fluorescence, time to reach 50% of maximal fluorensence and time to reach the established threshold.