

Sep 04, 2023

Salivary collection participant instructions using OMNIgene ORAL OM 501 device

DOI

<https://dx.doi.org/10.17504/protocols.io.n92ldm8qol5b/v1>

Elisa Menozzi^{1,2}, Roxana Mezabrovski^{1,2}, Sara Lucas^{1,2}, Jane Macnaughtan^{1,2}, Sofia Koletsi^{1,2}, Aleesa Nazeer^{1,2}, Nadine Loefflad^{1,2}

¹Department of Clinical and Movement Neurosciences, Queen Square Institute of Neurology, University College London (UCL), London, United Kingdom;

²Aligning Science Across Parkinson's (ASAP) Collaborative Research Network, Chevy Chase, MD, 20815



Elisa Menozzi

Create & collaborate more with a free account

Edit and publish protocols, collaborate in communities, share insights through comments, and track progress with run records.

Create free account

OPEN  ACCESS



DOI: <https://dx.doi.org/10.17504/protocols.io.n92ldm8qol5b/v1>

Protocol Citation: Elisa Menozzi, Roxana Mezabrovski, Sara Lucas, Jane Macnaughtan, Sofia Koletsi, Aleesa Nazeer, Nadine Loefflad 2023. Salivary collection participant instructions using OMNIgene ORAL OM 501 device. **protocols.io**
<https://dx.doi.org/10.17504/protocols.io.n92ldm8qol5b/v1>

License: This is an open access protocol distributed under the terms of the **Creative Commons Attribution License**, which permits unrestricted use, distribution, and reproduction in any medium, provided the original author and source are credited

Protocol status: Working

We use this protocol and it's working

Created: September 04, 2023

Last Modified: May 31, 2024

Protocol Integer ID: 87311

Keywords: ASAPCRN, salivary collection participant instruction, omnigene oral om 501 device, salivary sample, salivary samples from participant, dna genotek, oral om, om 501 device, dna, using omnigene, sample, protocol, instruction, adapted protocol, participant, device, protocol for collection

Funders Acknowledgements:

Aligning Science Across Parkinson's

Grant ID: ASAP-000420

Abstract

This is an adapted protocol for collection of salivary samples from participants. This protocol was adapted from DNA Genotek's instructions available at: <https://www.dnagenotek.com/us/support/collection-instructions/omnigene-oral/OM-501.html>



1



ASAL_saliva_collection_protocol.pdf 309.2KB