

Oct 16, 2023

Sample vitrification and cryo-EM data acquisition

DOI

<https://dx.doi.org/10.17504/protocols.io.kqdg39rreg25/v1>

Minghao Chen¹

¹Department of Molecular and Cell Biology, University of California, Berkeley, Berkeley, CA 94720, USA.



Minghao Chen

University of California, Berkeley

Create & collaborate more with a free account

Edit and publish protocols, collaborate in communities, share insights through comments, and track progress with run records.

Create free account

OPEN  ACCESS



DOI: <https://dx.doi.org/10.17504/protocols.io.kqdg39rreg25/v1>

Protocol Citation: Minghao Chen 2023. Sample vitrification and cryo-EM data acquisition. **protocols.io**

<https://dx.doi.org/10.17504/protocols.io.kqdg39rreg25/v1>

License: This is an open access protocol distributed under the terms of the **Creative Commons Attribution License**, which permits unrestricted use, distribution, and reproduction in any medium, provided the original author and source are credited

Protocol status: Working

We use this protocol and it's working

Created: June 05, 2023

Last Modified: May 31, 2024

Protocol Integer ID: 82920

Keywords: ASAPCRN, em data acquisition sample vitrification, sample vitrification, em data acquisition, data, cryo, sample

Funders Acknowledgements:

Aligning Science Across Parkinson's (ASAP) initiative

Grant ID: ASAP-000350

Abstract

Sample vitrification and cryo-EM data acquisition

Sample vitrification

- 1 Glow-discharge the grids at 25 mA for 30 sec with PELCO easiGlow system (Ted Pella)
- 2 Apply 3 μ l of protein solution to the grid:
QUANTIFOIL R1.2/1.3 mesh 300 (Electron Microscopy Sciences), or
QUANTIFOIL R2/1 mesh 300 (Electron Microscopy Sciences)
- 3 Vitrify the sample with a Vitrobot cryo-plunger (Thermo Fisher Scientific)
- 4 (Optional) Add 0.05%(w/v) of n-Octyl-Beta-D- Glucopyranoside as a surfactant before vitrification.

cryo-EM data acquisition at 300 kV Titan Krios microscope

- 5 Collect dataset at a magnification of 81,000x and a corresponding pixel size of 1.05 Å and a defocus range of -0.8 to -2.0 μ m. Image stacks contain 50 frames with a total dose of 50 e/Å².

cryo-EM data acquisition at 200 kV Talos Arctica microscope

- 6 Collect dataset at a magnification of 36,000x and a corresponding pixel size of 0.5575 Å and a defocus range of -0.8 to -2.0 μ m in a super-resolution correlated-double sampling mode. Image stacks contain 50 frames with a total dose of 50 e/Å².