

Oct 18, 2019

Sanger sequencing

DOI

dx.doi.org/10.17504/protocols.io.8fghtjw

Igem Dusseldorf¹

¹Heinrich-Heine Universität Düsseldorf



Igem Dusseldorf

Create & collaborate more with a free account

Edit and publish protocols, collaborate in communities, share insights through comments, and track progress with run records.

Create free account

OPEN  ACCESS



DOI: <https://dx.doi.org/10.17504/protocols.io.8fghtjw>

Protocol Citation: Igem Dusseldorf 2019. Sanger sequencing. protocols.io <https://dx.doi.org/10.17504/protocols.io.8fghtjw>

License: This is an open access protocol distributed under the terms of the [Creative Commons Attribution License](#), which permits unrestricted use, distribution, and reproduction in any medium, provided the original author and source are credited

Protocol status: Working

We use this protocol and it's working

Created: October 18, 2019

Last Modified: October 18, 2019

Protocol Integer ID: 28872

Troubleshooting



- 1 Label 1.5 mL tubes with barcodes from sequencing company
- 2 . Add template (400 to 500 ng), primer and water (10 uL total volume, details below)

		uL per Reac tion
	Tem plate	400 -500 ng
	10 uM Prim er	2.5
	Wat er	fill up with wate r
	Total Volu me	10.0

- 3 Place tubes inside a small plastic bag and drop off in sequencing company collection Box
- 4 Results are usually ready to download from the sequencing companys website