

Sep 23, 2020

Version 1

SapphireAmp PCR Master Mix -- CHEM 584 V.1

DOI

<https://dx.doi.org/10.17504/protocols.io.bmc7k2zn>

Ken Christensen¹

¹Brigham Young University

LAB PROTOCOLS



Ken Christensen

Brigham Young University

Create & collaborate more with a free account

Edit and publish protocols, collaborate in communities, share insights through comments, and track progress with run records.

Create free account

OPEN  ACCESS



DOI: <https://dx.doi.org/10.17504/protocols.io.bmc7k2zn>

Protocol Citation: Ken Christensen 2020. SapphireAmp PCR Master Mix -- CHEM 584. **protocols.io**
<https://dx.doi.org/10.17504/protocols.io.bmc7k2zn>

Manuscript citation:

https://www.takarabio.com/assets/documents/User%20Manual/RR350A_DS.v1902Da.pdf

License: This is an open access protocol distributed under the terms of the [Creative Commons Attribution License](#), which permits unrestricted use, distribution, and reproduction in any medium, provided the original author and source are credited

Protocol status: Working

We use this protocol and it's working

Created: September 14, 2020

Last Modified: September 23, 2020

Protocol Integer ID: 42111

Keywords: ideal for fast colony pcr screening, sapphireamp fast pcr master mix, fast colony pcr screening, sapphireamp pcr master mix, fast colony pcr amplification, fast pcr master mix, fast pcr, hot start pcr enzyme, pcr, direct loading of pcr product, agarose gel for electrophoresi, pcr product, sapphireamp, gel loading dye, electrophoresi, genomic dna template, kb from genomic dna template, optimized buffer, sample handling

Abstract

SapphireAmp Fast PCR Master Mix contains a hot start PCR enzyme, optimized buffer, dNTP mixture, gel loading dye (blue), and a density reagent as a 2X premix. SapphireAmp Fast PCR Master Mix is optimized for fast PCR and offers a rapid extension rate (10 sec. per kb). The inclusion of blue dye and a density reagent allows direct loading of PCR products on an agarose gel for electrophoresis. The master mix format simplifies workflows and sample handling; simply add primers, template, and water and then begin PCR. SapphireAmp Fast PCR Master Mix is ideal for fast colony PCR screening. Fast colony PCR amplification of a 5 kb insert can be completed in approximately 1 hr 15 min. Furthermore, it is possible to amplify fragments up to 6 kb from genomic DNA templates.

Guidelines

Storage: -20°C for long-term storage. 4°C for short-term storage (up to 3 months). (Note) If used frequently, store at 4°C ; the activity of the Master Mix may decrease with repeated freezing and thawing. Gently mix well before use and centrifuge briefly.

Application:

- DNA amplification by PCR
- Colony PCR

PCR Products: Since most PCR products amplified with SapphireAmp Fast PCR Master Mix have an A overhang added at 3'-termini, the obtained PCR product can be used directly for cloning into a T-vector. Additionally, it is possible to clone the product in a blunt-end vector after blunting and phosphorylation of the end.

Dye Migration During Electrophoresis: When 5 µl of the PCR sample is loaded on a 1% gel made with Agarose L03 [TAKARA] (Cat. #5003) and subjected to electrophoresis, the blue dye fronts are detected at positions corresponding to 1 kb and 3 - 5 kb. The absorption maxima for the dyes are ~ 260 nm and 620 nm, respectively. The dyes may be removed by isolating and purifying the DNA fragment from the gel or extracting DNA with NucleoSpin Gel and PCR Clean-Up (Cat. #740609.50/.250), if necessary

Setup Reaction

- 1 To a 12.5 μL aliquot of SapphireAmp PCR Master Mix, add template (4 μL cleared lysate for colony PCR or < 1 ng of purified DNA for typical PCR), forward and reverse primers to a final concentration of 100 nanomolar (nM) . Adjust final volume to 25 μL with nuclease free water or autoclaved water.

Sapphire Master Mix	12.5 ul (pre- aliquoted and stored in the freezer)
Template	4 ul of bacterial lysate or <1 ng DNA
Forward Primer	0.5 ul of 10 uM primer dilution
Reverse Primer	0.5 ul of 10 uM primer dilution
ddH ₂ O	to a final volume

of
25 ul

Run Reaction

2 followed by 30 cycles of 98°C, 5 sec; 55°C, 5 sec; and 72°C, 40 sec.

	Initial denature	98C	1 minu te
	Denature	94C	10 seco nds
	Anneal	55C	10 seco nds
	Extension	72C	30 seco nds
	Repeat steps 2-4		30- 40x
	Final extension	72	1 minu te
	Cool	4C	Until canc elled

A typical thermocycling program for a quick PCR for amplicons less than 1 kb. For longer amplicons, adjust the program to 10 seconds per kb for the extension and final extension times.