


Dec 16, 2022

Screening procedure to identify triazole-resistant *Aspergillus* spp. using agar plates

DOI

<https://dx.doi.org/10.17504/protocols.io.6qpvr492zgmk/v1>

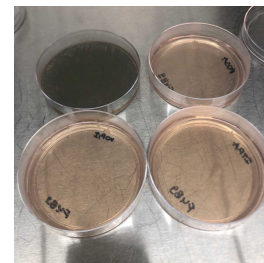

Juliê Alves¹

¹Universidade Federal de Ciências da Saúde de Porto Alegre



Juliê JA Alves

Universidade Federal de Ciências da Saúde de Porto Alegre

Create & collaborate more with a free account

Edit and publish protocols, collaborate in communities, share insights through comments, and track progress with run records.

Create free account

OPEN  ACCESS



DOI: <https://dx.doi.org/10.17504/protocols.io.6qpvr492zgmk/v1>

Protocol Citation: Juliê Alves 2022. Screening procedure to identify triazole-resistant *Aspergillus* spp. using agar plates. protocols.io <https://dx.doi.org/10.17504/protocols.io.6qpvr492zgmk/v1>

License: This is an open access protocol distributed under the terms of the **Creative Commons Attribution License**, which permits unrestricted use, distribution, and reproduction in any medium, provided the original author and source are credited

Protocol status: Working

We use this protocol and it's working

Created: December 15, 2022

Last Modified: December 16, 2022

Protocol Integer ID: 74082

Keywords: aspergillus spp, genus aspergillus spp, visual analysis of fungal growth, fungal growth, resistance to azole, posaconazole, itraconazole, voriconazole, triazole, detecting resistance, agar plates the methodology, using agar plate, screening procedure, azole, agar plate, resistant aspergillus spp

Abstract

The methodology of the screening procedure for detecting resistance to azoles in the genus *Aspergillus* spp. consists of an in house technique, with visual analysis of fungal growth on agar plates supplemented with azoles, Itraconazole at a concentration of 4 µg/ml, Voriconazole 2 µg/ml and Posaconazole 0.5 µg/ml. This protocol is based on the E.Def 10.1 from EUCAST

Materials

Reagents:

1L of sterile water:  1000 µL

RPMI 1640 powder –  10.4 g

D-Glucose 2% -  18 g

Bacteriological agar –  20 g

MOPS -  34.53 g (0.165 M)

Materials:

- 4 autoclavable vials (for each antifungal and control)
- 15ml Falcon type tube
- Pasteur pipette
- Pipettes and tips: 1000, 200, 100, 20, 10
- 1L beaker
- Funnel
- Glass stick
- 90 petri plates 60mmx15mm (If you use petri plates divided into 4 the medium will yield more)
- Swabs
- Eppendorfs

Before start

Separate all reagents and identify the plates



Preparation of antifungal solutions:

- 1 From the antifungal powder prepare stock solutions in DMSO, at concentrations at least 200 times higher than those tested on the agar plate
- 2 From the stock solution, prepare the working solutions at the following concentrations:
Itraconazole (ITZ) - 400 µg/ml,
Voriconazole (VRZ) - 200 µg/ml
Posaconazole (PSZ) - 50 µg/ml

Storage Temperature: -25 to -18° C (Freezer)

For calculations, use the formula:

Initial concentration x Initial volume = Final concentration x Final volume

	A	B	C	D	E	F
		Stock Solution Concentration	Stock Solution Volume	Volume of DMSO	Working Solution Concentration	Agar Final Concentration
		µg/mL	µL	µL	µg/mL	µg/mL
	ITZ	1600	2500	7500	400	4
	VRZ	400	5000	5000	200	2
	PSZ	400	1250	8750	50	0,5

Example: Preparation of 10 mL of working solution of each of the azoles from the stock solution

Preparation of the medium:

5m

- 3 Weigh all reagents:
RPMI 1640 powder –  10.4 g
D-Glucose 2% -  18 g
Bacteriological agar –  20 g
MOPS -  34.53 g (0.165 M)




- 4 In a 1L becker, Add  900 mL of distilled water and RPMI powder  10.4 g , glucose 2%  18 g , Bacto Agar  20 g , MOPS  34.53 g
- 5 Microwave  00:05:00 , little by little, stirring with a glass rod, until the solutes are completely dissolved
- 6 Confirm the pH of 7.0

5m



6.1 If necessary adjust with a solution of NaOH (10 M) to 7.0

7 Add more  100 mL of distilled water

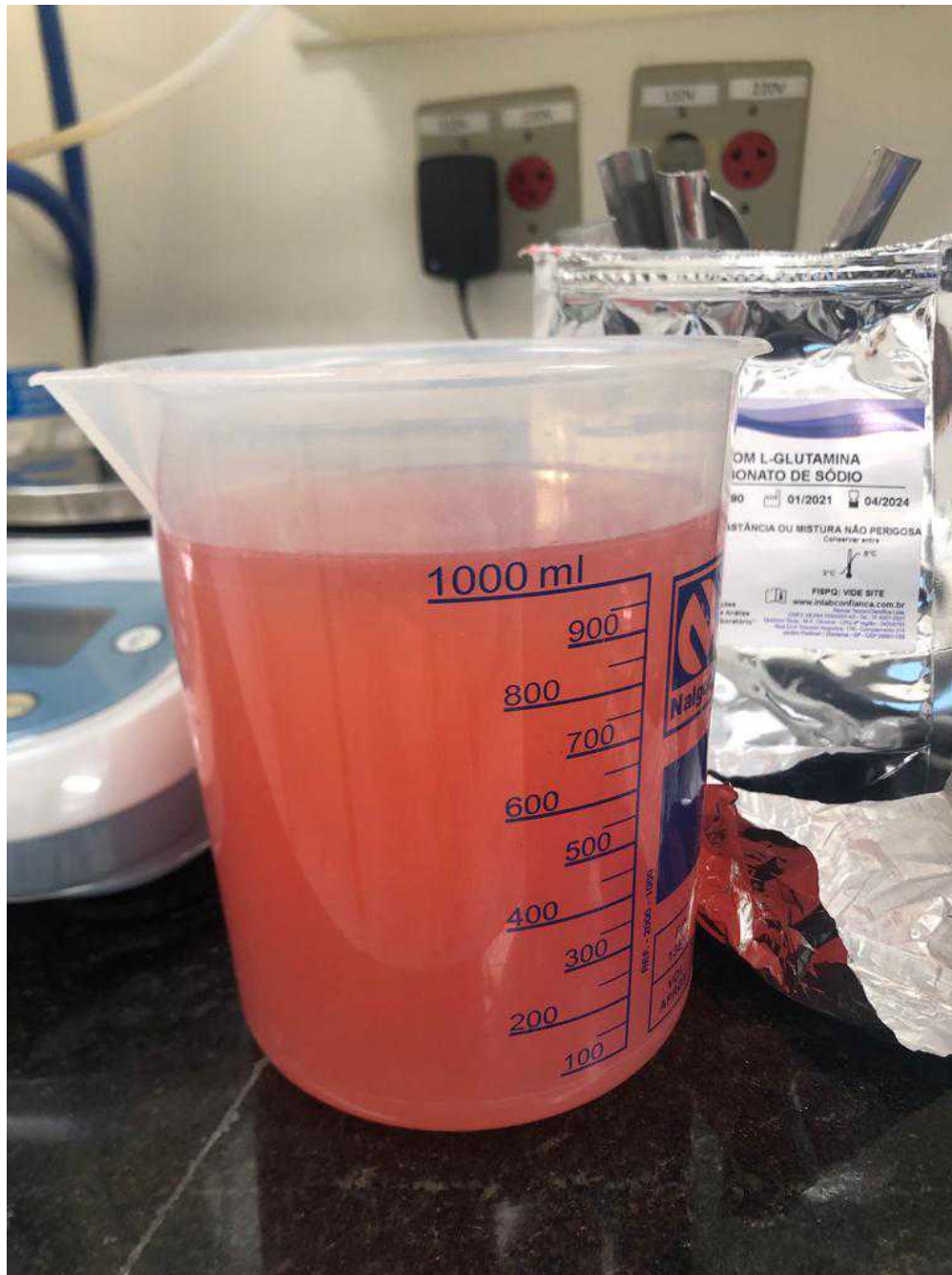

Incorrect beaker measurements, it was measured with the test tube.

8

Divide the medium into 4 Erlenmeyer flasks containing  250 mL each

9

Autoclave the medium at 121° for  00:15:00 .


15m



Preparation of the azole-containing plates:

- 10 Identify all the plates with the azole or "control"
- 11 Cool the agar down to 45°C
- 12 Add  2.5 mL of each azole working solution (100% final concentration) to each of the vials and ensure proper mixing
- 13 The remaining antifungal free vial will be used for the growth control
- 14 Pour the agar from each vial into the respective plates
- 15 After solidification of the plates (approximately  00:20:00) Storage the plates at 4°C for up to 6 months from the date of preparation.

20m

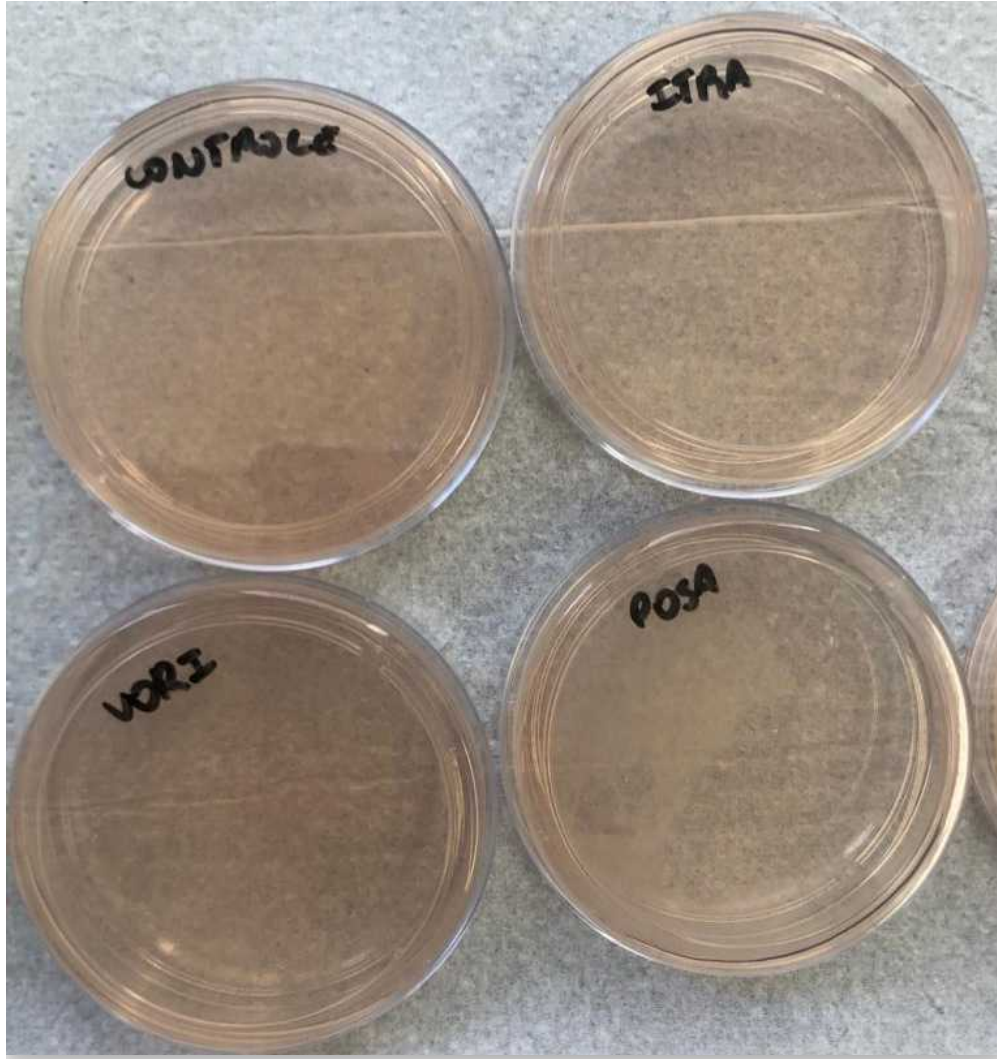

Preparation of the inoculum:

- 16 For inoculum suspensions use fresh and mature cultures incubated for at least 2–7 days
- 16.1 The isolates should be cultured on Sabouraud dextrose agar (with or without chloramphenicol) or other culture medium where the fungus is able to sporulate sufficiently and incubated at 34–37°C




Aspergillus fumigatus culture ready for the inoculum suspension

- 17 Make a diluent stock solution by adding 1 ml of Tween 20 to every 50 ml of distilled water
- 18 Working in the biological safety cabinet, add 3 ml of the water/Tween 20 solution to the 3-7 day slope. Tween 20 will help to put the *Aspergillus* conidia into suspension.
- 19 If necessary gently detach only the conidia from the culture with a handle until the solution becomes turbid
- 20 Put  1.5 mL of this conidial suspension in identified eppendorfs



- 20.1 This conidial suspension can be used in case of retest
- 21 Use the residual solution from the tube from steps 17-20 to perform the next steps
- 22 Add  2 mL of distilled water in a tube and transfer 1-2 drops of the residual solution with conidial
- 22.1 Or transfer 1-2 drops of the conidial suspension

- 22.2 In some cases you will need more drops, for *Aspergillus flavus* for example
- 23 Visually compare the suspension to the 0.5 McFarland Standard and make adjustments as described until the suspension visually matches the standard

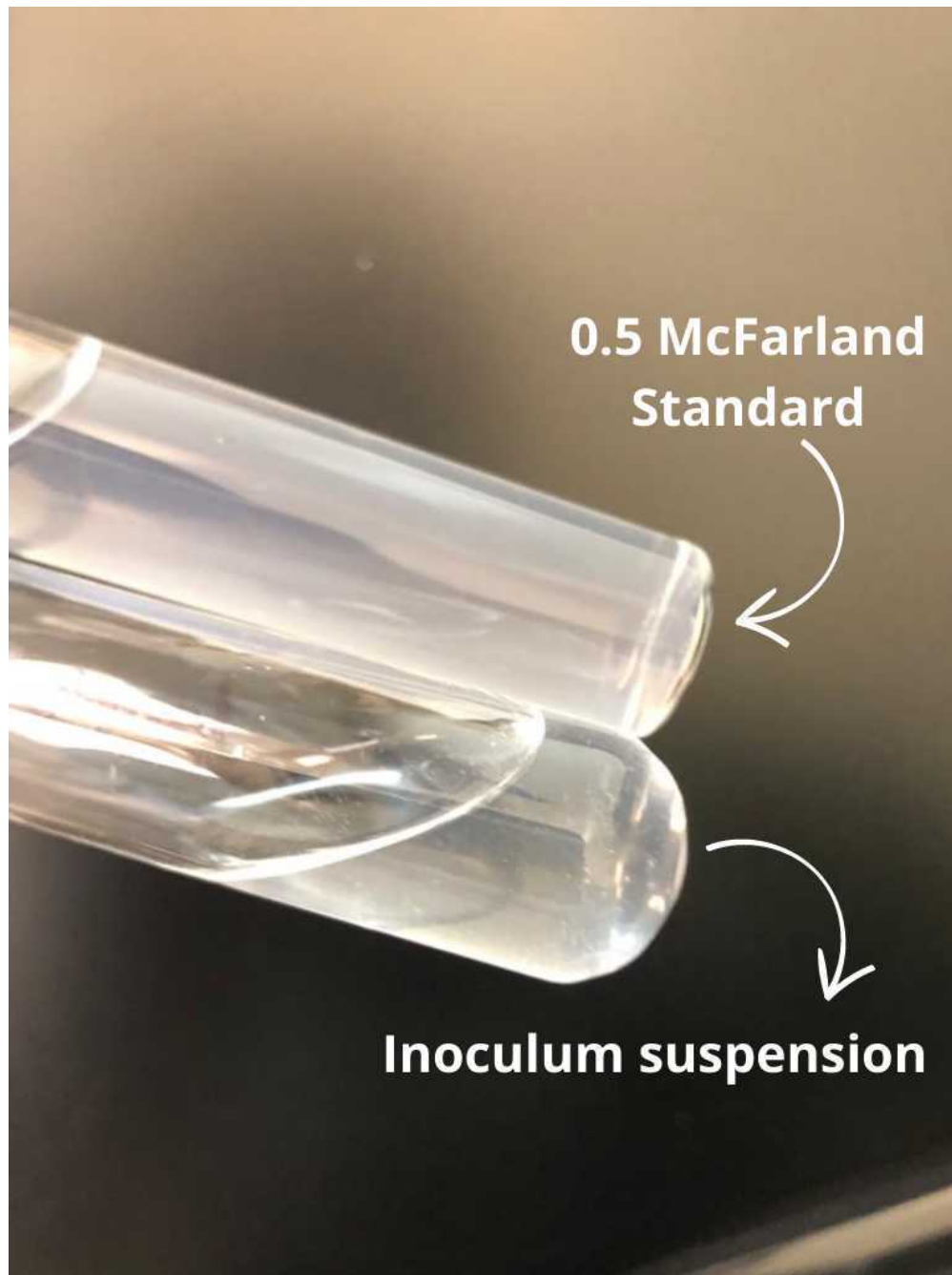



- 23.1 Adjustment is made by adding more distilled water if the suspension is very turbid and more conidia if the turbidity is very low.

Inoculation and incubation of agar plates:

2d

- 24 Dip a sterile cotton swab in the inoculum suspension and rub the entire surface of the control plate in a back-and-forth motion
- 25 Re-dip the cotton swab and rub the entire surface in the same way of each of the three remaining plates with the azoles
- 26 Incubate the plates at 35-37°C for  48:00:00 and observe growth.
- 26.1 Visual analysis score definition:
- 0: no visible growth
 - 1: weak/minimal growth (such as >5 tiny colonies or confluent weak growth where isolate was inoculated (covering $\leq$ half of the plate)
 - 2: clearly visible growth with hyphal extension, but not covering entire plate
 - 3: prominent uninhibited growth covering most of the plate


2d

Analyzing the results


- 27 The presence or absence of fungal growth on the surface of the plates will yield the following preliminary classification:
- **Azole susceptible isolate:** growth on the azole-free plate and absence of growth on the others plates
 - **Potentially an azole non-susceptible isolate:** growth on the azole-free plate and on any of the other supplemented plates.

Expected result

The control plate should always show growth (score definition = 3)
The azole containing plates will show growth only if the isolate is potentially resistant to the azole.

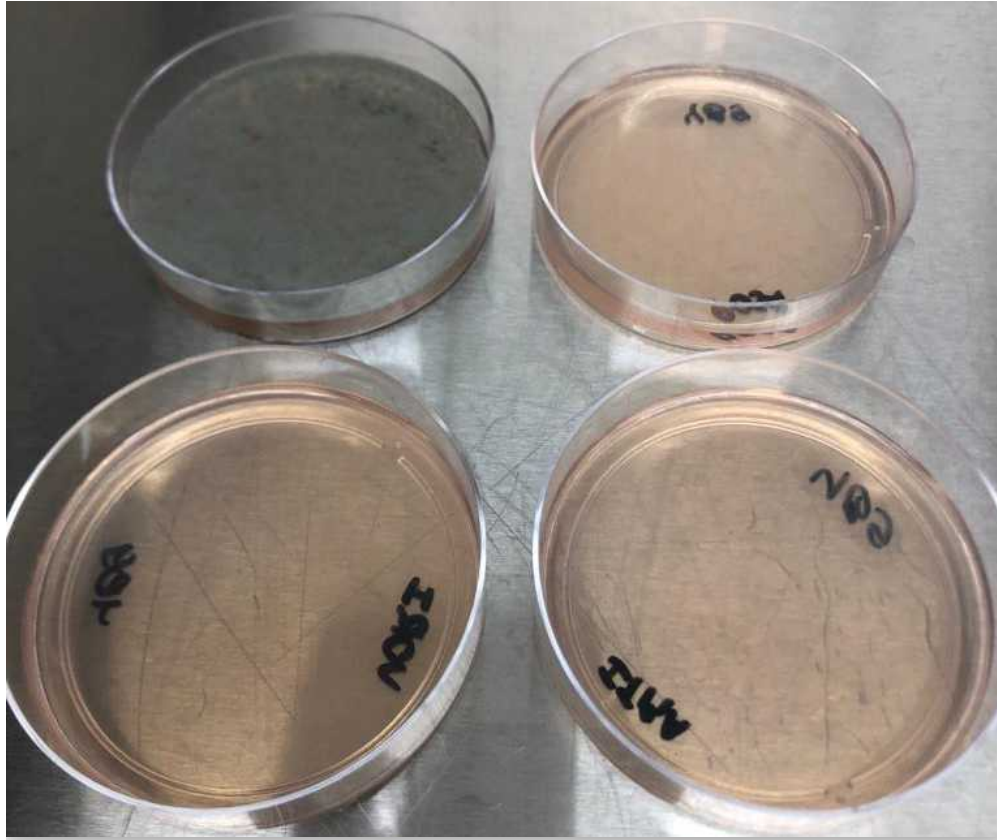

Azole susceptible isolate - Being the upper left control plate, without azole, upper right PSZ 0.5 $\mu\text{g/mL}$, lower left VRZ 2 $\mu\text{g/mL}$ and lower right ITZ 4 $\mu\text{g/mL}$.

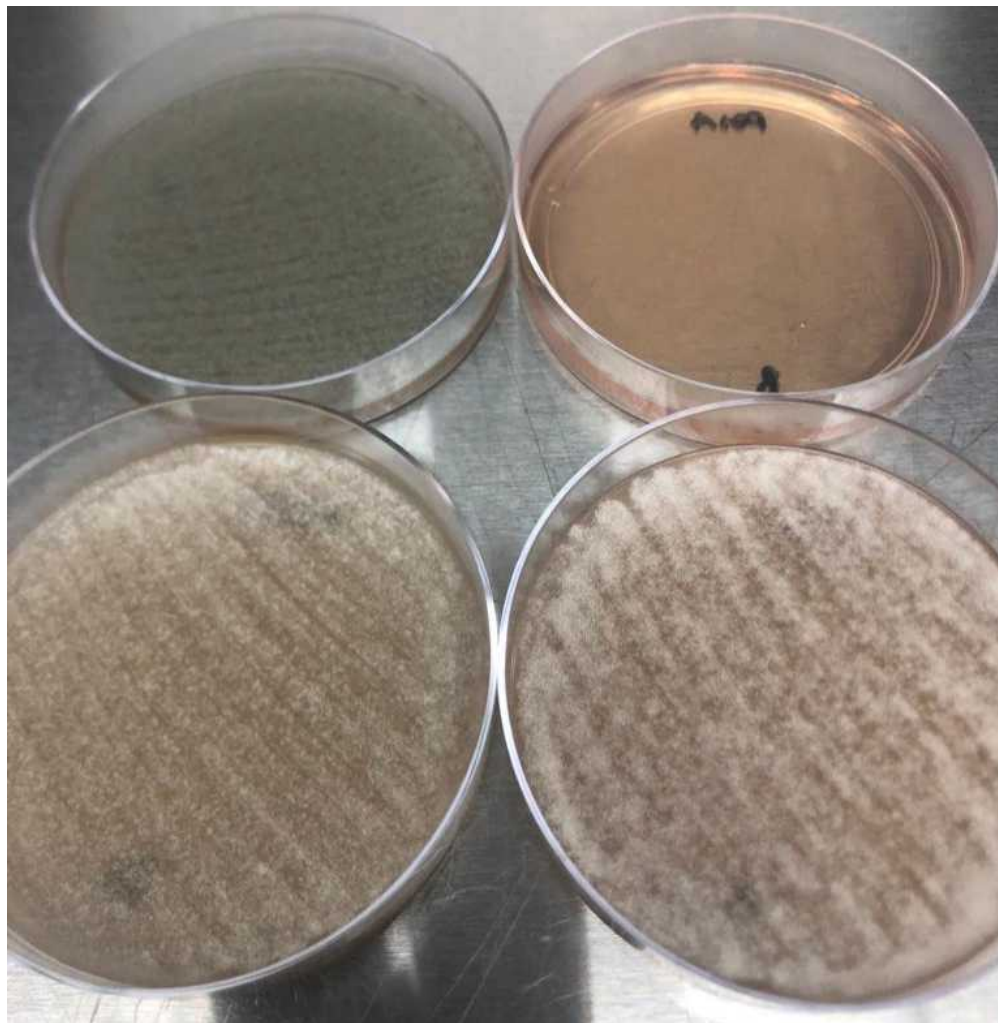

Potentially an ITZ and VRZ resistant isolate - Being the upper left control plate, without azole, upper right PSZ 0.5 µg/mL, lower left VRZ 2 µg/mL and lower right ITZ 4 µg/mL.