



Oct 14, 2024

SDS-PAGE and immunoblotting to assess whole-cell lysates and Endo-IPs and Lyso-IPs in hESCs and iNeurons

 In 1 collection

DOI

<https://dx.doi.org/10.17504/protocols.io.q26g71eq9gwz/v1>

Frances V Hundley^{1,2}, Miguel A. Gonzalez-Lozano^{1,2}, Harper JW^{1,2}

¹Harvard Medical School;

²Aligning Science Across Parkinson's (ASAP) Collaborative Research Network, Chevy Chase, MD 20815, USA



Frances V Hundley

Harvard Medical School

Create & collaborate more with a free account

Edit and publish protocols, collaborate in communities, share insights through comments, and track progress with run records.

Create free account

OPEN  ACCESS



DOI: <https://dx.doi.org/10.17504/protocols.io.q26g71eq9gwz/v1>

External link: <https://doi.org/10.1038/s41586-025-09059-y>

Protocol Citation: Frances V Hundley, Miguel A. Gonzalez-Lozano, Harper JW 2024. SDS-PAGE and immunoblotting to assess whole-cell lysates and Endo-IPs and Lyso-IPs in hESCs and iNeurons. **protocols.io**

<https://dx.doi.org/10.17504/protocols.io.q26g71eq9gwz/v1>



Manuscript citation:

Gonzalez-Lozano, M.A., Schmid, E.W., Miguel Whelan, E. *et al.* EndoMAP.v1 charts the structural landscape of human early endosome complexes. *Nature* **643**, 252–261 (2025). <https://doi.org/10.1038/s41586-025-09059-y>

License: This is an open access protocol distributed under the terms of the **Creative Commons Attribution License**, which permits unrestricted use, distribution, and reproduction in any medium, provided the original author and source are credited

Protocol status: Working

We use this protocol and it's working

Created: October 02, 2024

Last Modified: October 14, 2024

Protocol Integer ID: 108855

Keywords: ASAPCRN, negative control markers from other organelle, endolysosomal protein, cell lysate, other organelle, negative control marker, neuron, ips in hesc, protein, ips from hesc, general protocol for western blotting

Funders Acknowledgements:

Aligning Science Across Parkinson's

Grant ID: ASAP-025160

Disclaimer

DISCLAIMER – FOR INFORMATIONAL PURPOSES ONLY; USE AT YOUR OWN RISK

The protocol content here is for informational purposes only and does not constitute legal, medical, clinical, or safety advice, or otherwise; content added to **protocols.io** is not peer reviewed and may not have undergone a formal approval of any kind. Information presented in this protocol should not substitute for independent professional judgment, advice, diagnosis, or treatment. Any action you take or refrain from taking using or relying upon the information presented here is strictly at your own risk. You agree that neither the Company nor any of the authors, contributors, administrators, or anyone else associated with **protocols.io**, can be held responsible for your use of the information contained in or linked to this protocol or any of our Sites/Apps and Services.

Abstract

We present a general protocol for Western blotting to detect endolysosomal proteins and negative control markers from other organelles to assess whole-cell lysates and Endo-IPs and Lyso-IPs from hESCs and iNeurons.



SDS-PAGE and immunoblotting

5m

- 1 For whole-cell lysates, prepare samples following standard protocols, and final samples should be in LDS buffer with DTT or similar. For Endo-IP and/or Lyso-IP from hESCs and/or iNeurons, post-nuclear supernatant (PNS), flow through, and eluate samples should be in LDS buffer with DTT or similar. Incubate samples at 80 °C for 00:05:00 .
- 2 Load samples into a 4-20% Criterion TGX Stain-free precast gel (Bio-Rad) and separate by electrophoresis in 1xTris/Glycine/SDS buffer.
- 3 Scan gel for total protein using a ChemiDoc MP imager (Bio-Rad).
- 4 Transfer proteins to PVDF or nitrocellulose membranes by standard wet transfer in 20% methanol Tis/Glycine buffer.
- 5 Block membrane in blocking buffer (5% non-fat dry milk or 3% BSA in TBST) at Room temperature for 01:00:00 .
- 6 Incubate membrane in primary antibody solution (blocking solution plus primary antibody at 1:500-1:1,000, depending on the primary antibody) at 4 °C for 12:00:00 to 16:00:00 .
- 7 Wash membrane six times with TBST for 00:05:00 each wash.
- 8 Incubate membrane in secondary antibody solution (blocking solution plus secondary antibody conjugated to HRP at 1:5,000-1:10,000) at Room temperature for 01:00:00 .
- 9 Wash membrane four times with TBST for 00:05:00 each wash.
- 10 Apply Western Lightning Plus Chemiluminescence substrate (Revvity) to membrane and acquire blot images using a ChemiDoc MP imager.

5m

1h

1d 4h

5m

1h

5m



- 11 Process raw image files with Image Lab software (Bio-Rad).

Protocol references

Frances V. Hundley, Miguel A. Gonzalez-Lozano, Lena M. Gottschalk, Aslan N. K. Cook, Jiuchun Zhang, Joao A. Paulo, J. Wade Harper. Endo-IP and Lyso-IP Toolkit for Endolysosomal Profiling of Human Induced Neurons
bioRxiv 2024.09.24.614704; doi: <https://doi.org/10.1101/2024.09.24.614704>