

Sep 28, 2023

Version 1

Sea Water - Blue Treasure V.1

DOI

<https://dx.doi.org/10.17504/protocols.io.36wgq36b5lk5/v1>



Willian Barela Costa^{1,2}

¹FabLab Casa Firjan; ²@casafirjan



Willian Barela Costa

personal

Create & collaborate more with a free account

Edit and publish protocols, collaborate in communities, share insights through comments, and track progress with run records.

Create free account

OPEN  ACCESS



DOI: <https://dx.doi.org/10.17504/protocols.io.36wgq36b5lk5/v1>

Protocol Citation: Willian Barela Costa 2023. Sea Water - Blue Treasure. protocols.io

<https://dx.doi.org/10.17504/protocols.io.36wgq36b5lk5/v1>

License: This is an open access protocol distributed under the terms of the [Creative Commons Attribution License](#), which permits unrestricted use, distribution, and reproduction in any medium, provided the original author and source are credited



Protocol status: Working

We use this protocol and it's working

Created: September 27, 2023

Last Modified: September 28, 2023

Protocol Integer ID: 88418

Keywords: blue treasure sea water production for microalgae cultivation, blue treasure sea water production, microalgae cultivation, blue treasure supplier, salinity, sea water

Abstract

Sea water production for microalgae cultivation using the Blue Treasure supplier.

25 - 35 Salinity

[M] 27 mg/mL

Materials

Salt from Blue Treasure product

Distilled water



Salting process - 1 liter

- 1 [M] 0.027 g/mL → 🔬 Sample 27 g Sea Sal to 🧴 1 L Water distilled
- 2 Shake until diluted
- 3 Make sure the salinity is within the desired range using the salinity refractometer. In this protocol, the focus is 27 salinity.