

Jan 27, 2026

SEAP reporter assay for assessing NF- κ B inhibition by TGF- β 1/PD-L1 fusion protein

DOI

<https://dx.doi.org/10.17504/protocols.io.bp2l6e38rgqe/v1>

Marvin De los Santos¹

¹ChordexBio



Marvin De los Santos

ChordexBio

Create & collaborate more with a free account

Edit and publish protocols, collaborate in communities, share insights through comments, and track progress with run records.

Create free account

OPEN  ACCESS



DOI: <https://dx.doi.org/10.17504/protocols.io.bp2l6e38rgqe/v1>

Protocol Citation: Marvin De los Santos 2026. SEAP reporter assay for assessing NF- κ B inhibition by TGF- β 1/PD-L1 fusion protein. **protocols.io** <https://dx.doi.org/10.17504/protocols.io.bp2l6e38rgqe/v1>

Manuscript citation:

De Los Santos, M. I., & Bernal, S. D. (2024). U.S. Patent No. 11,912,751. Washington, DC: U.S. Patent and Trademark Office.

De Los Santos, Marvin I. (2019). Construction and Characterization of a Novel Bifunctional TGF- β 1/PD-L1 Fusion Protein (M.S. in Molecular Biology and Biotechnology Thesis). University of the Philippines Diliman. College of Science. National Institute of Molecular Biology and Biotechnology. UP Diliman CS Library. <https://cslib.science.upd.edu.ph/>

License: This is an open access protocol distributed under the terms of the [Creative Commons Attribution License](#), which permits unrestricted use, distribution, and reproduction in any medium, provided the original author and source are credited

Protocol status: Working

We use this protocol and it's working

Created: January 26, 2026

Last Modified: January 27, 2026

Protocol Integer ID: 241572

Keywords: NF- κ B signaling, SEAP reporter assay, secreted alkaline phosphatase, secreted embryonic alkaline phosphatase, HEK-Blue TLR2 cells, PD-1 receptor, TGF- β 1/PD-L1 fusion protein, Immunomodulation, Inflammatory signaling, Reporter gene assay, κ b inhibition by tgf, κ b activation, κ b pathway inhibition, κ b inhibition, κ b, I1 fusion protein this protocol, I1 fusion protein, seap reporter assay, tlr2 reporter hek 293 cell, human tlr2 reporter hek 293 cell, immunomodulatory protein, signaling pathway, treatment with the tgf, dependent inhibition of nf, seap secretion into the culture medium, tgf

Abstract

This protocol describes a secreted embryonic alkaline phosphatase (SEAP)-based reporter assay to evaluate NF- κ B pathway inhibition mediated by a TGF- β 1/PD-L1 fusion protein. Human TLR2 Reporter HEK 293 cells are stably modified to express the PD-1 receptor and used as a cellular model to assess downstream NF- κ B activation following Toll-like receptor 2 stimulation. NF- κ B signaling is induced using graded dilutions of Mycobacteria-containing Freund's Complete Adjuvant, and pathway activity is quantified via SEAP secretion into the culture medium. Treatment with the TGF- β 1/PD-L1 fusion protein is compared against antibody-blocked and untreated controls to determine receptor-dependent inhibition of NF- κ B activation. This assay provides a reproducible and modular framework for mechanistic validation of immunomodulatory proteins affecting inflammatory signaling pathways.

Protocol materials

 Imject[®]; Freund's Complete Adjuvant (FCA) **Thermo Fisher Catalog #77140**

 HEK Blue Detection Medium **InvivoGen**

Safety warnings

- ! This assay is sensitive to the presence of endotoxins. Before performing the experiment, ensure first that your culture is not contaminated and your reagents are endotoxin-free.

Before start

This protocol uses Human TLR2 Reporter HEK 293 Cells from InvivoGen, lipofected stably with pCMV3-PD1 to express the PD-1 receptor. Please contact the author regarding inquiries on how to reconstitute the PD-1 protein in these cells or different eukaryotic cell models.



Confirmation of receptor expression in PD1⁺ TLR2⁺ HEK 293 cells

44m

- 1 Perform immunofluorescence (IF) analysis using anti-TGFBR1 and anti-PD1 antibodies. Note that you can use FITC-conjugated antibodies (direct) or premix the primary antibodies with FITC-conjugated secondary antibodies in 1:1 ratio (indirect)

- 1.1 Harvest 1×10^6 cells and collect by centrifugation  500 x g, 25°C, 00:05:00

5m

- 1.2 Resuspend the cell pellet in  100 μ L of PBS, repeat  [go to step #1.1](#) once

- 1.3 Add  100 μ L of staining buffer and resuspend the cells well by pipetting

A	B
Reagent	Amount needed
Bovine serum albumin (BSA)	0.05% (w/v)
FITC-labelled antibody	1:100 (antibody:buffer ratio)
PBS	Fill to 100 μ L

Composition of staining buffer

- 1.4 Incubated at  4 °C for  00:30:00 in the dark with constant shaking. Note that over-exposure to light may photobleach FITC signal

30m

- 1.5 Wash the cells in  5 mL PBS-T, twice.

A	B
Reagent	Amount needed
Tween-20	0.01% (w/v)
PBS	5 mL

Composition of PBS-T buffer

- 1.6 Collect the cells by centrifugation  500 x g, 25°C, 00:03:00 and resuspend the cells in  25 µL of PBS with or without 0.5 µg/mL of Hoechst 33342. Incubate for  00:05:00 at room temperature
- 1.7 Wash the cells once with  5 mL PBS and resuspend to a final volume  100 µL of PBS
- 1.8 Aliquot desired amount and dispense to poly-L-lysine coated slides, incubate in the dark for at least  00:01:00 prior to visualization under a fluorescence microscope

8m

1m

SEAP reporter assay

2d

- 2 Seed 1×10^5 cells in a 96-well plate, each well should not exceed  200 µL
- 3 After  24:00:00, replace the medium with same volume of  HEK Blue Detection Medium **InvivoGen**
- 4 Challenge the cells with an increasing dilution of *Mycobacteria*-containing  Imject™ Freund's Complete Adjuvant (FCA) **Thermo Fisher Catalog #77140**

1d

	A	B
	Set-up	Dilution
	1	1:100
	2	1:1,000
	3	1:10,000
	4	1:50,000
	5	1:100,000
	Control	None (PBS only)

Dilution reactions for the experiment



Note: Consider the number of replicates in your assays

- 5 Divide the set-up into fusion protein-treated, antibody-blocked, and untreated reactions. For this protocol, the inhibition of NF- κ B activation by the TGF- β 1/PD-L1 (TP) fusion protein is determined. Note that you can use a different immunomodulating protein or molecule for this assay.

	A	B
	Reaction	Purpose
	TP-treated	Assess the direct effect of receptor binding of TP with receptors in NF- κ B activation
	Antibody-blocked	Identify the consequences of blocking the receptors of TP in NF- κ B activation
	Untreated	Serves as control

SEAP reporter assay set-up; TP = TGF- β 1/PD-L1 fusion protein

- 5.1 For TP-treated cells, add 50 ng/mL of fusion protein to the wells
- 5.2 For antibody-blocked cell, add 1 ug/mL of either anti-TGFB1 and anti-PD1 antibodies, incubate for  00:15:00 , then add 50 ng/mL of fusion protein to the wells
- 5.3 For untreated cells, add the same volume of PBS to the wells
- 5.4 Culture the cells as usual for  24:00:00
- 6 Check the culture plate for color development, image the plate and then read the optical density (OD) at 620-655 nm on a plate reader

1d

