

Mar 27, 2019 Version 3

Selection of stable transformants in *Ostreococcus tauri*, *Ostreococcus lucimarinus* and *Bathycoccus prasinos* V.3

 [Nature Methods](#)

DOI

dx.doi.org/10.17504/protocols.io.zj2f4qe

Francois-Yves bouget¹

¹Laboratoire d'Océanographie Microbienne. Banyuls sur Mer

Protist Research to Opti...



Francois Yves Bouget

Laboratoire d'Océanographie Microbienne

OPEN  ACCESS



DOI: dx.doi.org/10.17504/protocols.io.zj2f4qe

External link: <https://doi.org/10.1038/s41592-020-0796-x>

Protocol Citation: Francois-Yves bouget 2019. Selection of stable transformants in *Ostreococcus tauri*, *Ostreococcus lucimarinus* and *Bathycoccus prasinos*. **protocols.io** <https://dx.doi.org/10.17504/protocols.io.zj2f4qe>

Manuscript citation:

Faktorová D, Nisbet RER, Robledo JAF, Casacuberta E, Sudek L, Allen AE, Ares M, Aresté C, Balestreri C, Barbrook AC, Beardslee P, Bender S, Booth DS, Bouget F, Bowler C, Breglia SA, Brownlee C, Burger G, Cerutti H, Cesaroni R, Chiurillo MA, Clemente T, Coles DB, Collier JL, Cooney EC, Coyne K, Docampo R, Dupont CL, Edgcomb V, Einarsson E, Elustondo PA, Federici F, Freire-Beneitez V, Freyria NJ, Fukuda K, García PA, Girguis PR, Gomaa F, Gornik SG, Guo J, Hampl V, Hanawa Y, Haro-Contreras ER, Hehenberger E, Highfield A, Hirakawa Y, Hopes A, Howe CJ, Hu I, Ibañez J, Irwin NAT, Ishii Y, Janowicz NE, Jones AC, Kachale A, Fujimura-Kamada K, Kaur B, Kaye JZ, Kazana E, Keeling PJ, King N, Klobutcher LA, Lander N, Lassadi I, Li Z, Lin S, Lozano J, Luan F, Maruyama S, Matute T, Miceli C, Minagawa J, Moosburner M, Najle SR, Nanjappa D, Nimmo IC, Noble L, Vanclová AMGN, Nowacki M, Nuñez I, Pain A, Piersanti A, Pucciarelli S, Pyrih J, Rest JS, Rius M, Robertson D, Ruaud A, Ruiz-Trillo I, Sigg MA, Silver PA, Slamovits CH, Smith GJ, Sprecher BN, Stern R, Swart EC, Tsaousis AD, Tsylin L, Turkewitz A, Turnšek J, Valach M, Vergé V, Dassow Pv, Haar Tvd, Waller RF, Wang L, Wen X, Wheeler G, Woods A, Zhang H, Mock T, Worden AZ, Lukeš J, Genetic tool development in marine protists: emerging model organisms for experimental cell biology. *Nature Methods* 17(5). doi: [10.1038/s41592-020-0796-x](https://doi.org/10.1038/s41592-020-0796-x)



License: This is an open access protocol distributed under the terms of the **Creative Commons Attribution License**, which permits unrestricted use, distribution, and reproduction in any medium, provided the original author and source are credited

Protocol status: Working

We use this protocol and it's working

Created: March 27, 2019

Last Modified: March 27, 2019


Protocol Integer ID: 21850

Abstract

This protocol describes the selection and growth of stable transformants in semi solid agarose medium. Developed initially for *Ostreococcus tauri*, it also works for *Ostreococcus lucimarinus* and *Bathycoccus*.

Inclusion of cells in semi-solid agarose medium

- 1 Autoclave a solution of 2.1% low melting agarose in H₂O. Keep at 65-90 °C in a water bath.
- 2 For each transformation prepare 8 Petri dishes (55 mm diameter) and 8 × 15 ml tubes each containing 9 ml of ASW plus the required selection. (G418 at 1 mg/ml).
- 3 Add 1 ml of LMP agarose to the 9 ml in one of the tubes. Close the tube, and mix gently by inverting.
- 4 Add 0.5 ml of overnight transformed cells (see protocol on transient transformation), quickly mix and gently pour into the plate. pay attention to avoid bubbles . Repeat this process for all tubes.
- 5 Let the plates dry open in the **flowhood** for about one hour, so that the agarose solidifies.

 01:00:00
- 6 Close the plates and transfer them to large square Petri dishes. Add wet paper towel to keep the chamber humid. Seal the square plates with medical tape. Place the square plates in the incubator for about 10 days.

Selection of transformants

- 7 Transformant colonies should appear after 10 to 21 days. Pick colonies using with cut-off yellow tips. Suck out the green colony. Take care not to include any cells from neighbouring colonies.
- 8 Transfer the cells to 0,2 ml of ASW medium containing the selection (G418), in 96 wells microplate. Allow cultures to grow for 7 days in the culture incubator.
- 9 After one week transfer to 20 ml culture flasks and grow for 7 to 10 days. Stable integration into the genome by random insertion or homologous recombination can be detected using **PCR (see Lozano et al, Plant Journal 2014)**. When the transgene contains a luciferase reporter, a first screening can be performed directly by measuring luminescence in a microplate luminometer.
- 10