

Nov 23, 2022

Version 2

Senescence Cocktail V.2

DOI

dx.doi.org/10.17504/protocols.io.4r3l24xmxg1y/v2

Sakthikumar Mathivanan¹

¹University of Wisconsin System

Daniel's workspace



Daniel El Kodsí

Brigham and Women's Hospital and Harvard Medical School

Create & collaborate more with a free account

Edit and publish protocols, collaborate in communities, share insights through comments, and track progress with run records.

Create free account

OPEN  ACCESS



DOI: <https://dx.doi.org/10.17504/protocols.io.4r3l24xmxg1y/v2>

Document Citation: Sakthikumar Mathivanan 2022. Senescence Cocktail. **protocols.io**
<https://dx.doi.org/10.17504/protocols.io.4r3l24xmxg1y/v2> Version created by **Daniel El Kodsí**

License: This is an open access document distributed under the terms of the **Creative Commons Attribution License**, which permits unrestricted use, distribution, and reproduction in any medium, provided the original author and source are credited

Created: November 23, 2022



Last Modified: May 31, 2024

Document Integer ID: 73188

Keywords: ASAPCRN, senescence in fibroblast, senescence of ipsc, neurodegenerative disorders with human stem cell, aged fibroblast, senescence cocktail, features characteristic of aged fibroblast, senescence, axonal degeneration, related neurodegenerative disorder, phenotypes in neuron, human stem cell, derived neuron, stem cell, als patient ipsc, embryonic fibroblast, fibroblast, motor neuron, neuron, generation of reliable drug discovery platform

Abstract

Modeling age-related neurodegenerative disorders with human stem cells is difficult due to the embryonic nature of stem cell derived neurons. We developed a chemical cocktail to induce senescence of iPSC-derived neurons to address this challenge. We first screened small molecules that induce embryonic fibroblasts to exhibit features characteristic of aged fibroblasts. We then optimized a cocktail of small molecules that induced senescence in fibroblasts and cortical neurons without causing DNA damage. The utility of the "senescence cocktail" was validated in motor neurons derived from ALS patient iPSCs which exhibited protein aggregation and axonal degeneration substantially earlier than those without cocktail treatment. Our "senescence cocktail" will likely enhance the manifestation of disease-related phenotypes in neurons derived from iPSCs, enabling the generation of reliable drug discovery platforms.

Troubleshooting

SENESCENCE COCKTAIL- ZHANG LAB


Our senescence cocktail (SLO) is a combination of **SBI-0206965** (10mM) + **Lopinavir** (1mM) + **O151** (1mM).

Cell types tested: In our experiments, we tested the activity of SLO on Fibroblasts, H9 derived cortical neurons, hiPSC-TDP43 (Mutant and Isogenic) derived motor neurons.

Final timepoint tested: Neurons were treated with SLO on Day 7 post plating and performed analysis on Day 12.

Final Concentration: 1:1000 dilution (add the senescence cocktail directly to the fresh neuronal media and then feed to the cells).

Workflow:

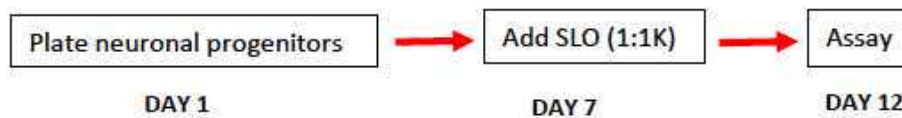

Workflow

User Optimization needed:

- 1) The user must optimize the concentration and timing for their experiments.
- 2) This depends on the cell type (neurons/ astrocytes) and timepoint (short term or long term).
- 3) We tested SLO on neuronal culture alone. If you are using neuron astrocyte co-culture, the concentration needs to be optimized.

Reagents:

- 1) SBI-0206965: Cat.no: S7885, Selleckchem
- 2) Lopinavir: Cat.no: S1380, Selleckchem
- 3) O151: Cat.no: 5153248, Hit2Lead (Chembridge)
- 4) Solvent- DMSO
- 5) Neuronal media- DMEM/ F12.

For more information please refer to our biorxiv preprint:

<https://www.biorxiv.org/content/10.1101/2021.07.11.451956v1>