

Jun 11, 2020

Version 2

Sentinel Flask Preparation for Cold Shipping Protocol of Human Islets V.2

DOI

<https://dx.doi.org/10.17504/protocols.io.bhdpj25n>



Integrated Islet Distribution Program¹

¹Integrated Islet Distribution Program, City of Hope

Integrated Islet Distribution Program

Tech. support email: iidp-email@coh.org



Integrated Islet Distribution Program

Integrated Islet Distribution Program, City of Hope

Create & collaborate more with a free account

Edit and publish protocols, collaborate in communities, share insights through comments, and track progress with run records.

Create free account

OPEN  ACCESS



DOI: <https://dx.doi.org/10.17504/protocols.io.bhdpj25n>

External link: <https://dx.doi.org/10.17504/protocols.io.bac5iay6>; <https://iidp.coh.org/Investigators/Policies-Standard-Operating-Procedures>

Protocol Citation: Integrated Islet Distribution Program 2020. Sentinel Flask Preparation for Cold Shipping Protocol of Human Islets. protocols.io <https://dx.doi.org/10.17504/protocols.io.bhdpj25n> Version created by Integrated Islet Distribution Program

License: This is an open access protocol distributed under the terms of the Creative Commons Attribution License, which permits unrestricted use, distribution, and reproduction in any medium, provided the original author and source are credited

Protocol status: Working

We use this protocol and it's working

Created: June 10, 2020

Last Modified: June 11, 2020

Protocol Integer ID: 38031

Keywords: Human Islets, culture, sentinel flask, IIDP, QA, Quality Assurance, Islet Equivalent, Prodo Labs, sentinel flask of human islet, research in the integrated islet distribution program, integrated islet distribution program, representative sample of the islet preparation, flask of islet, islet preparation, islet integrity at the production center, islet shipment, islet quality, sentinel flask preparation for cold shipping protocol, islet integrity, human islet, islet, sentinel flask preparation, cold shipping protocol, cold shipping condition, cold shipping, national institute of diabetes, cultured in the same shipping media, diabetes, flask,

Abstract

To establish a standardized method for the National Institute of Diabetes and Digestive and Kidney Diseases (NIDDK) sponsored research in the Integrated Islet Distribution Program (IIDP) for preparing a sentinel flask of human islets in order to monitor sterility and islet integrity at the production center. This is a representative sample of the islet preparation that has been distributed to investigators and is held at the distributing center as a quality control sample. The islets are cultured in the same shipping media used for islet shipment, for 18 hours at 4-8°C, simulating the cold shipping conditions followed by 2 days at 37°C culture. The flask of islets are verification in the event of a dispute over sterility or islet quality.

Note

Integrated Islet Distribution Program (IIDP) (RRID:SCR_014387)

Attachments



Prodo Media-Preparat...

441KB



GemCell Human Serum

...
81KB



Attachment 1-Solutio...

176KB



Attachment 2-Sentine...

209KB

Guidelines

- ***Integrated Islet Distribution Program (IIDP) (RRID:SCR_014387)***: The IIDP is a program commissioned and funded by the NIDDK to provide quality human islets to the diabetes research community to advance scientific discoveries and translational medicine. The IIDP consists of the NIDDK Project Scientist and Program Official, the External Evaluation Committee and the CC at City of Hope (COH). The IIDP CC integrates an interactive group of academic laboratories including the subcontracted IIDP centers.
- ***IIDP Coordinating Center (CC)***: Joyce Niland, Ph.D., IIDP Principal Investigator leads CC staff to coordinate the activities of the IIDP and assists the participating centers and investigators in the distribution of human islets.
- ***Islet Equivalent (IEQ)***: An IEQ is defined as an islet with a diameter of 150 microns. The number of IEQ in each size class is calculated by multiplying the number of islets by a conversion factor for each micron size.
- ***Approved Investigators***: Researchers who have requested islets from the IIDP for basic science studies and whose research protocols have been reviewed and approved by the IIDP.
- ***Sentinel Flask***: A representative sample of the islet preparation that has been distributed to investigators that is held at the distributing center as a quality assurance sample. This sample should be held in the 30 mL shipping flask at the center and should contain at least 100 IEQ cultured in the same shipping media used for islet shipment. *One flask must be prepared for each purity batch of islets distributed.*

Materials

MATERIALS

- ⊗ Human AB Serum (ABS) HI **Gemini Bio-Products Catalog #100-512**; Heat Inactivated
- ⊗ PIM(G)[®] (5 mL Glutamine/Glutathione) **Prodo Laboratories, Inc Catalog #PIM(G)[®]**
- ⊗ PIM(T)[®] **Prodo Laboratories, Inc Catalog # PIM(T)[®]**
- ⊗ Corning[™] Ciprofloxacin Hydrochloride **Fisher Scientific Catalog #MT61277RG** (Corning[™] 61277RG)

Equipment	
Invitrogen [™] EVOS [™] XL Core Imaging System or equivalent	NAME
Inverted Microscope	TYPE
Light Microscope	BRAND
AMEX1000	SKU
https://www.fishersci.com/shop/products/evos-xl-core-imaging-system/12562751#?keyword=Inverted+Microscopes	LINK

Equipment

Corning 4492 or equivalent

NAME

Sterile Polystyrene Serological Pipets

TYPE

Corning™ Stripette™ Wide-Tip Disposable Polystyren

BRAND

07-200-619

SKU

[https://www.fishersci.com/shop/products/stripette-disposable-plastic-wrapped-polystyrene-serological-pipettes-2/07200619?](https://www.fishersci.com/shop/products/stripette-disposable-plastic-wrapped-polystyrene-serological-pipettes-2/07200619?searchHijack=true&searchTerm=07200619&searchType=RAPID&matchedCatNo=07200619)

LI
N
K

[searchHijack=true&searchTerm=07200619&searchType=RAPID&matchedCatNo=07200619](https://www.fishersci.com/shop/products/stripette-disposable-plastic-wrapped-polystyrene-serological-pipettes-2/07200619?searchHijack=true&searchTerm=07200619&searchType=RAPID&matchedCatNo=07200619)

Sterile, individually wrapped, and certified nonpyrogenic and DNase-/RNase-free
Accuracy within $\pm 2\%$ at full volume

SPECIFICATIONS



Equipment

Gilson™ PIPETMAN Classic™ Pipets- F123601 or equivalent

NAME

Adjustable pipettor -P200

TYPE

Gilson

BRAND

F123601G

SKU

<https://www.fishersci.com/shop/products/gilson-pipetman-classic-pipets-model-p200-vol-50-200-l/f123601g>

LI
N
K

50 to 200 μ L, ± 0.5 , $\pm 1\mu$ L

SPECIFICATIONS



Equipment

Fisherbrand™ Large-Orifice Pipet Tips, 1 to 200µL or equivalent	NAME
Genomic/Wide Orifice Pipet Tips	TYPE
Fisherbrand	BRAND
02-707-134	SKU
https://www.fishersci.com/shop/products/fisherbrand-large-orifice-pipet-tips-6/02707134	LINK
1 to 200µL	SPECIFICATIONS



Equipment	
Drummond™ Fixed-Volume Microdispensers or equivalent	NAME
Drummond 3000385	TYPE
Drummond™	BRAND
21176F	SKU
https://www.fishersci.com/shop/products/drummond-fixed-volume-microdispensers-16/21176f?searchHijack=true&searchTerm=21176F&searchType=RAPID&matchedCatNo=21176F	LINK
Volumetric Range 100/200UL with borosilicate glass bores	SPECIFICATIONS
	

Equipment

Bottle Top Filters with PES Membrane or equivalent

NAME

0.2 µm Bottle Top Filter

TYPE

Thermo Scientific™ Nalgene™ Rapid-Flow™ Sterile Di

BRAND

09-741-09

SKU

<https://www.fishersci.com/shop/products/nalgene-rapid-flow-sterile-disposable-bottle-top-filters-pes-membrane/0974109>

LINK

0.2 µm membranes for sterile filtration; PES is the best membrane for cell culture fluids; lowest protein binding to maintain protein balance, lowest extractables to maintain media purity

SPECIFICATIONS



Equipment

Thermo Scientific™ Nalgene™ Square PETG Media Bottles with Closure	NAME
30 mL PETG Media Bottles with Closure	TYPE
Thermo Scientific™ Nalgene™	BRAND
03-311-1V	SKU
https://www.fishersci.com/shop/products/nalgene-square-petg-media-bottles-closure/033111V#?keyword=true	LINK

30 mL PETG Media Bottles

SPECIFICATIONS



Safety warnings

⚠ Please see attached SDS (Safety Data Sheet) for hazards and safety warnings.

Ciprofloxacin Hydrochloride

Precautionary statements:

- P280 - Wear protective gloves and eye/face protection
- P305 + P351 + P338 - IF IN EYES: Rinse cautiously with water for several minutes. Remove contact lenses, if present and easy to do. Continue rinsing.
- P337 + P313 - If eye irritation persists: Get medical advice/attention.
- P273 - Avoid release to the environment.

GemCell™ U.S. Origin Human Serum AB

GemCell™ human serum AB is collected from healthy male donors of the AB serotype at FDA-licensed facilities in the United States.

Hazardous Components:

- Biohazard contains human source material. Handle as though capable of transmitting infectious agents.
- Toxicity: Not Established.

Target Organs/Systems: Product could possibly irritate the skin, eyes and respiratory system. Do not ingest this product.

Before start

Prodo Labs, Inc. Protocols and Website: <http://www.prodolabs.com>

Citation

Ricordi, C. (1992). Pancreatic Islet Cell Transplantation. Austin: R.G. Landes Company, 1992:137-138..

Citation

Ricordi, C. (Ed). (1995). Methods in Cell Transplantation. Austin: R.G. Landes Company, 1995: Section G..

SUPPLEMENTAL MATERIALS

1 The following supplies are necessary for the preparation of Sentinel Test Flasks for human islet distribution.

- Islet preparation for distribution
- Wide mouth pipettes and pipettor
- 30 mL shipping bottles - 1 per islet purity batch broadcasted.
- Shipping media that is to be used for distribution as described below.

MEDIA PREPARATION

2 Shipping media that is to be used for distribution is prepared as described below.

- The Prodo Labs PIM(T) should be stored between 2° and 8°C upon receipt but is stable at room temperature.
- The (heat inactivated) Gemini AB serum and the PIM(G) vials should be stored at -5° to -20°C.

3 The Ciprofloxacin powder can be stored on the shelf but filter sterilized suspension aliquots should be stored at -5° to -20°C.

Pre-Preparation of Ciprofloxacin Powder for Addition to Media

- Remove 0.5 gm (500mg) of ciprofloxacin hydrochloride from the bottle and QS to 50mL with distilled water. This will give a stock concentration of 10mg/ml.
- Mix with a stir bar and stirring plate until totally dissolved.
- Filter sterilize the solution using a 0.2µM filter.
- Aliquot into sterile tubes, 5mL samples, label, and freeze for later use.
- The expiration date of the solution is indicated on the Certificate of Analysis and/or the bottle. Document expiration date as date of CoA.
- Diluted solution is good for 1 year frozen (if less than CoA expiration date) and 1 month thawed.

4 Prepare 1 bottle of PIM(T) media prior to the collection by adding the following:

- Thaw and add 5 mL of PIM(G).
- Add 12.5 mL of AB serum (2.5%v/v).
- Add 0.5 mL of prepared sterile ciprofloxacin aliquot.
- Once all additives have been added to the bottle of PIM(T), it is now referred to as PIM(T) complete.

**Note**

Note: If a prepared media bottle is to be used from a previous isolation, it must have been filter sterilized at the end of the previous use. The media will expire within 30 days, once it has been fully supplemented.

SENTINEL FLASK PREPARATION

- 5 Assemble all items described in the *Material and Media Preparation Sections*.
- 6 At the time of preparation for islet distribution, the cultured islets are consolidated and counted. Calculate the volume of suspension that is needed from the islet preparation to provide at least 100 IEQs.

The formula will be:

$$\text{Total IEQ} / (\text{Total mL} * 1000 \mu\text{L}) = x \text{ IEQ} / \mu\text{L} \text{ then } 100 \text{ IEQ} / x \text{ IEQ} / \mu\text{L} = y \mu\text{L}$$

- 7 Thoroughly resuspend the preparation with a large mouth pipet and remove the calculated aliquot. Place in the 30mL shipping flask. Top with appropriate amount of media using the same batch of PIM(T) complete that will be used for distribution. *Repeat for each purity batch that is being distributed.*
- 8 Store this/these Sentinel QA Flask(s) between 4°C and 8°C for approximately 18 hours after which the flask(s) should be moved to a 37°C incubator for an additional 2 days.

Note

Note: This culture procedure should simulate the average shipping time followed by culture by an investigator. A lab refrigerator should be appropriate for the first 18 hours of cold temperature culture.

- 9 Observe under a microscope daily for visual islet integrity, quality, and sterility. Record results on *Attachment 2 - Sentinel Flask Documentation Sheet*.
- 10 Alert the IIDP and the recipients of the distributed islets if there is a problem with the sterility of the preparation. In which case, send a copy of the *Attachment 1-Sentinel Flask Documentation Sheet* to IIDP.





Note

The sentinel flask is mainly for the documentation of the preparation by the distribution center in circumstances where islet sterility, quality or integrity of the distributed islets is questioned by a recipient. Photographs may be required to document the visual state of the sentinel flask if there is a challenge.

Citations

Ricordi, C.. Pancreatic Islet Cell Transplantation.

Ricordi, C. (Ed).. Methods in Cell Transplantation.